# QuikSip™ BT-Aspirator

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## Operating Manual

Please read carefully before use, and follow all operating and safety instructions!

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## Safety Instructions

This aspirating instrument is not designed for universal use. This instrument may be involved with hazardous materials, operations, and equipment. This Manual does not purport to address all of the safety problems associated with its use. It is the responsibility of whomever uses this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

Please follow the safety instructions of the cell-culture™ Aspiration Controller.

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## Application

This instrument is designed for aspirating aqueous solutions.

Valve block (H), filling tube (J), and valves (G, J) withstand steam sterilization at 121 °C. The dispensing cartridge (E) and pump assembly (B) are not resistant to steam sterilization.

The cell-culture™ Aspiration Controller (K) is not resistant to steam sterilization. The included tube and the adapters are resistant to steam sterilization.

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## Operating range

Aqueous solutions of inorganic and organic substances such as biological buffers, liquids for cell culture, etc.

Please see our catalog for a listing of chemical compatibility with components of this instrument. Please ask for further information.

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## Operating Exclusions

Never use this instrument with liquids which attack PP, PE or EPDM. If used in combination with the cell-culture™ Aspiration Controller never use with liquids, which additionally attack silicone and PVC.

Liquids that should not be aspirated include, but are not limited to:

- non-polar solvents like hydrocarbons and halogenated hydrocarbons
- concentrated or oxidizing acids.

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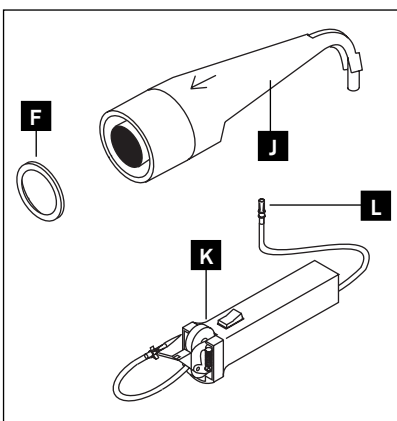
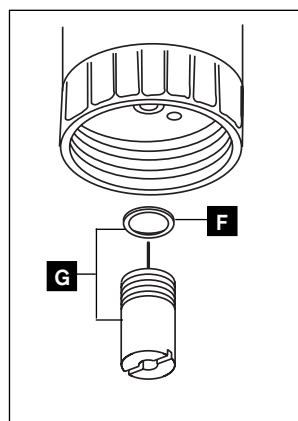
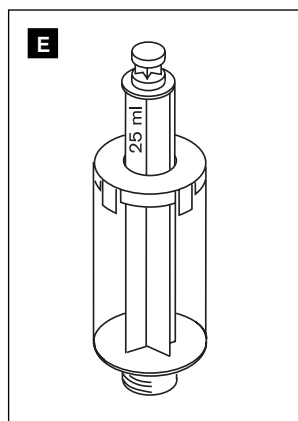
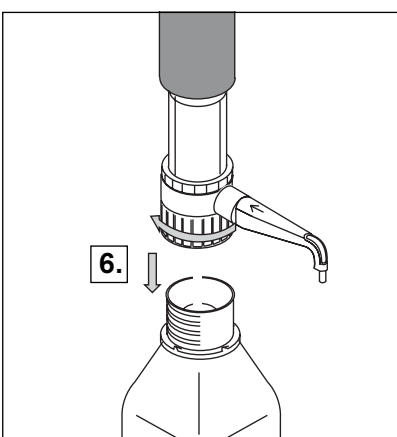
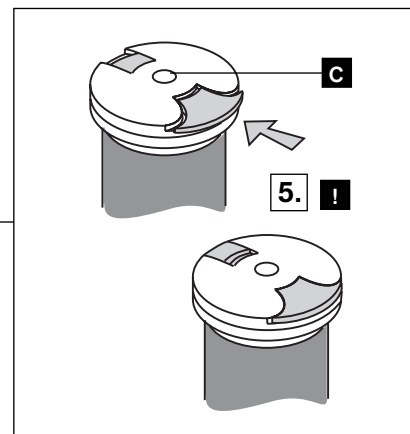
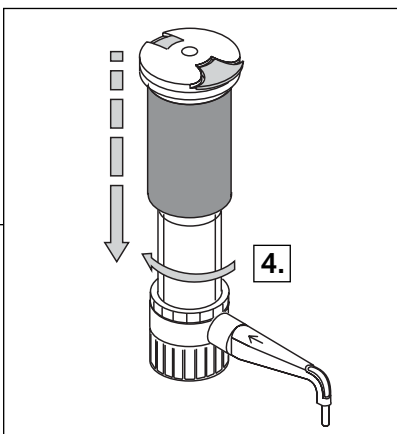
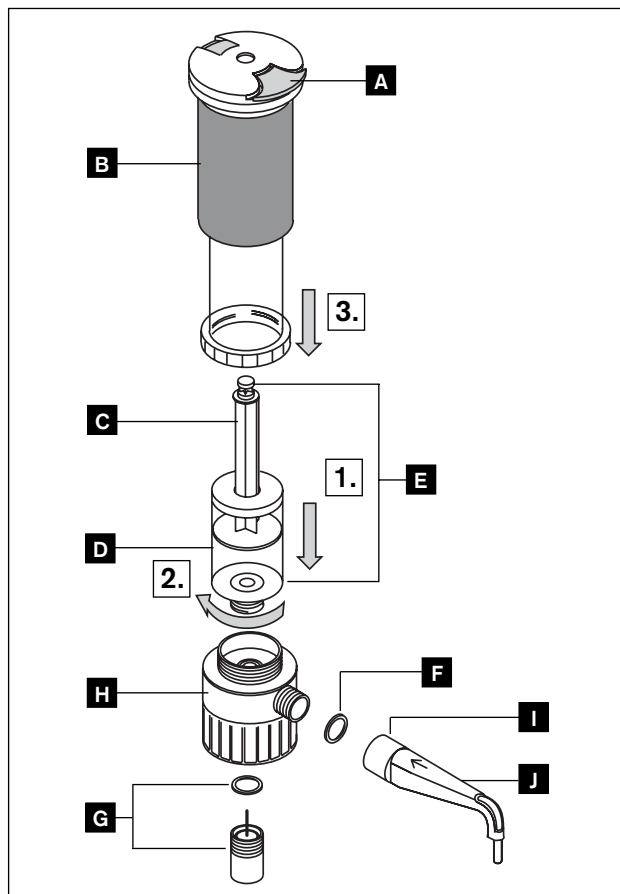
**Components**

- A Piston lock
- B Pump assembly, PC
- C\*** Piston, PE
- D\* Dispensing cylinder, PP
- E\* Dispensing cartridge, PP/PE
- F\* Seals, EPDM
- G\* Discharge valve, PP/EPDM
- H\* Valve block, PP
- I Locking nut, PP
- J\* Filling tube, PP, with filling valve, PP/EPDM
- K\* cell-culture™
- L\* adapter for cell-culture™, PP

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### Warning!

Wear protective clothing, goggles and gloves. Follow all Safety Instructions and observe the Operating Exclusions (⇒ pages 4-7).

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### Note:

Follow all instructions on page 8/9 for proper assembly.

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To control the vacuum, connect the cell-culture™ Aspiration Controller with the QuikSip™ BT-Aspirator using the included silicone tube. The link is effected by the included adapter.

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### Attention!

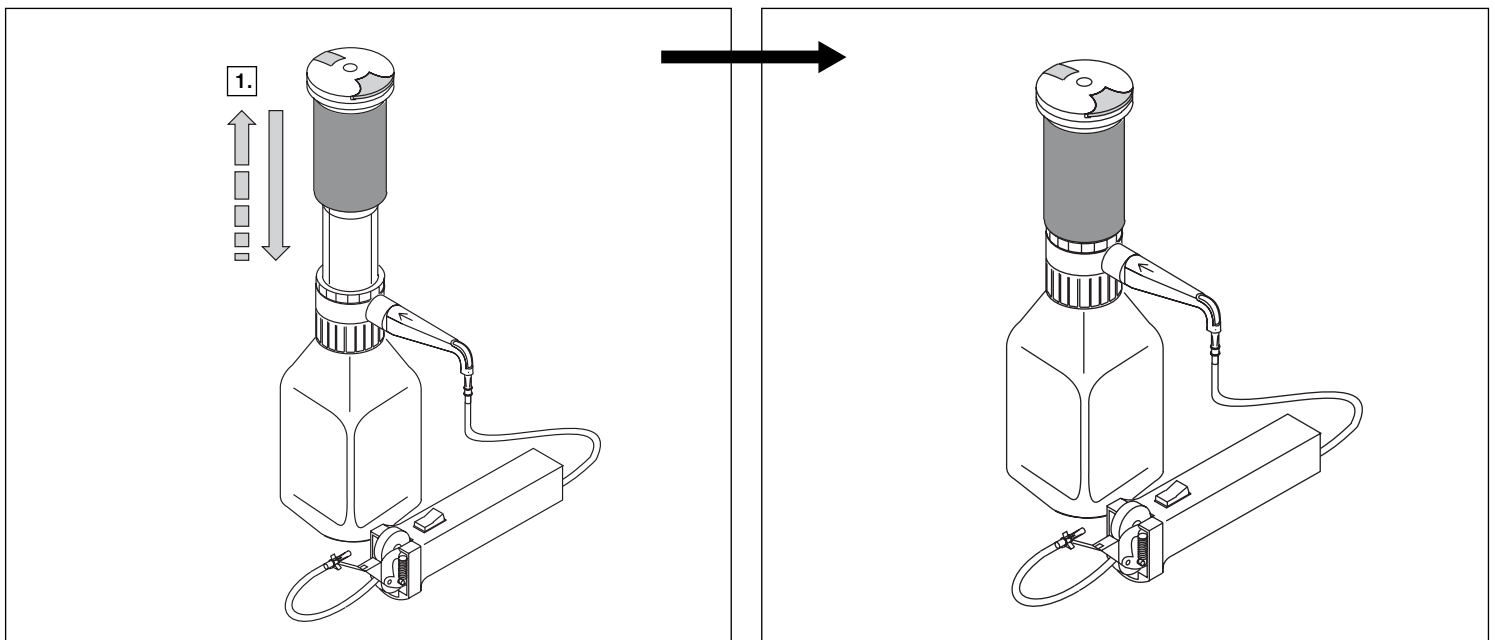
Use a bottle stand when using smaller or other potentially unstable bottles to help protect against tipping. Observe all Safety Instructions (⇒ p. 4-5).

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### Attention!

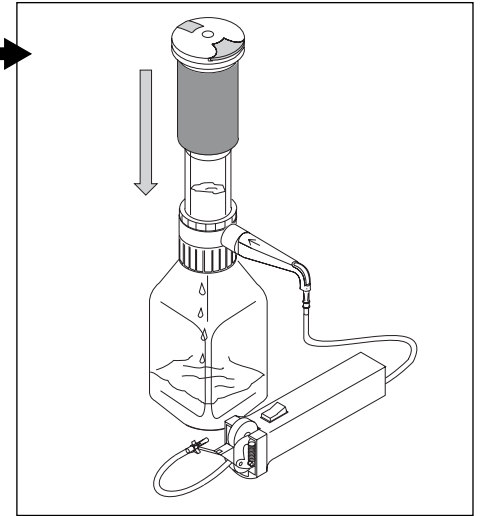
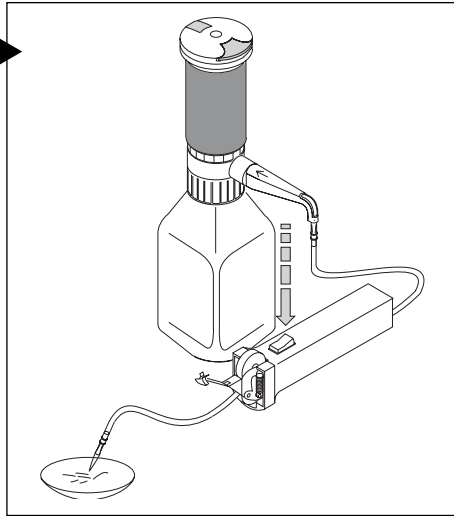
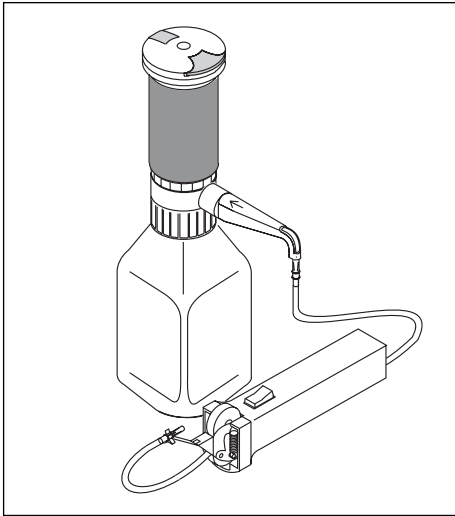
Follow all Safety Instructions and observe the Operating Exclusions (⇒ pages 4-7).

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Once the pump assembly has been pushed down, aspiration is accomplished by depressing the button on the cell-culture™. Aspiration rate is proportional to the pressure on the button. Releasing the button ceases aspiration.

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## Replacing the dispensing cartridge

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### Attention!

The dispensing cartridge is a wearing part. Replace the cartridge when the piston is difficult to move or is not airtight. To prevent chemical injury, the dispensing cartridge has to be first cleaned (⇒ page 15) then emptied prior to replacement.

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### Note:

Follow the diagrams on pages 8/9 for replacing the dispensing cartridge.

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## Cleaning

The instrument must be cleaned in the following situations to assure correct operation:

- immediately if the piston becomes difficult to move
- prior to long term storage
- before replacing the valves
- before replacing the dispensing unit

Depending upon the frequency of use, the dispensing cartridge should be replaced regularly.

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### Warning!

Follow all Safety Instructions (⇒ page 4-5).

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For cleaning the cell-culture™ Aspiration Controller please follow the instructions in the corresponding cell-culture™ manual.

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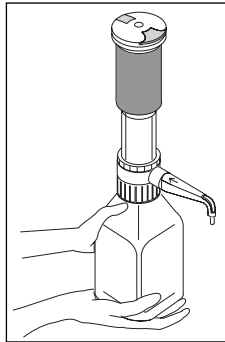
## Normal cleaning procedure

### Warning!

Carry the mounted instrument only as shown in the figure.

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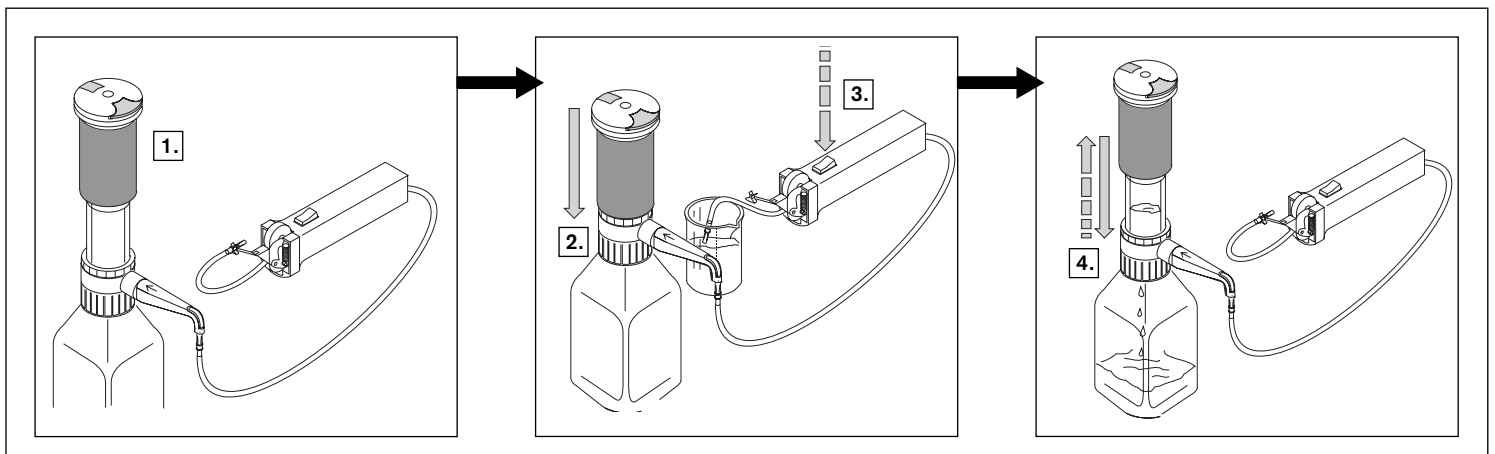
**B** suitable solvent, e.g., distilled water

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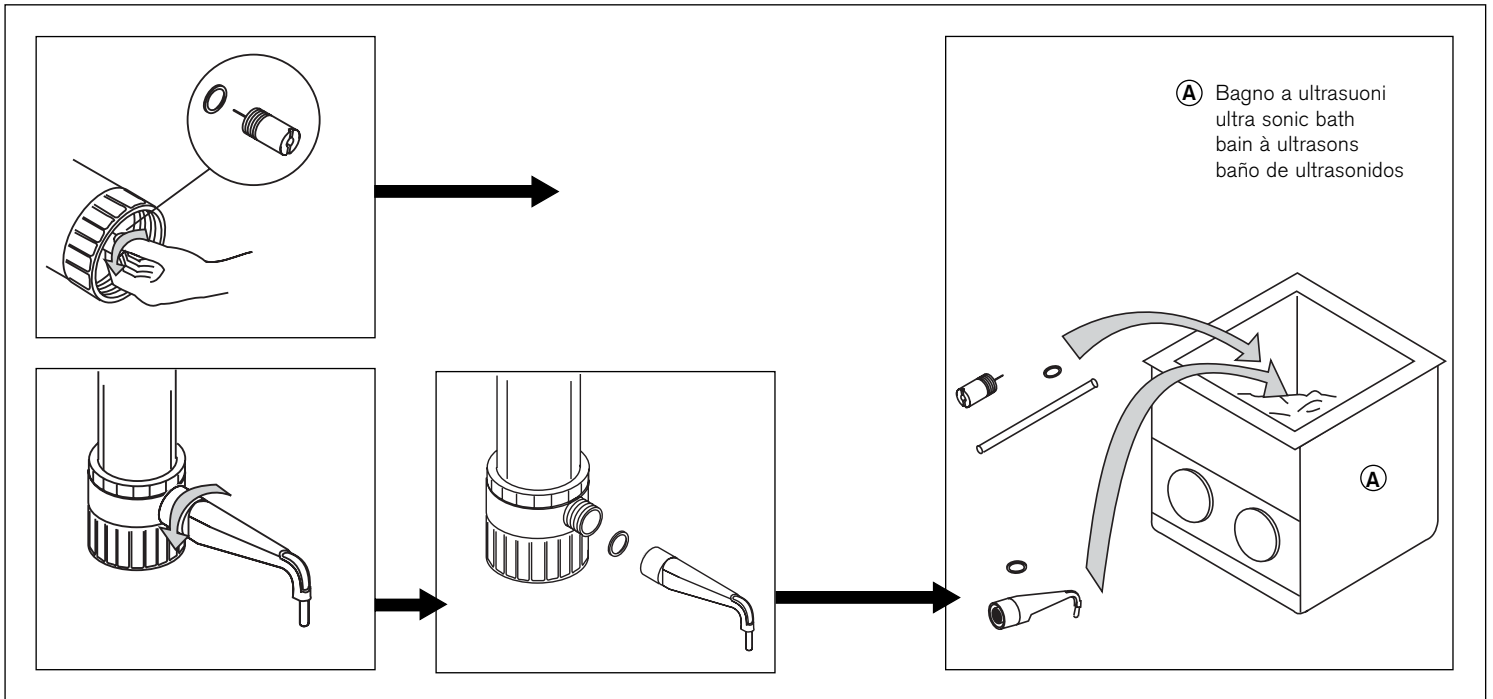
## Troubleshooting

Problem	Possible cause	Corrective action
<b>Will not aspirate</b>	<ul style="list-style-type: none"> <li>- Dispensing cartridge not locked</li> <li>- Dispensing cartridge leaks</li> <li>- Filling valve leaks or is sticking</li> <li>- Connection between filling tube and tube of the cell-culture™ leaks</li> <li>- Tube of cell-culture™ blocked or sticking</li> </ul>	<ul style="list-style-type: none"> <li>- Lock dispensing cartridge (⇒ page 8).</li> <li>- Mount new dispensing cartridge (⇒ page 13).</li> <li>- Clean filling valve (⇒ page 16) or replace valve and seal.</li> <li>- Stick adapter firmly onto the filling tube and pull tube firmly onto the adapter.</li> <li>- Clean tube or replace.</li> </ul>
<b>Piston and dispensing cartridge difficult to move</b>	<ul style="list-style-type: none"> <li>- Crystal deposition</li> <li>- Improper liquid aspirated</li> </ul>	<ul style="list-style-type: none"> <li>- Replace dispensing cartridge (⇒ page 13).</li> <li>- Check, if liquid is proper to aspirate.</li> </ul>
<b>Liquid behind the piston of the dispensing cartridge</b>	<ul style="list-style-type: none"> <li>- Dispensing cartridge leaks</li> </ul>	<ul style="list-style-type: none"> <li>- Replace dispensing cartridge (⇒ page 13).</li> </ul>
<b>Aspirated liquid can't be delivered into the bottle</b>	<ul style="list-style-type: none"> <li>- Discharge valve sticking or damaged</li> </ul>	<ul style="list-style-type: none"> <li>- Clean valve (⇒ page 16) or replace.</li> </ul>

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The liquid volume, theoretically to be aspirated with one stroke, is 25 ml. The tube length, the viscosity of the aspirated liquid and other losses however could result in reduction of the aspirated volume.

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**Attention!**

For safety reasons only clean/decontaminated instruments can be checked/repared.

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- Therefore: Clean and decontaminate the instrument carefully.
- Complete the "Declaration on Absence of Health Hazards" (ask your supplier or manufacturer for forms. The forms can also be downloaded on **www.brand.de**) and send the instrument to the manufacturer or supplier and describe:
  - the nature of the problem
  - the liquids disensed.

Return shipment has to be effected at the risk and cost of the sender.

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**Warranty**

We shall not be liable for the consequences of improper handling, use, servicing or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original parts have been used.

Subject to technical modification without notice. Errors excepted.

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