accu-jet[®]-Pipettierhelfer accu-jet[®] Pipette Controller

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The device named below fulfills the relevant fundamental requirements of the EC directives and standards listed. In the case of unauthorised modifications to the device, the declaration becomes invalid.

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Read carefully before use!

Safety Instructions

This instrument may be involved with hazardous materials. This Operating Manual does not purport to address all of the safety problems associated with its use. It is the responsibility of whoever uses this instrument to consult and establish appropriate safety and health practices and to determine the applicability of regulatory limitations prior to use.

1. Follow the general instructions for hazard prevention and safety regulations, e.g., wear protective clothing, eye protection and gloves.
2. Observe the reagent manufacturer's information.
3. Every user must read and understand this Operating Manual before using the instrument.
4. Never use the instrument in an atmosphere with danger of explosion.
5. Always work in a way which neither endangers the user, nor any other person.
6. Use the instrument only for pipetting, and within the recommended limits of its chemical resistance and mechanical properties.
7. Never use force when using this instrument.
8. In case of trouble, immediately stop pipetting. Clean and repair the instrument according to the

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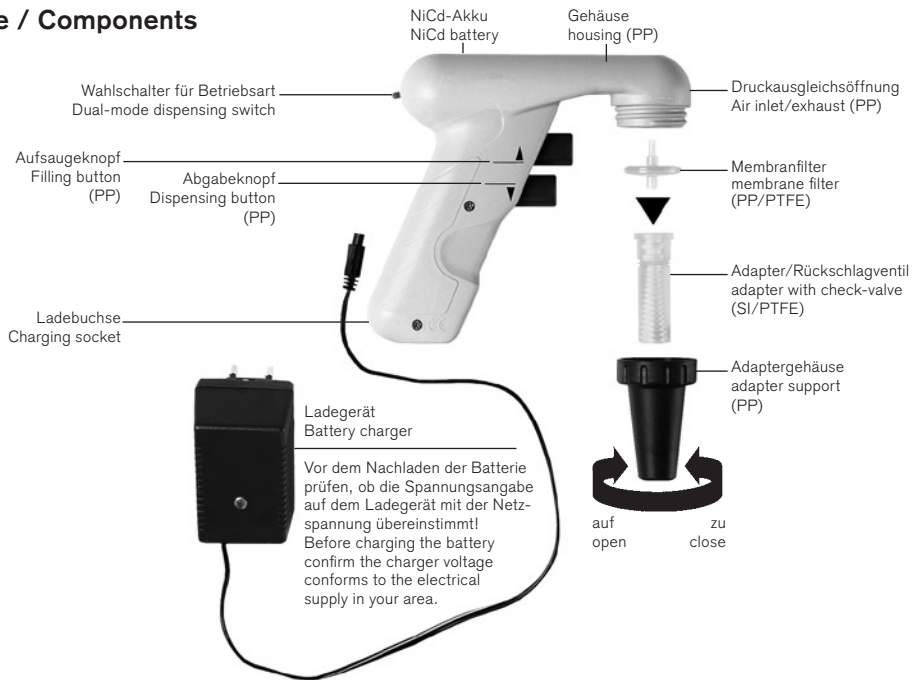
instructions before any further use of the instrument (➤ pages 13-18) or contact an authorized dealer or the manufacturer.

9. Only use original manufacturer's accessories and spare parts.
10. Only the original recharger can be used for recharging of the NiCd battery.
11. The recharger has to be protected against moisture and must only be used for this instrument.
12. Only authorized electricians are allowed to carry out repairs or interventions into the instrument.
13. Only open the instrument for replacing and disposal of the NiCd battery.
14. The built-in battery must not be replaced by another battery, especially not by primary cells. Only use original manufacturer's battery!
15. Dispose only completely discharged batteries. Please dispose the NiCd battery according to the relevant provisions!

Warning!

Improper use of the instrument or the battery (short-circuit, mechanical damage, overheating etc.) may cause the explosion of the battery.

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Storage conditions

Store instrument and accessories in a cool, dry place.

Storage temperature:

- 20 °C to + 50 °C

Starting operation

Before starting to operate, confirm the charger voltage conforms to the electrical supply in your area and recharge the battery, (> pages 16-17).

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Application

The instrument is designed to assist the filling and dispensing of graduated and volumetric pipettes of glass or plastic used for measuring liquids in the volume range of 0.1 mL to 100 mL.

It is not designed for Pasteur pipettes.

If the instrument is used correctly, the pipetted liquid will only contact the pipette.

Restrictions of use

Fumes of liquids come in contact with silicone, PTFE, PP, EPDM and PF (phenol-formaldehyde mass). Therefore the instrument must not be used with fuming liquids that are incompatible with materials above.

Warning!

Never use or recharge the instrument in an atmosphere with danger of explosion. Highly flammable liquids (e. g. ether, acetone and other liquids with a flash point below 0 °C) must not be pipetted.

Temperaturbeschränkungen

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+ 10 °C bis + 35 °C

Restrictions of temperature

Working- and recharging temperature:

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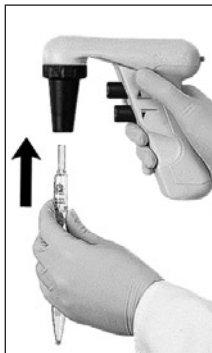
Pipetting

Attach the pipette

Hold the pipette as near to its upper end as possible, and carefully insert it into the adapter until it fits tightly.

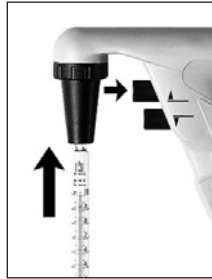
Attention!

Be sure that the pipette fits tightly into the adapter. Never use force. Thin pipettes are particularly liable to break. Avoid the risk of injury!

Once the pipette has been securely attached, always hold the instrument in a vertical position, tip down.

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Filling

Immerse the pipette tip into the liquid. Slowly press the upper pipetting button. The speed of aspiration is dependent on the amount the pipetting button is depressed. Fill the pipette so that the meniscus is slightly above the mark desired.

Attention!

Take care to avoid aspiration of liquid into the instrument, which can impair instrument performance. If liquid is aspirated, clean the pipette adapter/check valve, and replace the membrane filter, if necessary (> pages 13-16).

Adjust the volume

Use suitable lint-free tissue to wipe the pipette tip. Slowly press the lower button until the meniscus is adjusted exactly to the desired volume.

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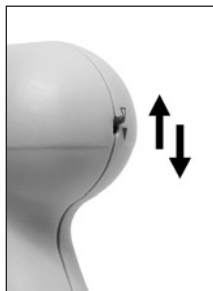
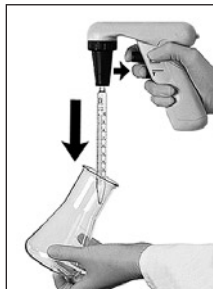
Blow-out and gravity-driven delivery

Slowly press the lower pipetting button. The harder the pipetting button is pressed, the faster the pipette is emptied.

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Storing the instrument

Attach the instrument to the wall support for storage.

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Note:

To mount the wall support remove the protective film from the snap tapes and tightly press the wall support on the desired place. The surface has to be smooth, clean and grease-free. Clean the surface with Isopropyl alcohol / deionized water (1:1). The snap tapes allow for easy detaching of the wall support for cleaning purposes. Alternatively the wall support can be fixed with two screws.

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Cleaning

Occasionally wipe the housing (PP) with a damp cloth. When properly used, the instrument is maintenance-free.

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Replacing the filter and cleaning the non-return valve

Attention!

Wear protective gloves and eye protection. Avoid the risk of contamination!

If the pipette has been overfilled or the suction capacity is impaired, proceed as described below:

1. Pull out the pipette.
2. Unscrew the adapter support.
3. Pull out the membrane filter.

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4. Remove the adapter from the adapter support by pushing it upwards.
Do not dismantle further!

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5. For cleaning the adapter with built-in non-return valve, carefully rinse it by using a wash bottle. Blow out the liquid and allow to dry completely.

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6. Place the new membrane filter with the thicker end facing downwards into the adapter (with the smaller conical side up).

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7. The entire pipette adapter incl. membrane filter is autoclavable at 121°C.

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Note:

Not every membrane filter is suitable. Only use genuine manufacturer's recommended accessories. Each instrument will be supplied with a 0.2 µm membrane filter (red color code). The membrane filter is autoclavable up to 5 times at 121 °C, 2 bar absolute (i.e. 1 bar above atmospheric pressure) for 20 minutes. It is the user's responsibility to ensure effective autoclaving.

The filter with 0.2 µm pore size (for cellculture) is sufficient to meet class B tolerances.

For class A accuracy it is necessary to use a 3 µm membrane filter. This filter has a larger pore size and thus less influence on the time of gravity-driven delivery.

(Spare parts ➤ page 27).

8. Reassemble the instrument in reverse order.
9. For leak test see next page (➤ page 16).

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Leak Test

When the meniscus has been set, liquid should not drip out of the pipette until the valves are activated. Should the pipette drip, disassemble the pipette adapter and reassemble it carefully.

Recharging the battery

One full charge of the NiCd battery allows 8 hours of non-stop pipetting. Decreasing motor power indicates the need for recharging.

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Warning!

Use original battery charger only!
Use of a different charger could damage the battery and would void the warranty. Confirm the voltage of the charger matches the voltage of your electrical outlet. The power supplied to the charger should differ by no more than $\pm 10\%$.

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Recharging procedure

1. **Temperature range for recharging:**
+ 10 °C to + 35 °C
2. A light-emitting diode (LED) will glow during recharging.
3. For the first six hours, the battery is charged in a high-current mode. Afterwards, the charger switches to a pulsed charge for long term battery life (LED will flash; 5 sec. on, 20 sec. off).
4. After 16 hours the battery will be fully charged.

Note:

- The instrument can be used during recharging.
- If the recharging procedure is not successful (> page 24).

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Replacing the battery

Note:

Only use original manufacturer's accessories
(> page 27). The battery must not be replaced
by primary cells or other rechargeable batteries.

Warning!

Use of the wrong batteries or improper use of the
instrument (e.g., short-circuit, mechanical damage,
overheating etc.) may cause the batteries to
explode.

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
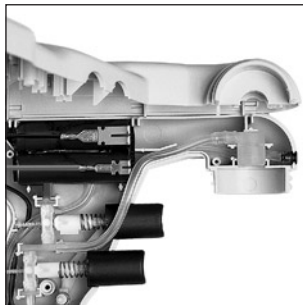
Pierce through the sealing wax and
remove the screws.

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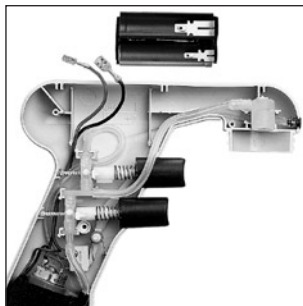
Pull off the pole plugs and remove the battery.

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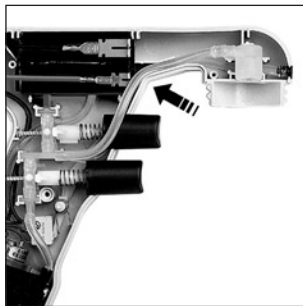
Insert the new battery and push on the pole plugs as far as they will go. Push the cables to the bottom of the housing.

Ensure correct position of the hose.

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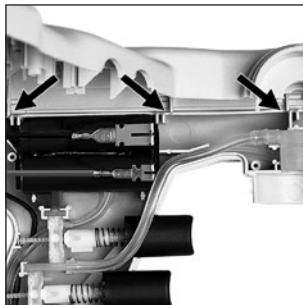Align the parts of the housing; watch for correct position of the three pins on the edge.

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When closing the two housing parts, push the hose inside, past the edges.

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To check if the motor switches on, operate both pipetting buttons once. Insert and tighten the screws and mount the adapter support.

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Disposal

Dispose NiCd-battery only completely discharged and according to the relevant provisions. For disposal of the instrument, loosen the screws, remove the batteries and dispose them separately.

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Trouble Shooting

Trouble	Possible Cause	Action to be taken
Suction capacity impaired	<ul style="list-style-type: none"> - Non-return valve closed - Filter or non-return valve dirty - Batteries discharged 	<ul style="list-style-type: none"> - Press the lower button down (blow-out) (> page 11) - Clean and/or replace the membrane filter (> pages 13-16) - Recharge batteries (> p. 16-17)
Filling only possible after fully pressing the lower button	<ul style="list-style-type: none"> - Filter or non-return valve dirty 	<ul style="list-style-type: none"> - Clean and/or replace the membrane filter (> p. 13-16)
Pipette drips	<ul style="list-style-type: none"> - Filter not properly positioned - Adapter or pipette damaged 	<ul style="list-style-type: none"> - Insert filter properly (> p. 14) - Replace adapter or pipette (> p. 13-16)
Pipette not held properly	<ul style="list-style-type: none"> - Adapter dirty or damaged 	<ul style="list-style-type: none"> - Clean the adapter, allow it to dry and/or replace it (> p. 13-16)
Recharging procedure not successful; LED-display at the charger doesn't light up	<ul style="list-style-type: none"> - Instrument or charger defective 	<ul style="list-style-type: none"> - Send the instrument in for repair or replace the charger (> p. 25)
Recharging procedure not successful; LED-display at the charger does light up	<ul style="list-style-type: none"> - Batteries defective 	<ul style="list-style-type: none"> - Replace the batteries (> p. 18-22)

If recommendations in the Trouble Shooting do not solve current problems, contact your authorized dealer or the manufacturer directly.

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Return for Repair

Attention:

For safety reasons only clean/decontaminated instruments can be checked/repared.

- Therefore: Clean and decontaminate the instrument carefully.
- Complete the “Declaration on Absence of Health Hazards” (ask your supplier or manufacturer for forms. The forms can also be downloaded on **www.brand.de**) and send the instrument to the manufacturer or supplier and describe:
 - the nature of the problem
 - the pipetted liquids.

Return shipment has to be effected at the risk and the cost of the sender.

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(www.brand.de/deutsch/agb/agb-d.pdf)

Warranty

As provided by law, any and all warranties are null and void if the product has been misused, modified or repaired by unauthorized personnel, if the defects are caused by negligence (instruction manual, maintenance) or by normal wear and tear. Use only original manufacturer's accessory/spare parts.

For conditions and extend of warranty refer to our General Conditions of Sale.

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Best.-Nr. 264 04

Ordering Data

accu-jet® pipette controller with charger,

incl. wall support and
2 spare membrane filters (0.2 µm)

	accu-jet® pipette controller			
	grey	blue	yellow	mint
with recharger for	Cat. No.	Cat. No.	Cat. No.	Cat. No.
Europe (continental) 230 V/50 Hz	264 00	264 30	264 31	264 32
UK/Ireland 230 V/50 Hz	264 01	264 40	264 41	264 42
USA 120 V/60 Hz	264 02	264 60	264 61	264 62
Australia 230 V/50 Hz	264 03	264 50	264 51	264 52
Japan 100 V/50 Hz	264 05	264 06	264 07	264 08

accu-jet® pipette controller without charger,

incl. wall support and
2 spare membrane filters (0.2 µm) **Cat. No. 264 04**

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NiCd-Akkublock **Best.-Nr. 265 31**

Spare Parts

**Silicone adapter
with non-return valve** **Cat. No. 265 08**

Adapter support **Cat. No. 265 12**

Spare membrane filter 0.2 µm, sterile
(red color code)
1 pc. in blister pack **Cat. No. 265 30**

Spare membrane filter 0.2 µm, non-sterile
10 pc. in PE-bag **Cat. No. 265 35**

Spare membrane filter 3 µm, non-sterile
1 pc. in blister pack **Cat. No. 260 52**

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This sign certifies that the product meets the requirements of the EEC directive and has been tested according to the specified test methods.