EASYCAL™ 5

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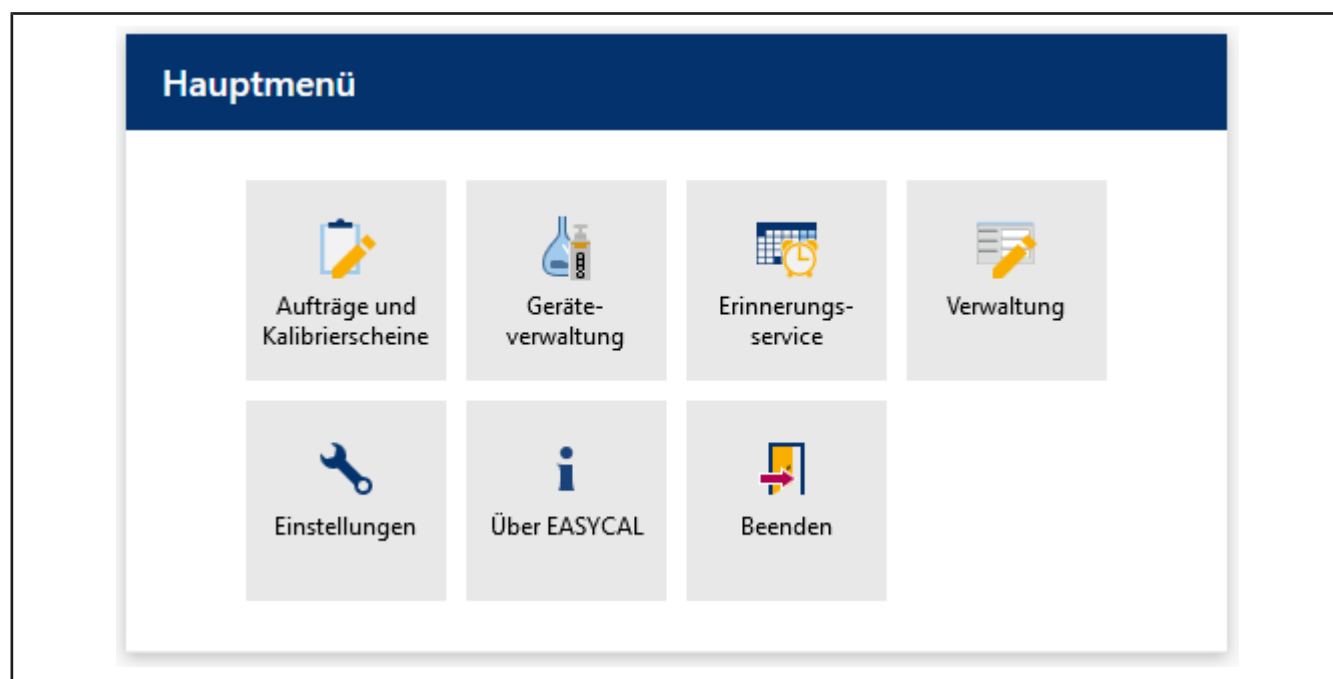
info@brand.de

www.brand.de

More operating manuals and translations?

Please refer to <http://www.brand.de/om> or use the following Quick Response Code:

The original operating manual is in German. Other languages are translations of the original operating manual.

Subject to technical changes, errors, and misprints.

Contents

- 1 Preliminary remarks 35**
- 2 Installation..... 36**
- 3 First steps in EASYCAL™ 5..... 45**
 - 3.1 Tutorial..... 45
 - 3.1.1 Set up balance as test equipment..... 45
 - 3.1.2 Creating order 45
 - 3.1.3 Determine test equipment 50
 - 3.1.4 Performing the test 50
 - 3.1.5 Complete test..... 58

1 Preliminary remarks

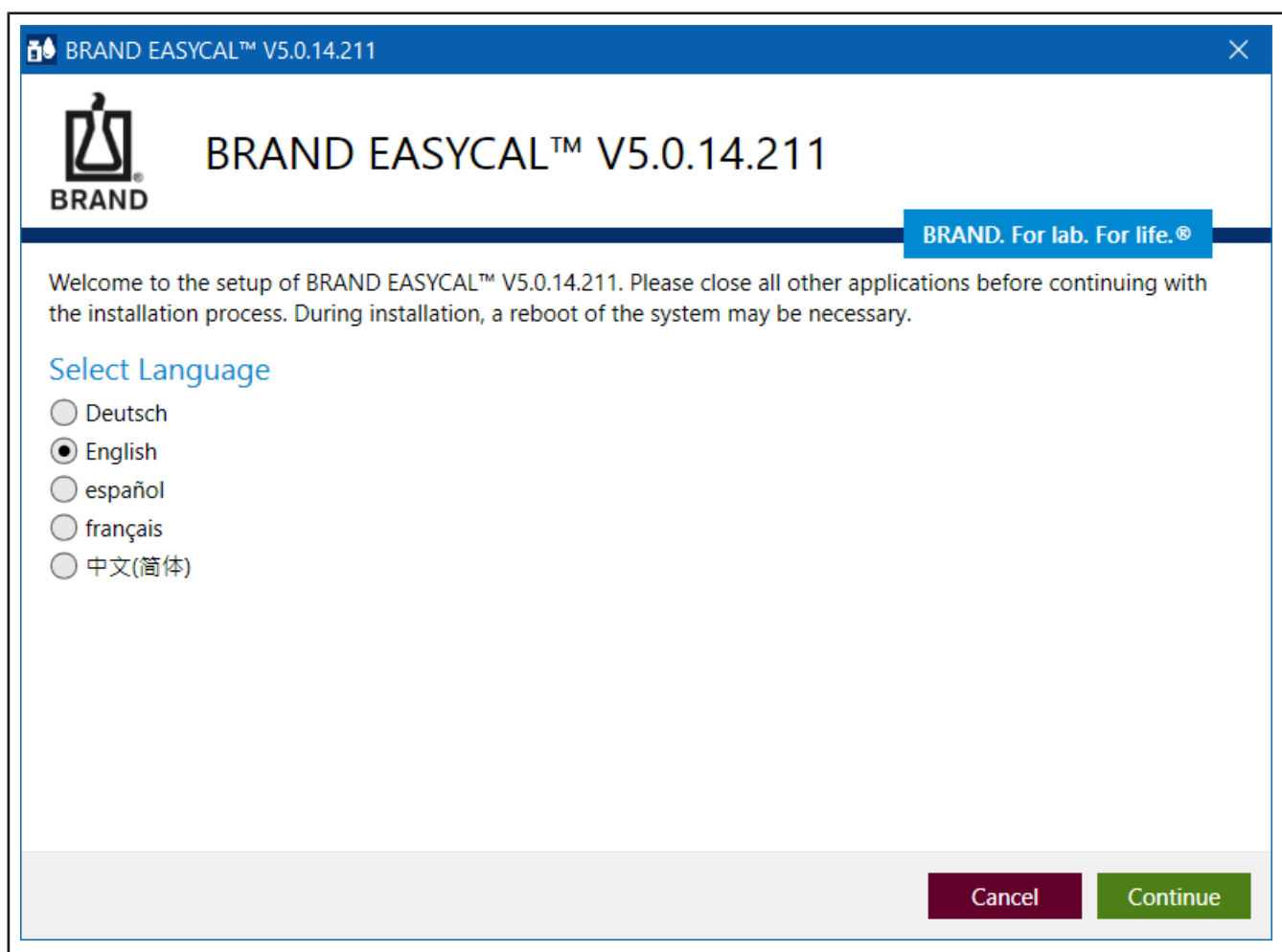
This tutorial is intended to make it easier for you to use EASYCAL™ 5 as part of the 30-day test period. Using an exemplary process, you will learn how to fill different areas of EASYCAL™ 5, how to carry out the first device test and how to print out a certificate.

Please note that in the course of this introduction, sample data will be generated in the database, which cannot be deleted later. We recommend coordinating this with the regulatory requirements of your quality management system.

2 Installation


Prerequisite:

- ✓ You have already received the EASYCAL™ 5 installation file, EasyCal-Setup.exe, from the BRAND home-page and saved it locally.
 - ✓ You have also received a test license (ticket) by E-mail.
 - ✓ You have local admin rights and an existing internet connection.
- a. Start the installation by opening the EasyCal-Setup.exe and follow the instructions in the installation wizard.


➔ EASYCAL™ 5 and all software components required for operation are installed on your PC.

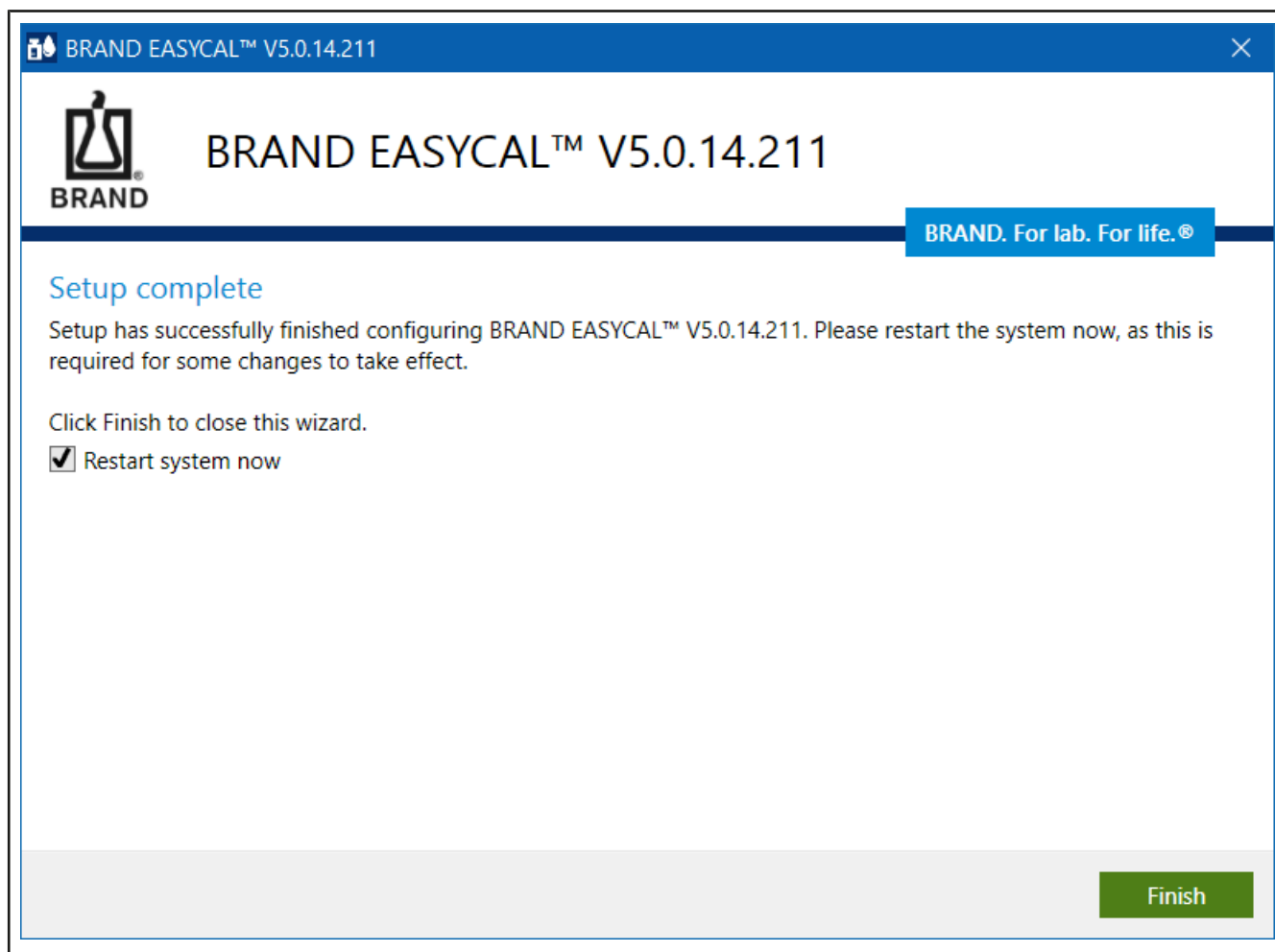- b. After completing the installation, start EASYCAL™ 5 by ticking the box and clicking on Complete.

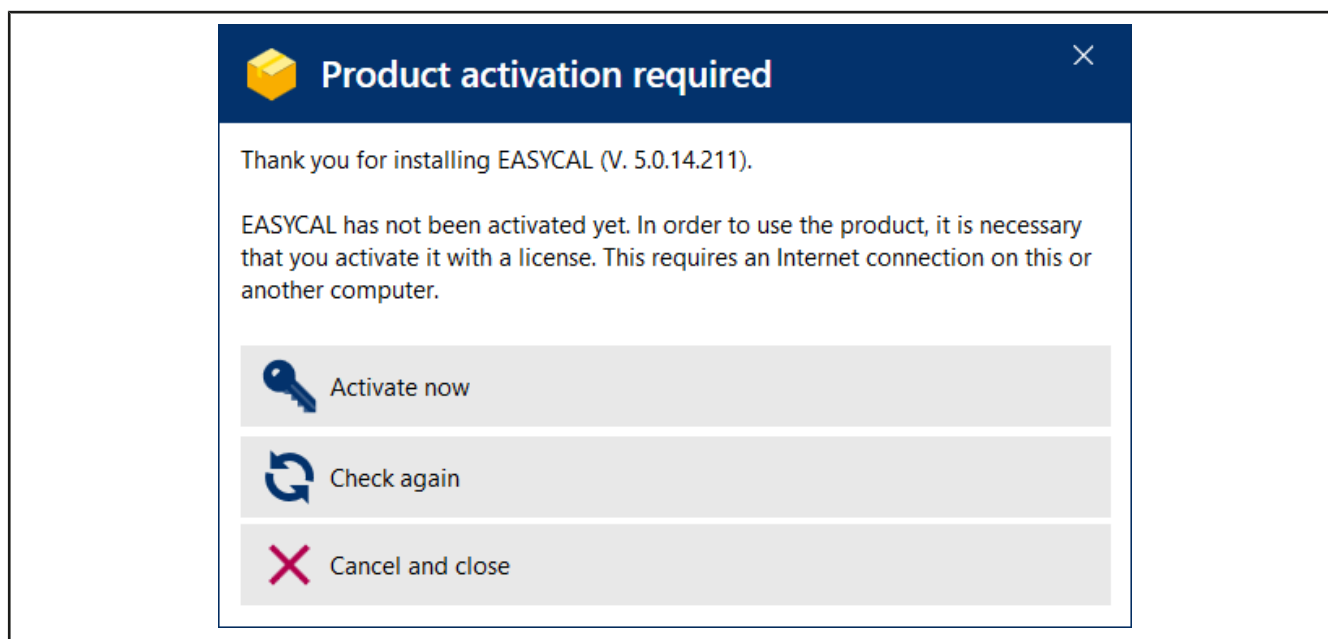
- c. You will be asked to activate EASYCAL™ 5.

- d. Click on Activate now.

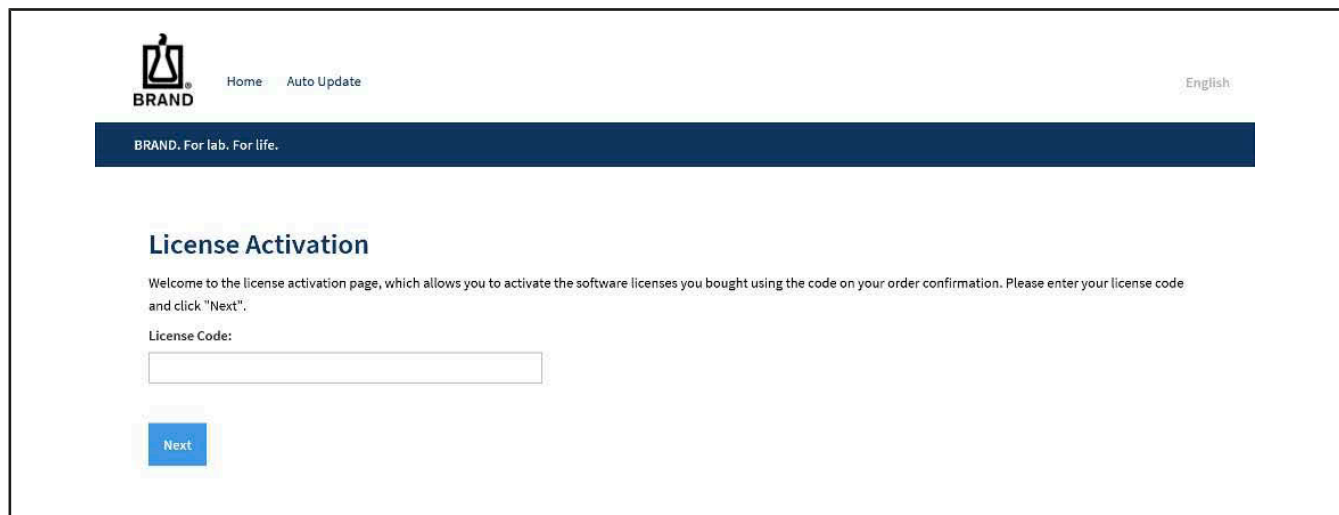➔ Your standard web browser opens with the page <https://license.brand.de/index.php>.

BRAND Home Auto Update English

BRAND. For lab. For life.

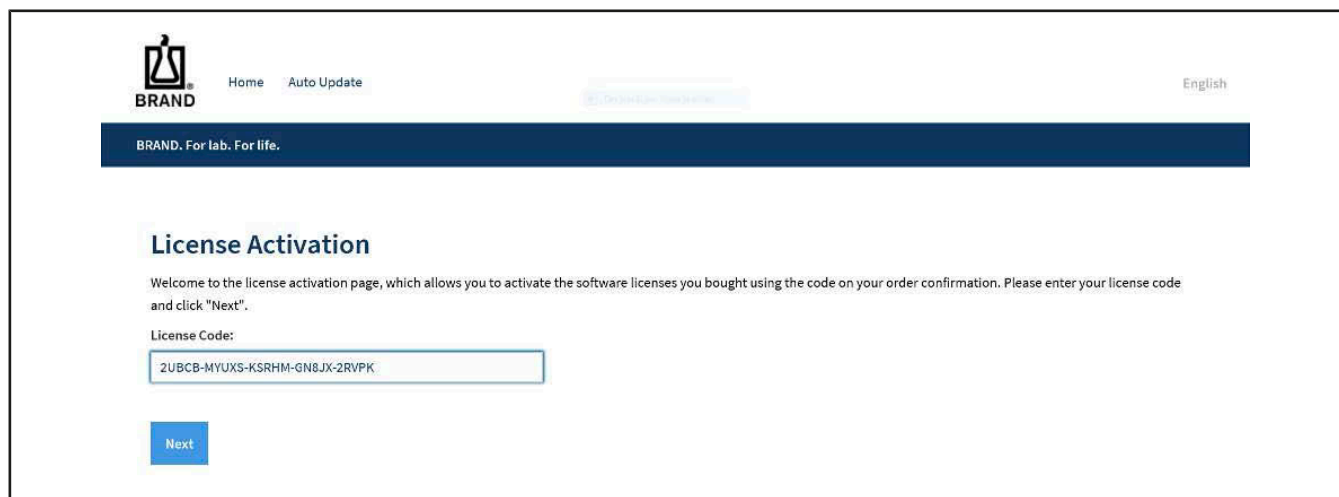
License Activation

Welcome to the license activation page, which allows you to activate the software licenses you bought using the code on your order confirmation. Please enter your license code and click "Next".

License Code:

Next

e. Enter your ticket (= license key) here.

BRAND Home Auto Update English

BRAND. For lab. For life.

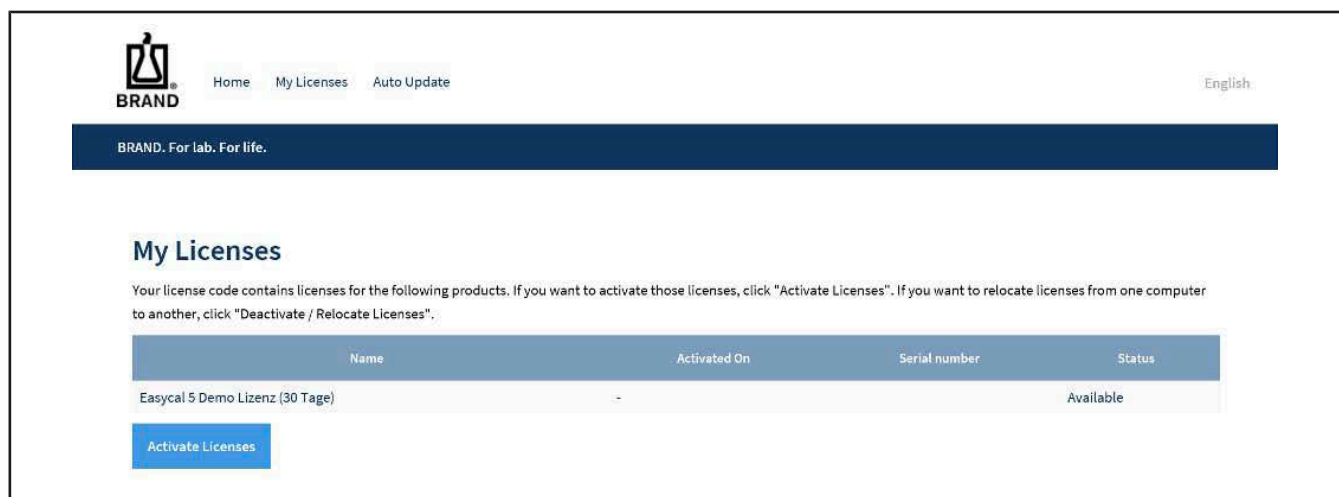
License Activation

Welcome to the license activation page, which allows you to activate the software licenses you bought using the code on your order confirmation. Please enter your license code and click "Next".

License Code:

Next

f. Click on Continue.

BRAND Home My Licenses Auto Update English

BRAND. For lab. For life.

My Licenses

Your license code contains licenses for the following products. If you want to activate those licenses, click "Activate Licenses". If you want to relocate licenses from one computer to another, click "Deactivate / Relocate Licenses".

Name	Activated On	Serial number	Status
Easycal 5 Demo Lizenz (30 Tage)	-		Available

Activate Licenses

g. Click on Activate License.

Activate Licenses

To activate your licenses:

1. Select the licenses you want to activate.
2. Select a license file to load the licenses into.
3. Click "Activate Selected Licenses Now".

☒	Name	Activated On	Serial number	Status
☒	EasyCal 5 Demo Lizenz (30 Tage)			Available

Select license file

select / create automatically

[Manual license transfer](#)

[Activate Selected Licenses Now](#)

h. If you do not change any of the settings, click on Activate Selected Licenses Now.

→ The online license transfer is carried out.

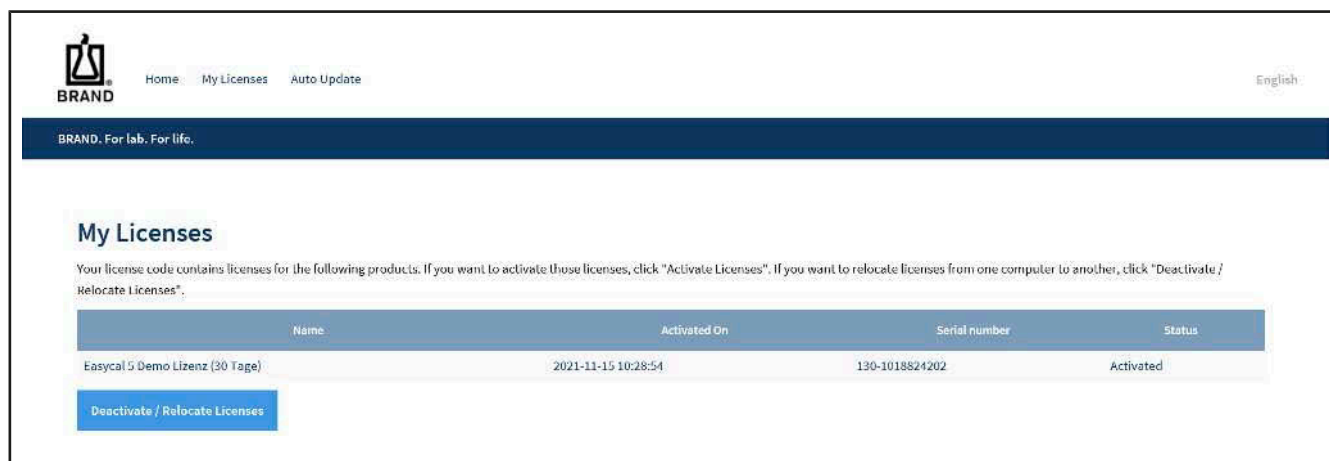
Online License Transfer

Starting license transfer.
 Downloading license template.
 Registering license template.
 Creating license request.
 Downloading license update.
 Importing license update into target.
 Creating receipt.
 Uploading receipt.

☒ License transfer completed successfully!


[OK](#)

- i. Click on OK.


- j. The license is now activated. You can close the browser window.

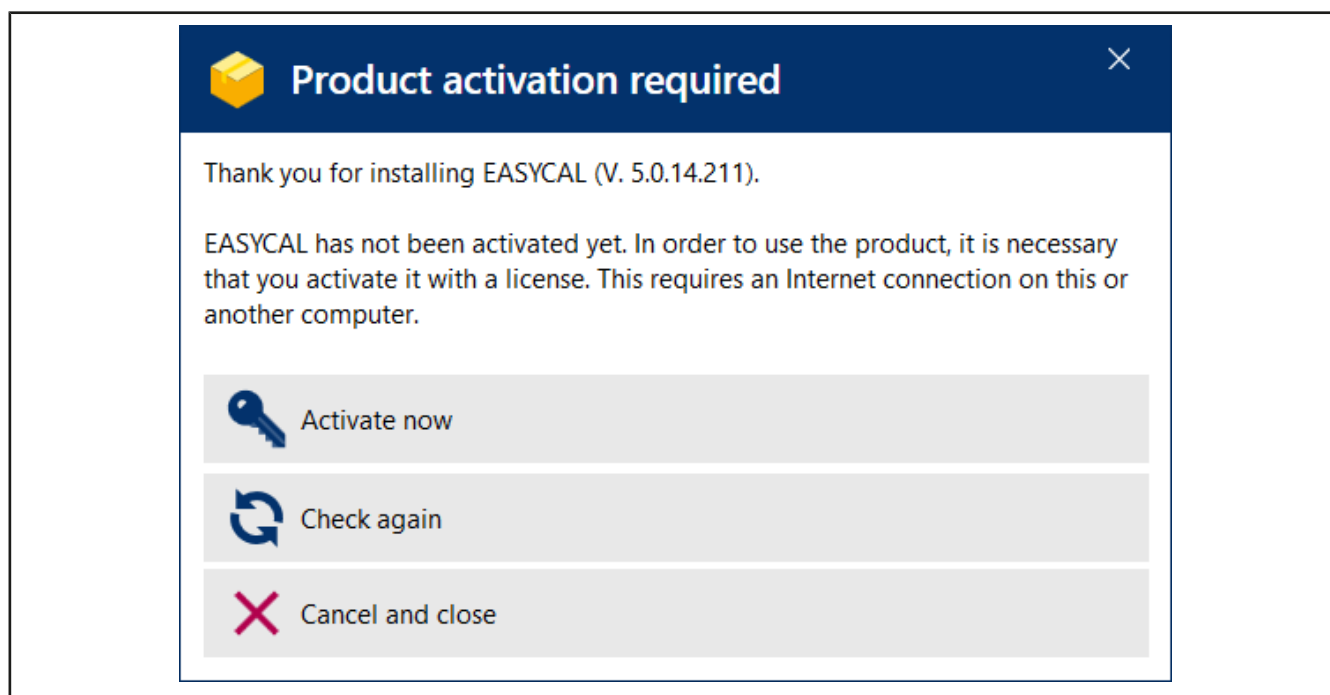
- k. Go to the product activation window.

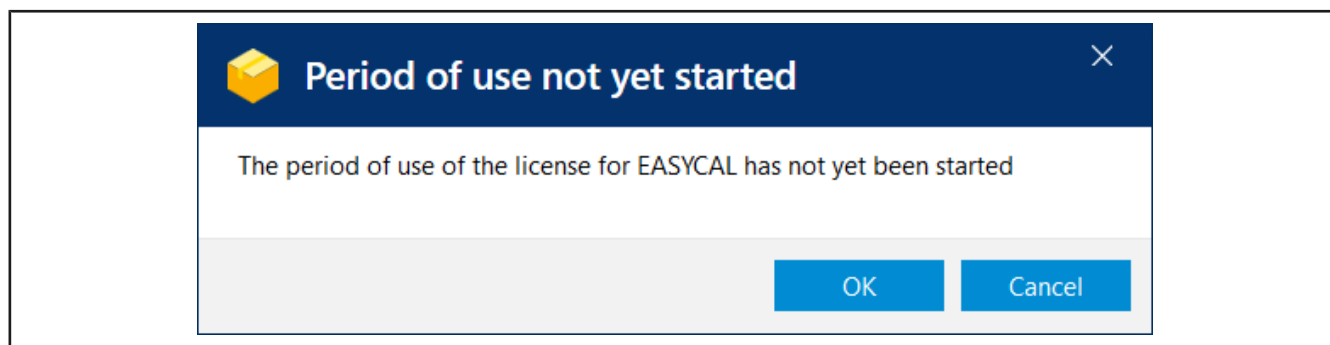
- l. Click "Check Again".

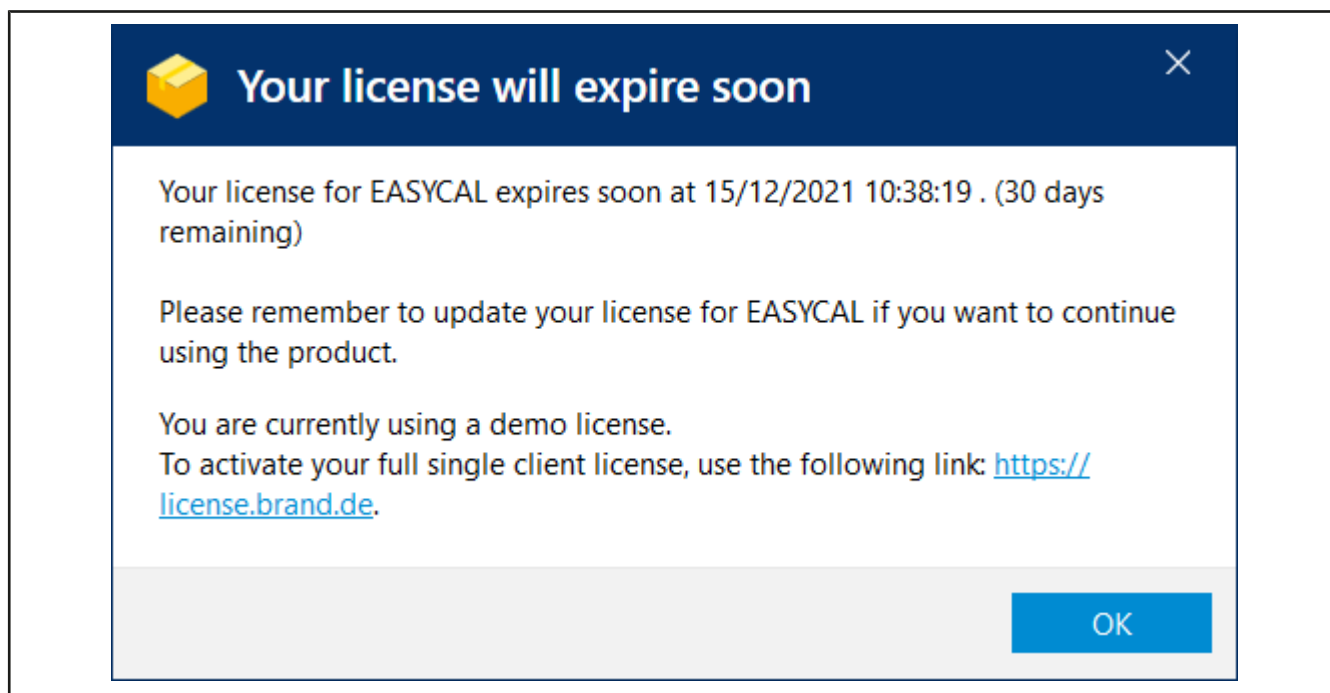
- m. Click on OK to start the usage period. An information window appears.

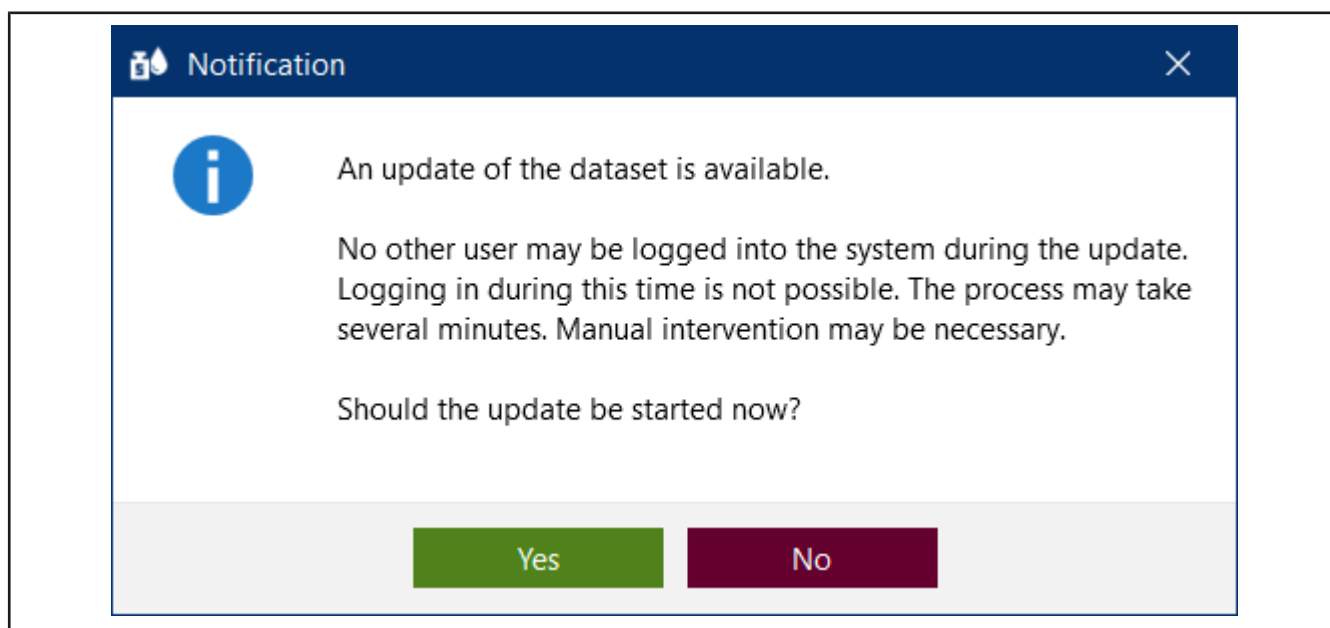
- n. Click on OK.
- o. At the same time, EASYCAL™ 5 is started and a notification appears.

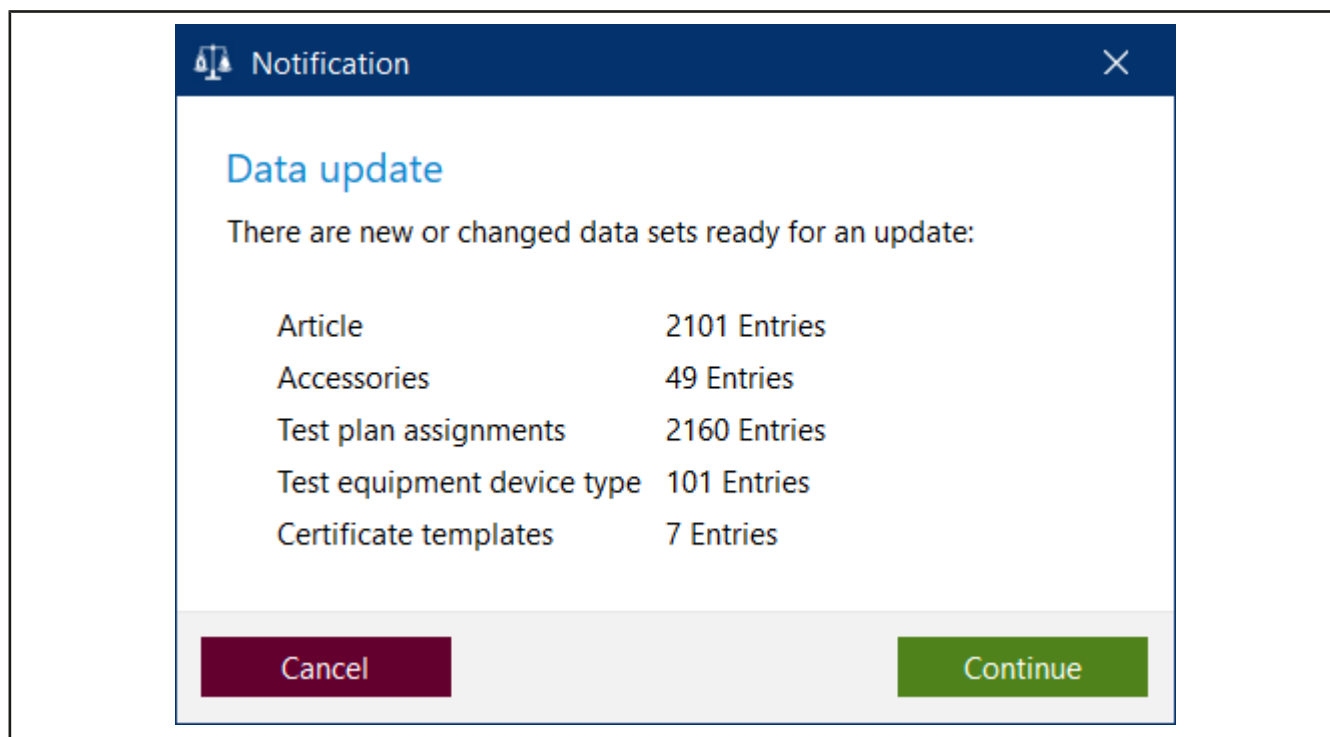
- p. Acknowledge the notification by clicking Yes so that the database is initially filled.
- q. A notification will appear after a short time.

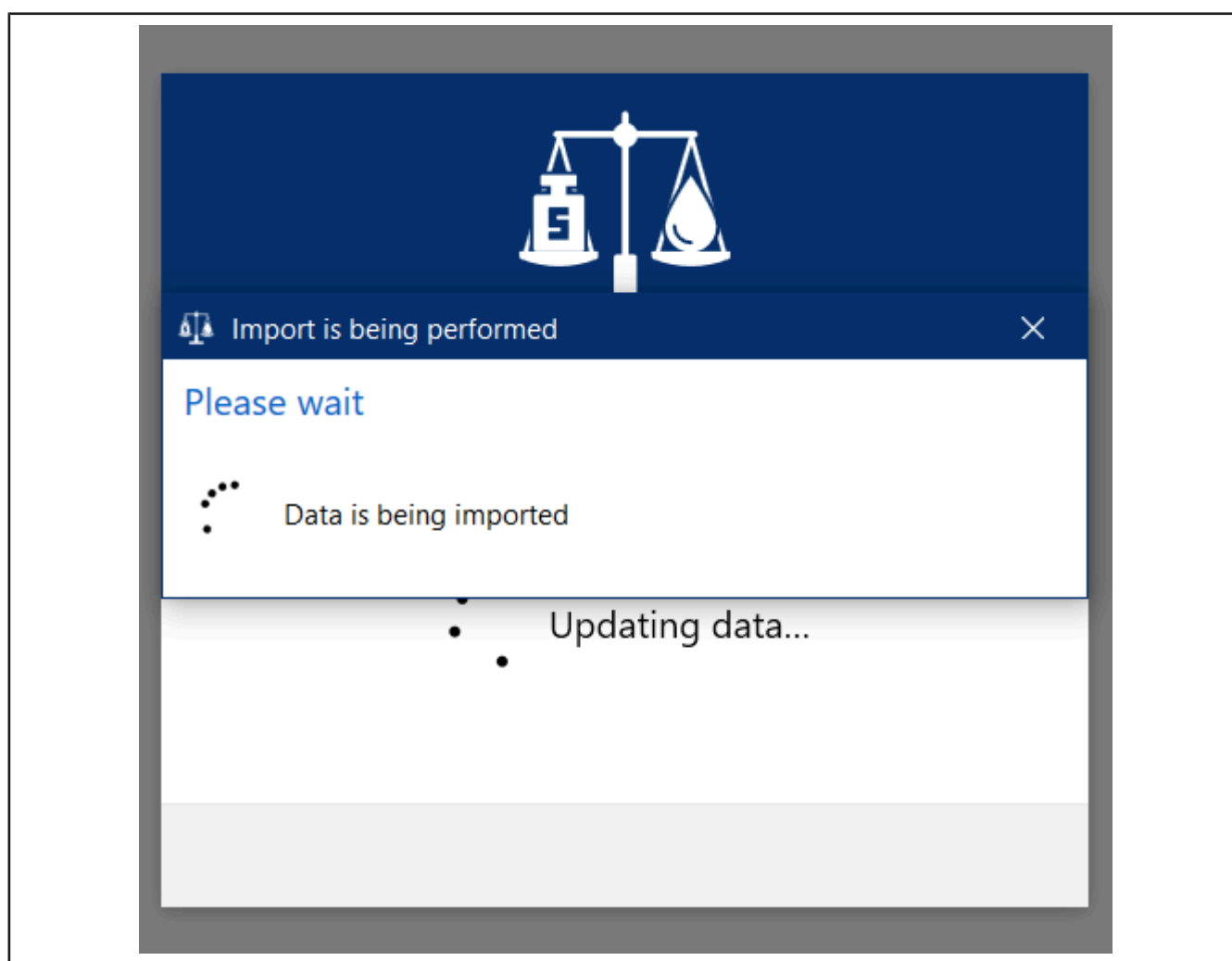
r. Close your entries by clicking Continue.

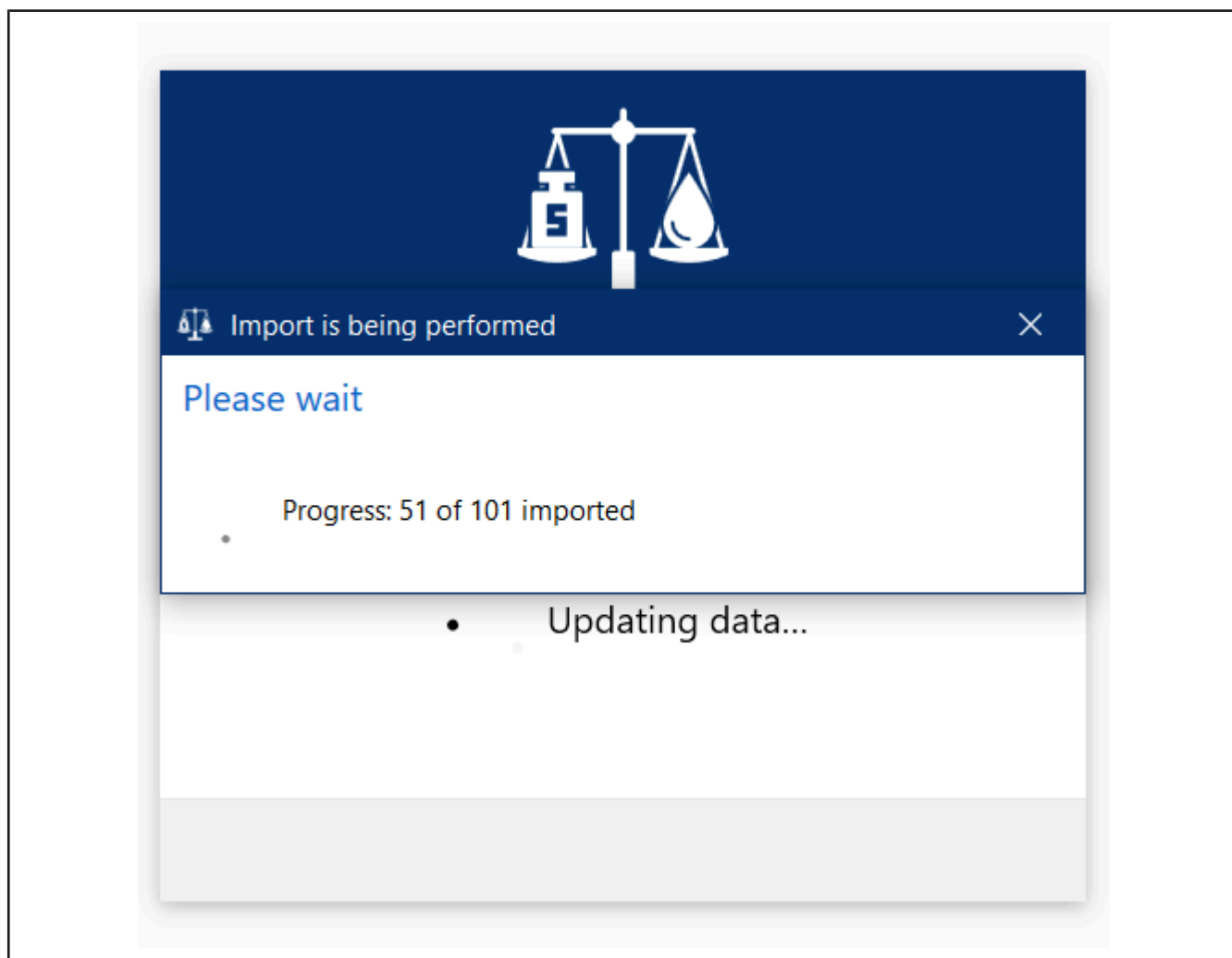- s. The new data is imported and the progress is displayed.

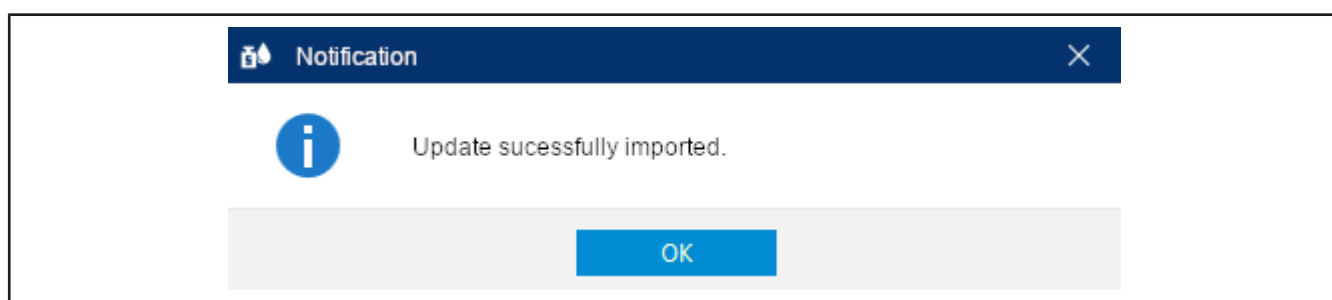
- After a while, a notification will appear that the update is complete.

- t. Acknowledge these by clicking on OK.

- EASYCAL™ 5 will start. Various messages will appear to show the progress.

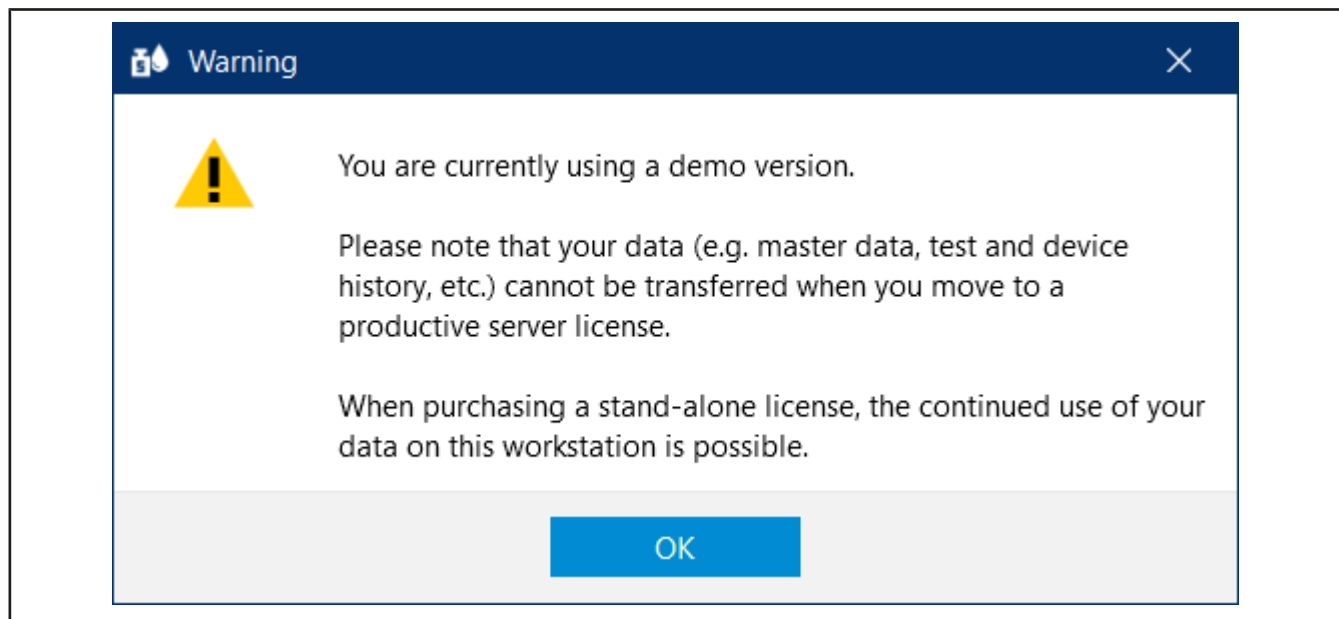- A warning appears that you are using a demo version. Read the notes.

- u. After you acknowledged the warning you can now test EASYCAL™ 5.
- You will be automatically logged in as "EASYCAL Admin". Note that EASYCAL™ 5 will always start in single user mode after installation. You can change this later in the "Global Settings", but shouldn't do so at this point in time.

3 First steps in EASYCAL™ 5.

3.1 Tutorial

This example shows the process of a test – from the creation of an order to the test to the printing of the certificate. The section refers to master data available in the demo license and serves as a first introduction to the program.

Login

If the program EASYCAL™ 5 is started, a log-in dialog appears. Here, the user logs in with user name and password.

3.1.1 Set up balance as test equipment

In order set up a balance as a piece of test equipment, proceed as follows. The set-up is always done in the order: Set up test equipment in the master data, set up test equipment at the workstation, and select test equipment for the orders.

Prerequisite:

- ✓ You have connected a balance to the workstation.
 - ✓ You know the connection parameters of the balance.
 - ✓ You know to which COM port the balance is connected at the workstation.
- a. Name balance (see Administration > Master data maintenance > Test equipment > Test equipment).
 - b. Set up location for test equipment (if not already done).
 - c. Configure the balance as a test equipment (see Test equipment device type). In doing so, you specify both the function and the connection of the balance to the workstation (see Measuring function, Communication).
- Exit this menu. Change to the Settings menu.
- d. Register the balance as test equipment at the workstation (see Settings > Workstation settings > Test equipment).
 - e. Select balance and the connection for the workstation. You can also test the connection to the balance.
- Exit this menu. Change to the Orders and calibration slips menu.
- f. You can select the connected balance in the order when selecting the test equipment. In the tutorial (see Determine test equipment, p. 50), in the user manual under Select test equipment for the test.

3.1.2 Creating order

- a. Navigate from Main menu to Orders and calibration slips.

Orders and calibration slips

Test type filter Not specified
Search
from 1/11/2022 to 1/26/2022

Order overview
Calibration slips

Order I	Test typ	Order number	Article	Status
2	SPEC	2022-01-25-15-2	EC-12484 - Burette BLAUBRAND®, straight stopcock, Schilling, class AS; 0.02 - 10 ml	Created
1	STD	123456	EC-704106 - Transferpette® Fixed volume; 5 µl	Finished

Back
Individual device test
Create order
Open

b. Click on Create order.

Select order type

Batch test

Use when testing random samples (of any size). Batch certificates are issued.

Individual test

Use for 100% test or test of a single device.

Each device receives its own certificate.

Back

c. Click on Individual test.

d. For Test type, select the value STD – Standard values.

Test types organize your test procedures. You cannot create additional test types; instead, you must choose from the test types already stored. The test plans are linked to the test type as well as the certificate templates. In the user administration, the user must be given the respective access right for the corresponding test type. Because you are currently logged in as Global Admin, you have access to all data.

Individual test - Create order

Test type: **STD - standard values** | Number of devices: **1** | Customer: [dropdown]

Order no.: [input] | Order no.: [input]

Article code: [input] | OC-Number: [input]

Article: [input]

☐ Prepare calibration slips

☐ Pre- and post-test required

Notes on the order: [text area]

End customer: [dropdown]

Volume ranges: [input]

Test plan: [dropdown]

Back | **Start test**

e. Click on the button at Item code in order to display already created articles.

Individual test - Create order

Test type: **STD - standard values** | Number of devices: **1** | Customer: [dropdown]

Order no.: [input] | Order no.: [input]

Article code: [input] | OC-Number: [input]

Article: [input]

☐ Prepare calibration slips

☐ Pre- and post-test required

Notes on the order: [text area]

End customer: [dropdown]

Volume ranges: [input]

Test plan: [dropdown]

Back | **Start test**

Select article

Search: [input]

Article code	Vei	Manufacturer	Product	Product variation	Volume range	Status
10_copy	1	BRAND GMBH I + CO KG	Dispensette® S Organic	Fixed volume, Special fixed volume	0.05 - 5 ml	Released
10002	0	BRAND GMBH I + CO KG	Burette	SILBERBRAND, ohne I lahn, class B, Economy	0.1 - 25 ml	Released
10012	0	BRAND GMBH I + CO KG	Compact Burette	SILBERBRAND, class B, Economy	0.1 - 25 ml	Released
100710	6	Eppendorf AG	Research® plus	Fixed volume, medium grey	10 µl	Released
100711	7	Eppendorf AG	Research® plus	Fixed volume, light grey	20 µl	Released
100711_copy	0	Eppendorf AG	Research® plus	Fixed volume, light grey	20 µl	Released
100212	6	Eppendorf AG	Research® plus	Fixed volume, yellow	10 µl	Released
100212_copy	0	Eppendorf AG	Research® plus	Fixed volume, yellow	10 µl	Released
100213	6	Eppendorf AG	Research® plus	Fixed volume, yellow	20 µl	Released
100214	7	Eppendorf AG	Comforpette®	Fixed volume	20 µl	Released
100215	6	Eppendorf AG	Research® plus	Fixed volume, yellow	50 µl	Released
100216	6	Eppendorf AG	Research® plus	Fixed volume, yellow	100 µl	Released
100217	6	Eppendorf AG	Research® plus	Fixed volume, yellow	200 µl	Released
100218	6	Eppendorf AG	Research® plus	Fixed volume, blue	200 µl	Released
100219	6	Eppendorf AG	Research® plus	Fixed volume, blue	250 µl	Released
100220	6	Eppendorf AG	Research® plus	Fixed volume, blue	500 µl	Released
100221	6	Eppendorf AG	Research® plus	Fixed volume, blue	1000 µl	Released
100310	6	Eppendorf AG	Research® plus	Adjustable volume, dark grey	0.1 - 2.5 µl	Released
100311	8	Eppendorf AG	Research® plus	Adjustable volume, dark grey	0.5 - 10 µl	Released

© Daniel Iking | Last status: The application is prepared

f. Enter Transferpette® S 1000 in the search field. As you type, you will see the article selection change. This behavior is the same in all search fields of EASYCAL™ 5. We recommend always using the search function first because EASYCAL™ 5 comes with an extensive master database.

The prefix EC- indicates that it is a prefabricated and unchangeable data set. If you would like to change

these records, copy the record, and edit and save it under a different name. For more information, see “Help” in EASYCAL™ 5.

Individual test - Create order

Test type: **STD - standard values** | Number of devices: **1** | Customer:

Order no.: | Order no.: | OC-Number:

Article code: | Article:

Volume ranges:

Test plan:

Select article

Search:

Article code	Vei	Manufacturer	Product	Product variation	Volume range	Status
700762	2	Treff AG	TreffLab-Transferpette®	Fixed volume	1000 µl	Released
700700	2	Treff AG	TreffLab-Transferpette®	Adjustable volume	100 - 1000	Released
700784	2	Treff AG	TreffLab-Transferpette®	Adjustable volume	1000 - 1000	Released
700791	0	Treff AG	TreffLab-Transferpette®	Adjustable volume	100 - 1000	Released
700793	0	Treff AG	TreffLab-Transferpette®	Adjustable volume	1000 - 1000	Released
700965	0	BRAND GMBH + CO KG	Transferpette® α	1.0 ml, 1-Stop	1000 µl	Released
700966	0	BRAND GMBH + CO KG	Transferpette® α	100 - 1000 µl, 1-Stop	100 - 1000	Released
704162	3	BRAND GMBH + CO KG	Transferpette®	Fixed volume	1000 µl	Released
704180	4	BRAND GMBH + CO KG	Transferpette®	Adjustable volume	100 - 1000	Released
704762	5	BRAND GMBH + CO KG	Transferpette® S	Fixed volume	1000 µl	Released
704780	6	BRAND GMBH + CO KG	Transferpette® S	Adjustable volume	95 - 1000	Released
704780-4711	0	BRAND GMBH + CO KG	Transferpette® S	Adjustable volume	100 - 1000	Released
704780-4712	0	BRAND GMBH + CO KG	Transferpette® S	Adjustable volume	100 - 1000	Released
704780-4713	0	BRAND GMBH + CO KG	Transferpette® S	Adjustable volume	100 - 1000	Released
704784	4	BRAND GMBH + CO KG	Transferpette® S	Adjustable volume	1000 - 1000	Released
704941	2	BRAND GMBH + CO KG	Transferpette® S	1.0 ml, 1-Stop	1000 µl	Released
704945	2	BRAND GMBH + CO KG	Transferpette® S	100 - 1000 µl, 1-Stop	100 - 1000	Released
704980	5	BRAND GMBH + CO KG	Transferpette® α	Adjustable volume	100 - 1000	Released
704984	5	BRAND GMBH + CO KG	Transferpette® α	Adjustable volume	1000 - 1000	Released

Back

Daniel Ivrig | Last status: The application is prepared

- g. Select the article code EC-704780. Confirm your selection with a double click.
- h. For Test plan, select the appropriate test plan. As a rule, you will find 1–2 entries here.
- i. The general conditions for the test as well as the defect limits of the assigned article are stored in the test plan.

Individual test - Create order

Test type: **STD - standard values** | Number of devices: **1** | Customer:

Order no.: | Order no.: | OC-Number:

Article code: **704780** | Article: **Transferpette® S**
Adjustable volume
BRAND GMBH + CO KG

Volume ranges: **1000µl (95 - 1000)**

Test plan: **4437.0 - EC-STD-ISO-PiPip-var-LHDiv**

Notes on the order:

End customer:

Back | **Start test**

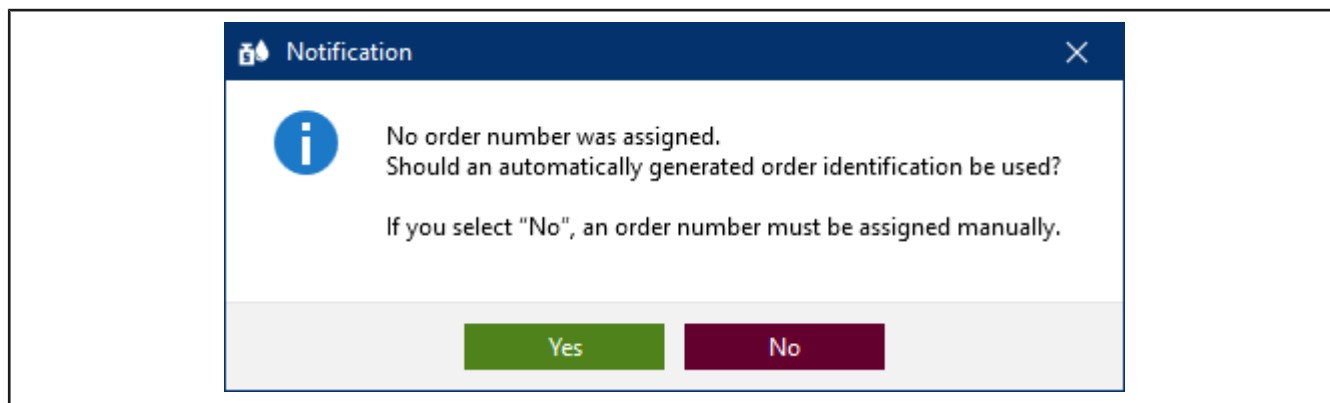
- j. Click on Start test.

A notification that no order number has been assigned appears.


If there are specifications in your company/quality management process for organizing tests, select no, and assign an order number.

If there are no specifications for organizing tests, select Yes and EASYCAL™ 5 will assign a unique order number

We recommend deciding on a uniform approach at an early stage.

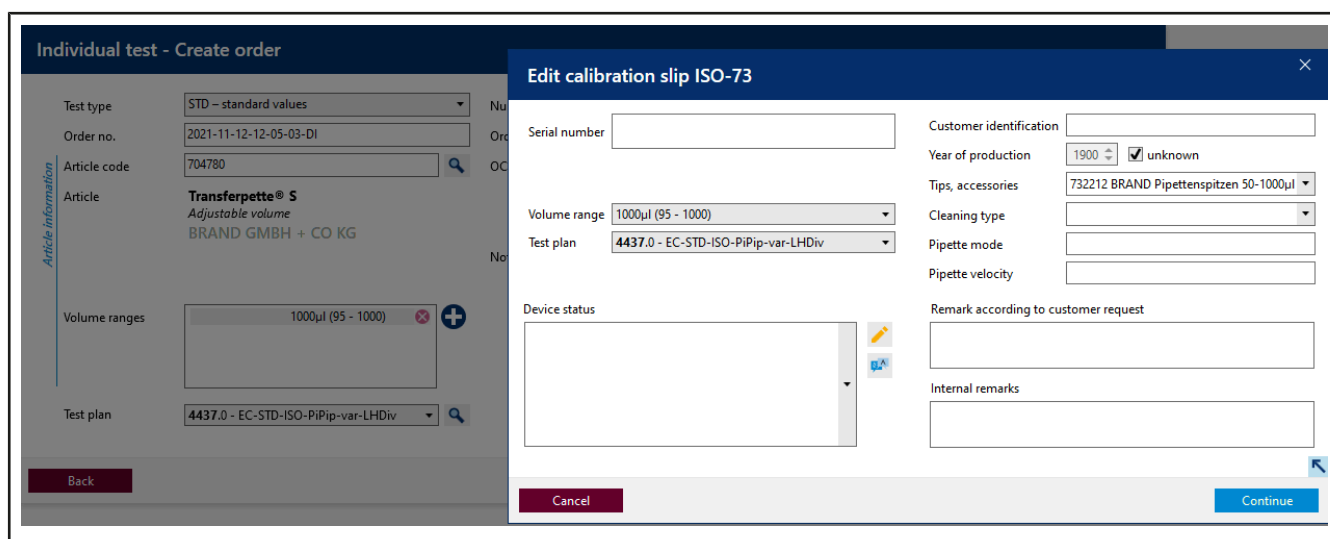k. Continue the tutorial by clicking on Yes. The job number is created based on the current time stamp.

l. You can enter a serial number for the identification of the device. This is necessary for real tests. However, it is not necessary for the tutorial.

m. Click on Continue. The test order has been created.

The window for starting the actual test will open.

3.1.3 Determine test equipment

BRAND EASYCAL™ 5

2021-11-12-12-05-03-DI / ISO-73 - - 704780 BRAND GMBH + CO KG Transferpette® S Adjustable volume 95 - 1000µl

Test equipment > 2. TV > 1. TV > NV

✓ / ✗ A / B / C Measurement + i

Air temperature	Thermometer		351000-71
Temperature of the medium	Thermometer		351000-71
Relative humidity	Hygrometer		600029-10
Air pressure	Barometer		600029-10
Weight	Single-channel balance	Calibrate now	20210708-RZ-01
Weight	Single-channel balance	Calibrate now	20210708-RZ-01
Weight	Single-channel balance	Calibrate now	20210708-RZ-01

Back

Previous device Previous characteristic Next characteristic Additional device

Device evaluation Complete test Print preview

Prerequisite:

- ✓ The order is created.
- a. The necessary parameters as well as the test equipment type and the test equipment (test equipment number) are preselected for the test equipment. Because only one piece of test equipment is created, we leave it at this selection.

3.1.4 Performing the test

- a. Continue by clicking on Next characteristic.

You see the central input screen of the measurement functionality of EASYCAL™ 5.

BRAND - EASYCAL™ 5

2021-11-12-12-05-03-DI / ISO-73 - - 704780 BRAND GMBH + CO KG Transferpette® S Adjustable volume 95 - 1000µl

Ambient data

Request ambient data

Target value: 100 µl

Number: 0 / 10

Test equipment: 2. TV, 1. TV, NV

Graph of the measured values

Statistics of device

Statistics of order

Back, Previous device, Previous characteristic, Next characteristic, Additional device, Device evaluation, Complete test, Print preview

- b. Enter fictitious weight values for Neue Messung. Start with 0, 1, and click on “Apply measured value”. A window in which you must enter or complete the measured values appears.

The selection of the separator (comma or dot) is done via the language selection in the software.

Enter measured values

Temperature difference	0.00 °C	0.5 °C	
Weight	0.1 g		☒ Manual
Air temperature	0 °C	(20°C - 25°C)	☒ Manual
Temperature of the medium	0 °C	(20°C - 25°C)	☒ Manual
Relative humidity	0 %rh	(40%rh - 65%rh)	☒ Manual
Air pressure	0 mbar	(960mbar - 1025mbar)	☒ Manual

Cancel Apply

- c. Enter fictitious values for the environment parameters as follows:

Enter measured values

Temperature difference	0.00 °C	0.5 °C	
Weight	0.1	g	☒ Manual
Air temperature	21	°C	(20°C - 25°C) ☒ Manual
Temperature of the medium	21	°C	(20°C - 25°C) ☒ Manual
Relative humidity	40	%rh	(40%rh - 65%rh) ☒ Manual
Air pressure	961	mbar	(960mbar - 1025mbar) ☒ Manual

Cancel
Apply

- d. Click on Apply. The first measurement is hereby completed. The measured value and the volume calculated from it appear in the table.

Test equipment > 2. TV > 1. TV > NV				
Number	Time	Measured value	Volume [µl]	Sta
1	11/12/2021 2:06:42 PM	0.10000 g	100.3	✓

- e. Now enter the value 0.1 g as the measured value eight times in the same way. The table is filled up.

2021-11-12-14-10-58-DI / ISO-74 - - 704780 BRAND GMBH + CO KG Transferpette® S Adjustable volume 95 - 1000µl

Ambient data: 21 °C, 21 °C, 40 %rh, 961 mbar

Request ambient data

Target value: 100 µl

Number: 9 / 10

Letzte Messung: 0.1 g, 1003.01211312 µl/g, 100.3 µl, 2:13:39 PM

Neue Messung: g, µl/g, µl

Apply measured value

Test equipment disconnected

Fast measurement, Automatic printing, Automatic tare, Allow all measured values, Check measured values before entering

Discard entry, Undo discard

Test equipment > 2. TV > 1. TV > NV

Number	Time	Measured value	Volume [µl]	Sta
2	11/12/2021 2:13:27 PM	0.10000 g	100.3	✓
3	11/12/2021 2:13:29 PM	0.10000 g	100.3	✓
4	11/12/2021 2:13:30 PM	0.10000 g	100.3	✓
5	11/12/2021 2:13:32 PM	0.10000 g	100.3	✓
6	11/12/2021 2:13:34 PM	0.10000 g	100.3	✓
7	11/12/2021 2:13:35 PM	0.10000 g	100.3	✓
8	11/12/2021 2:13:37 PM	0.10000 g	100.3	✓
9	11/12/2021 2:13:39 PM	0.10000 g	100.3	✓

Graph of the measured values

Statistics of device

100.3 ± 0

Accuracy: 0 % / 8 % ✓

Coefficient of variation: 0 % / 3 % ✓

Mean: 100.3 µl

Standard deviation: 0 µl

Expanded measurement uncertainty: 0 µl

Statistics of order

Accuracy: 0 %

Mean: 100.3 µl

Standard deviation of the mean: 0 µl

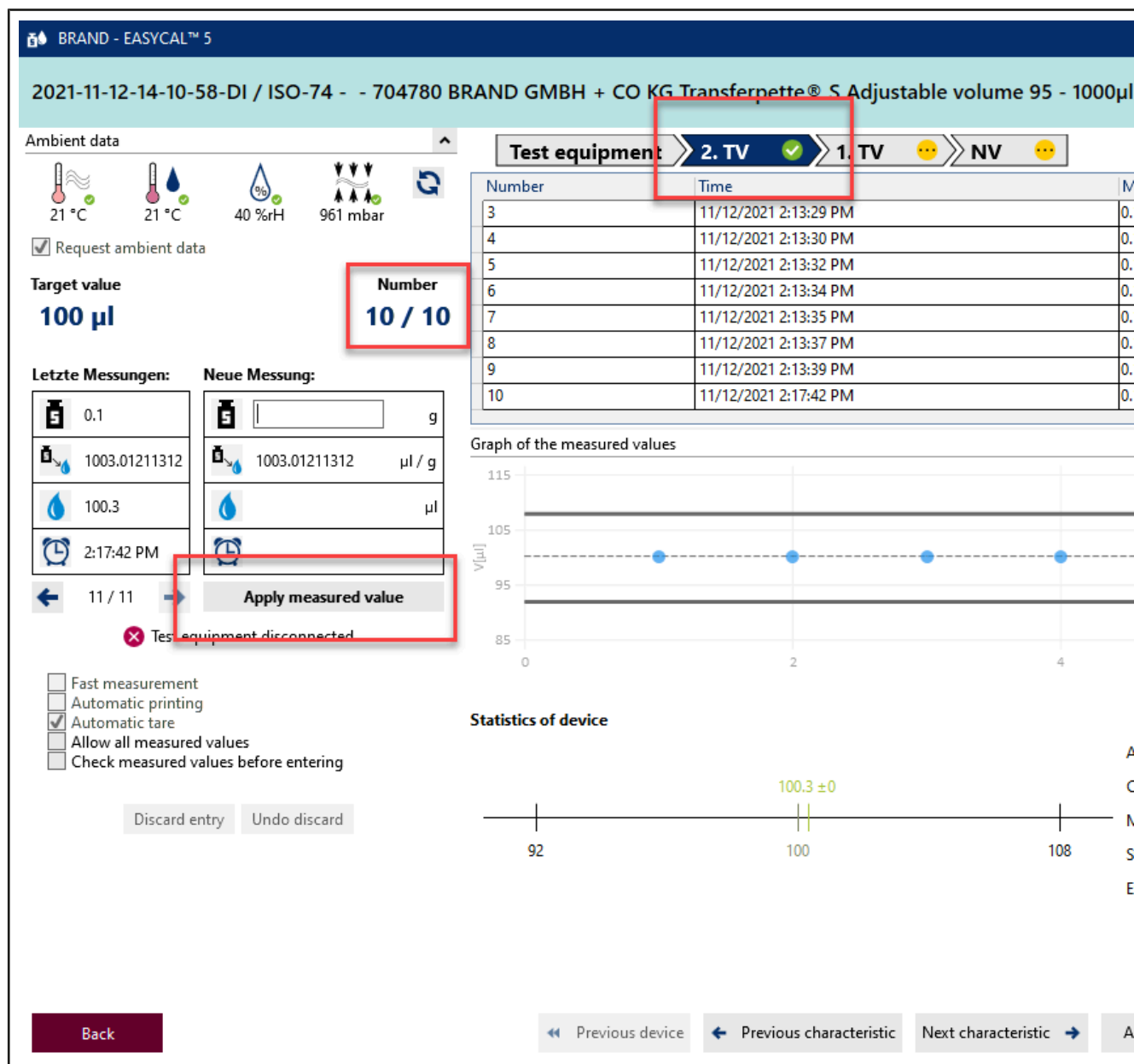
Coefficient of variation of accuracy: 0 %

Mean value CV: 0 %


CPK: 0

Back, Previous device, Previous characteristic, Next characteristic, Additional device, Device evaluation, Complete test, Print preview

- f. The light blue color in the section above the table shows that the order is being processed. Click the icon in order show and hide the extended certificate information.
- g. Enter 0.1 again as the measured value, and pay attention to the progression display above the measured value table.

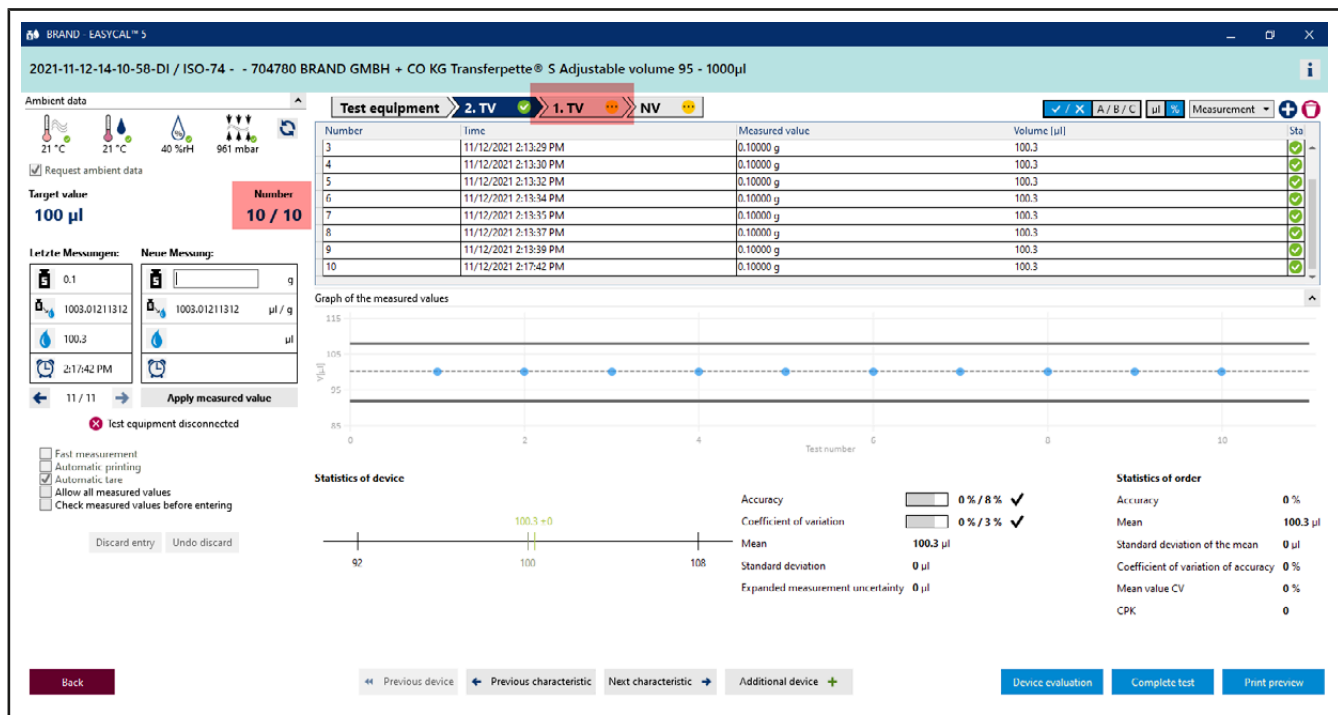- h. With input of the 10th measured value, all required measurements for characteristic 2 . TV have been performed. The characteristic gets a green check. You can enter further measured values; however, this is not required.


i. Click on 1. TV in the progress bar in order to jump to the next characteristic.

j. For Neue Messung, enter the following values in sequence:


0.498
0.499
0.500

Repeat twice.

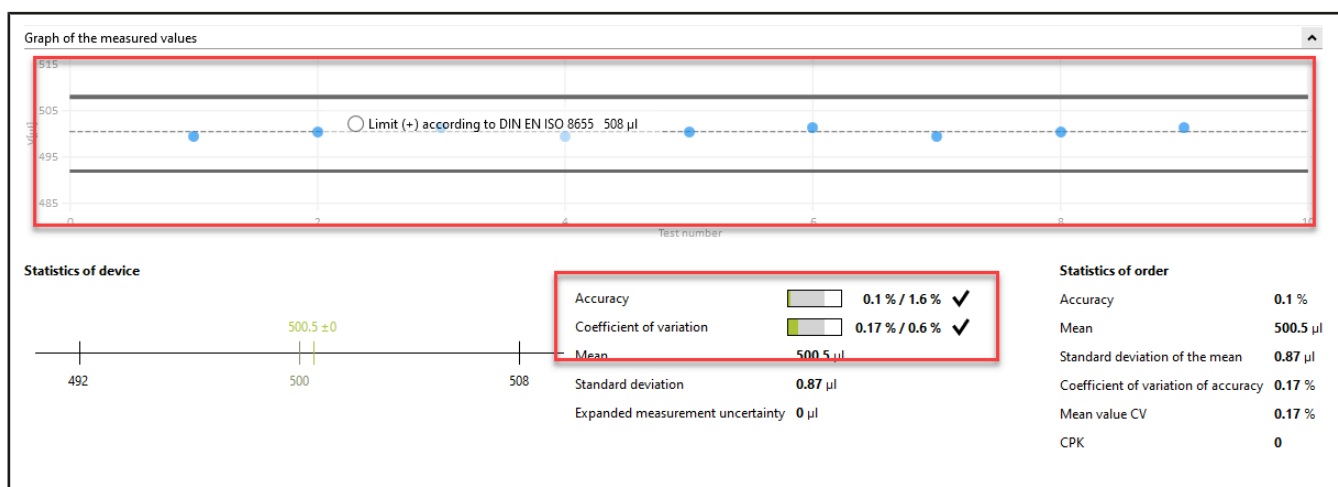
➔ You can see the position of the individual measured values in relation to defined error limits (solid lines) and the calculated mean value in the display Graph of the measured values. Under Statistics of device, you can see the calculated actual values for accuracy and coefficient of variation compared with the defined limits.

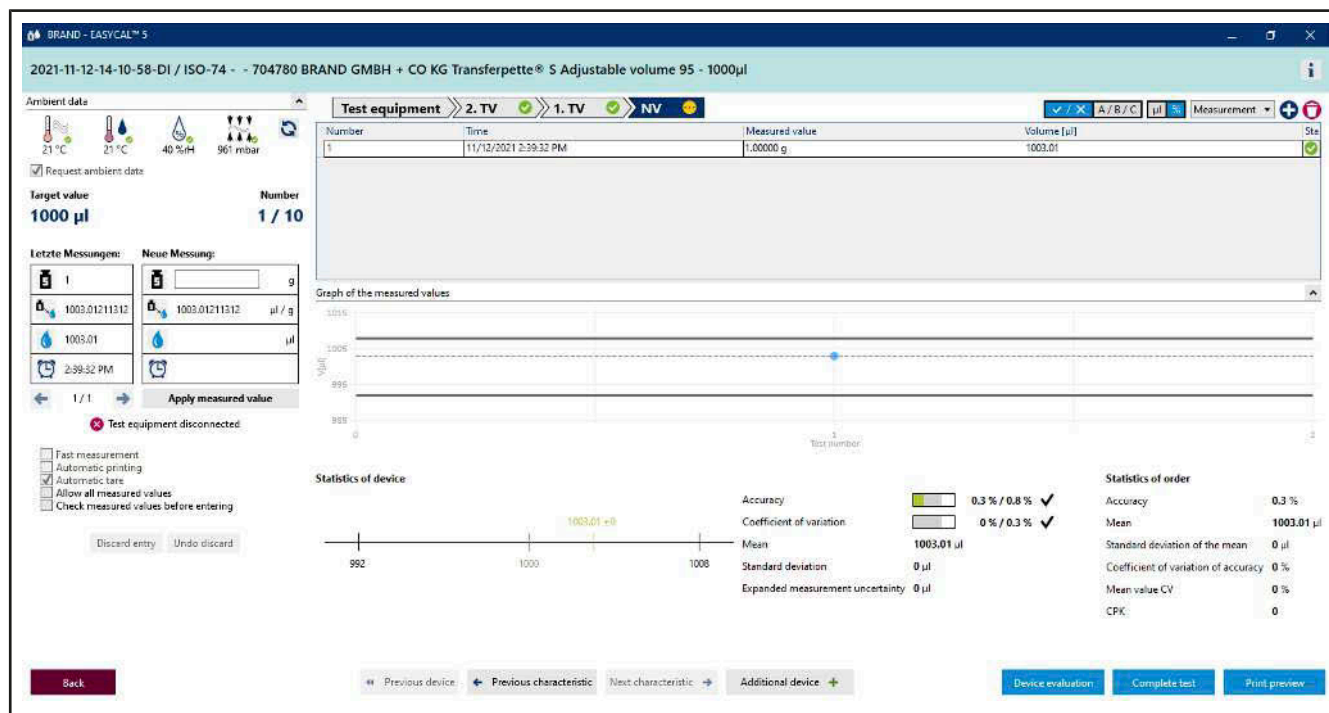- k. Enter 0.498 again as the measured value, confirm the environmental parameters, and click NV in the progress display. Then enter 1 as the measured value.

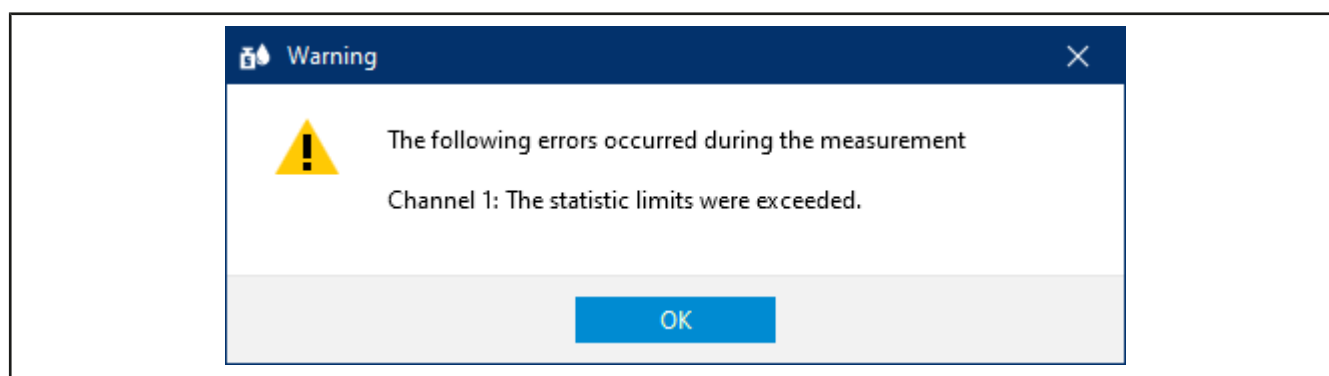- l. We assume that you have made a typing error. To do this, enter 1.1 as the measured value, and confirm the environmental parameters. The following warning appears:

- m. Acknowledge this warning by clicking OK.

You can see that the different areas of the input screen react according to the new measured value. You cannot delete this measured value; however, you can discard it. The value remains visible but is no longer used for statistics.


The screenshot shows the EASYCAL 5 software interface. At the top, it displays the date and time: 2021-11-12-14-10-58-DI / ISO-74 - 704780 BRAND GMBH + CO KG Transferpette® S Adjustable volume 95 - 1000µl. Below this, there are tabs for 'Test equipment' (2. TV, 1. TV, NV) and a 'Measurement' dropdown menu. A table lists measurements with columns for Number, Time, Measured value, and Volume [µl]. The first measurement is 1, 11/12/2021 2:39:37 PM, 1.00000 g, 1003.01 µl. The second measurement is 2, 11/12/2021 2:40:16 PM, 1.10000 g, 1103.31 µl. Below the table, there is a 'Graph of the measured values' showing a line graph with two data points. To the left of the graph, there are checkboxes for 'Fast measurement', 'Automatic printing', 'Automatic tare', 'Allow all measured values', and 'Check measured values before entering'. A 'Discard entry' button is highlighted with a red box. To the right of the graph, there are 'Statistics of device' and 'Statistics of order' sections. The 'Statistics of device' section shows a mean value of 1053.16 µl and a standard deviation of 71 µl. The 'Statistics of order' section shows an accuracy of 5.3 % and a coefficient of variation of 6.73 %.

- n. Click on the number of the measurement in the table, and then click on “Discard entry”.

A message appears.


The screenshot shows a 'Discard measurements' dialog box. The dialog box has a title bar with a close button (X). The main text says 'Specify a reason to discard the selected measurements.' Below this text is a drop-down menu with the option 'identified as outliers' selected. At the bottom of the dialog box, there are two buttons: 'Cancel' and 'Continue'.

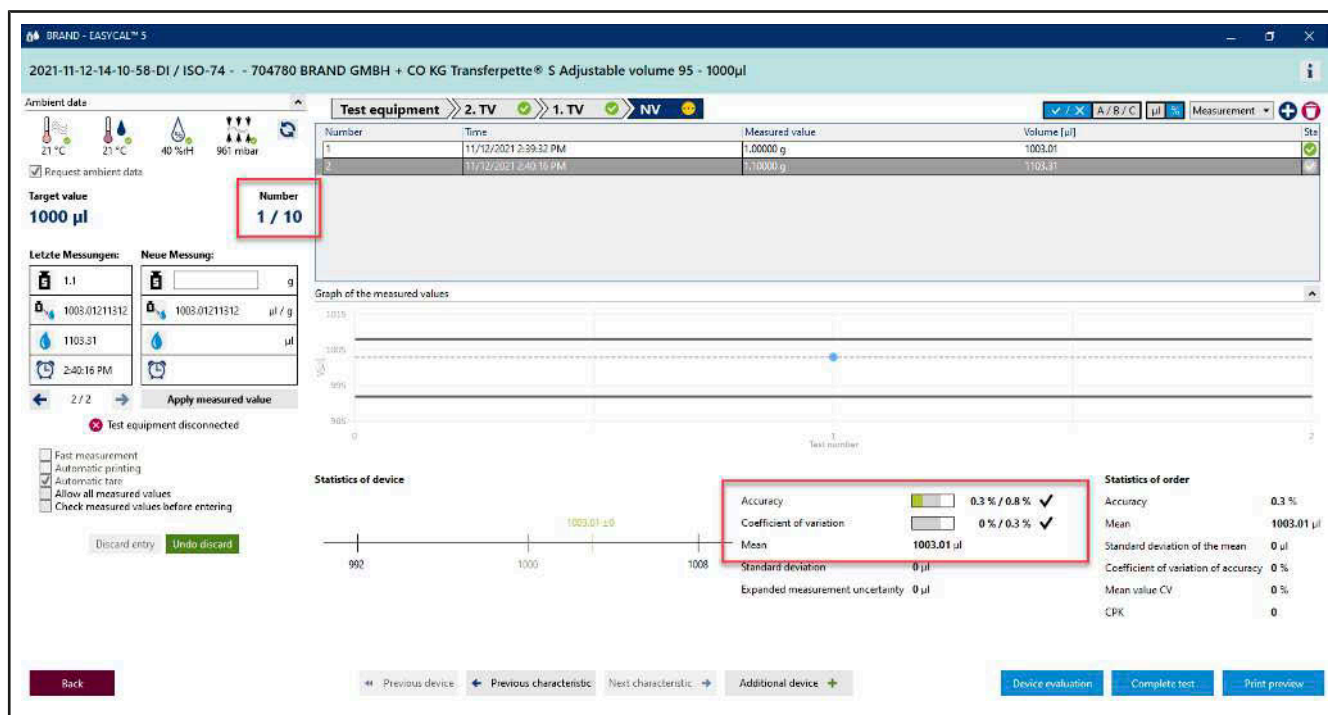- o. Specify a reason (e.g., Identified as outliers.). You can also select a reason from the drop-down menu. These texts can be maintained in the master data. Confirm your selection by clicking on Continue.

p. The measured value 1.1 is not considered further in the evaluation.

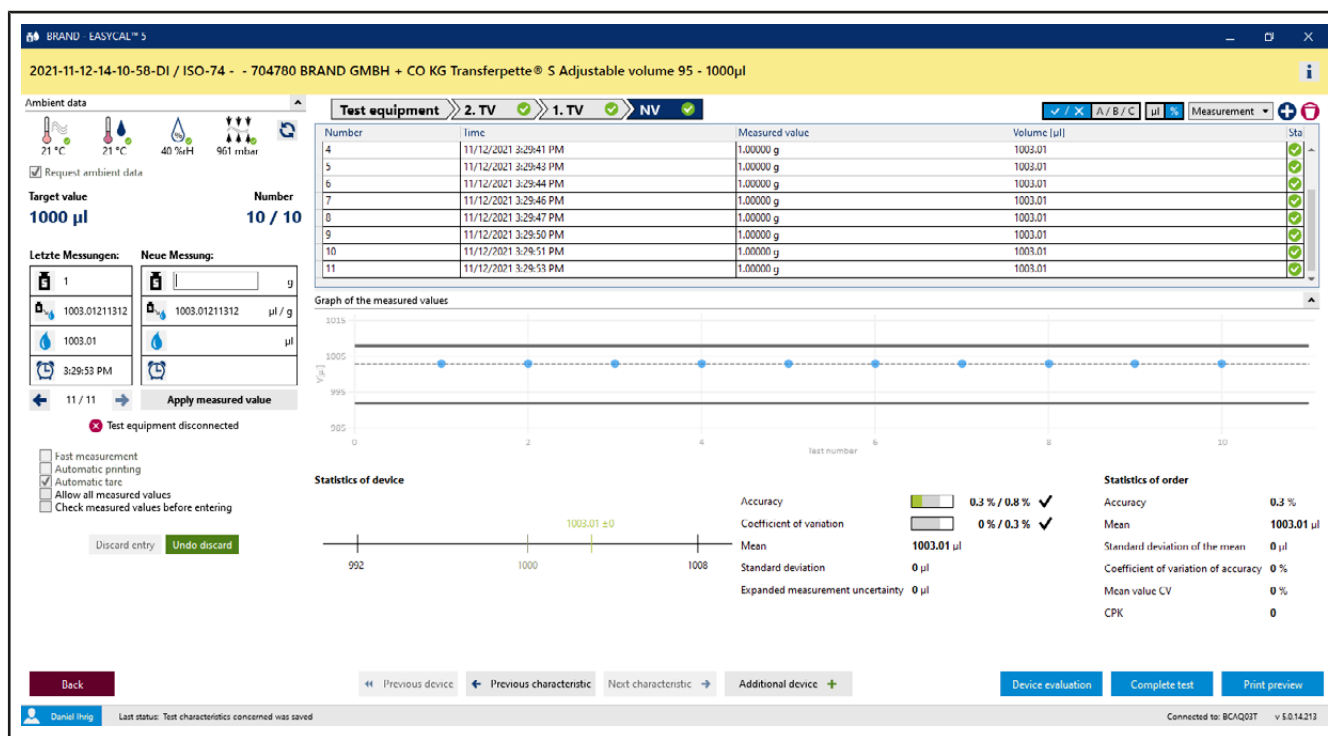q. Continue entering measured values until 10/10 is displayed for “Number”. When the last required measured value is entered, the color in the section above the table changes from light blue to yellow. This signals to you that the order is complete.

Furthermore, all characteristics in the progress bar are marked with a green check. The discarded measured value is highlighted in pink.

3.1.5 Complete test

The screenshot shows the EASYCAL™ 5 software interface. At the top, the title bar reads 'BRAND EASYCAL™ 5'. Below it, the header displays '2021-11-12-14-10-58-DI / ISO-74 - 704780 BRAND GMBH + CO KG Transferpette® S Adjustable volume 95 - 1000µl'. The main area is divided into several sections:

- Ambient data:** Shows temperature (21 °C), humidity (40 %rH), and pressure (961 mbar).
- Test equipment:** Displays '2. TV', '1. TV', and 'NV' with status indicators.
- Target value:** Set to '1000 µl'.
- Number:** Set to '10 / 10'.
- Letzte Messungen:** Shows the last measurement with a value of 1003.01 µl.
- Neue Messung:** Shows the new measurement with a value of 1003.01 µl.
- Table of measured values:** A table with columns for Number, Time, Measured value, and Volume [µl]. It lists 10 measurements, all with a measured value of 1003.01 µl.
- Graph of the measured values:** A line graph showing the measured values over time, with a mean value of 1003.01 µl.
- Statistics of device:** Shows accuracy (0.3 % / 0.8 %), coefficient of variation (0 % / 0.3 %), mean (1003.01 µl), standard deviation (0 µl), and expanded measurement uncertainty (0 µl).
- Statistics of order:** Shows accuracy (0.3 %), mean (1003.01 µl), standard deviation of the mean (0 µl), coefficient of variation of accuracy (0 %), mean value CV (0 %), and CPK (0).
- Buttons:** Includes 'Back', 'Previous device', 'Previous characteristic', 'Next characteristic', 'Additional device', 'Device evaluation', 'Complete test' (highlighted in red), and 'Print preview'.

a. Click on Complete test.

The screenshot shows the EASYCAL™ 5 software interface after clicking 'Complete test'. The title bar reads 'BRAND EASYCAL™ 5'. The header displays 'Order 2021-11-12-14-10-58-DI'. The main area is divided into several sections:

- Test type:** STD - standard values.
- Article:** 704780 - Transferpette® S Adjustable volume: 95 - 1000 µl.
- Order no.:** 2021-11-12-14-10-58-DI.
- Test plan:** 4437.0 - EC-STD-ISO-PIPip-var-LHDIV.
- Order type:** Individual test.
- Tester:** D. Ihrig.
- Lot size:** 1.
- Random sample size:** 1 / 1.
- Status:** Finished.
- Order no.:** [Empty field]
- Customer:** [Empty field]
- OC Number:** [Empty field]
- End customer:** [Empty field]
- Notes on the order:** [Empty field]
- Communication with the customer:** [Empty field]
- Test characteristic:** A table with columns for Test characteristic, Description, Target value, Mean, Tol target, Tol OK, Standard devi, A % target, A %, A % OK, CV % target, CV %, CV % OK, QS, CPK, Min CPK, and CPK OK. It lists 3 test characteristics: 1. TV, 1. TV, and NV.
- Device/calibration slip overview:** A table with columns for Calibration slip type, Internal no., Calibration slip no., Serial number, Article code, Conducted tests, Status, Device status, Tester, and Linked calibration slip. It lists 1 individual test with a status of 'Finished'.
- Buttons:** Includes 'Evaluation/release', 'Print internal log' (highlighted in red), 'Print certificates', 'New calibration slip', 'Back', 'Back to the order overview', and 'Save'.

b. You will now see the order with a summary of the measurements for the respective test characteristics. Click on Print internal log.

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Back

c. You will see the certificate editor in view mode. Here you can print or export the stored report.

d. Exit the certificate editor by clicking Back.

Order 2021-11-12-14-10-58-DI

Test type: STD - standard values Article: 704780 - Transferpette S Adjustable volume: 95 - 1000 µl

Order no.: 2021-11-12-14-10-58-DI Test plan: 4437.0 - EC-STD-ISO-PIP-var-LHDIV

Order type: Individual test Tester: D. Ihrig

Lot size: 1

Random sample size: 1 / 1

Status: Finished

Order no.: Customer: OC-Number: End customer:

Notes on the order: Communication with the customer:

Test characteristic:

Test characteristic	Description	Target va	Mean	Tol target	Tol OK	Standard devi	A% targ	A% OK	CV% targ	CV% OK	CV% OK OS	CPK	Min CPK	CPK OK
1	2. TV	100 µl	100.3 µl	OK	0 µl	0.2 µl	0.1 %	0.1 %	0.3 %	0.0 %	0.0 %	0	0	OK
2	1. TV	500 µl	500.4 µl	OK	0.80 µl	1.6 %	0.1 %	0.1 %	0.6 %	0.10 %	0.10 %	0	0	OK
3	NV	1000 µl	1000.01 µl	OK	0 µl	0.0 %	0.3 %	0.2 %	0.3 %	0.0 %	0.0 %	0	0	OK

Device/calibration slip overview

Calibration slip type	Internal no.	Calibration slip no.	Serial number	Article code	Conducted tests	Status	Device status	Tester	Linked calibration slip
Individual test	13014	ISO / 74	704780	30	Finished			D. Ihrig	

Evaluation/release Print internal log Print certificates New calibration slip

Back Back to the order overview Save

e. Click on Evaluation/release.

Evaluation/release

Current status

Finished

New status

Released

Remarks

☒ Apply to all

Cancel

Save

f. In the window that opens, select Released, and confirm the entry with Saving....

BRAND EASYCAL™ 5

Order 2021-11-12-14-10-58-DI

Test type

STD – standard values

Article

704780 - Transferpette® 5 /adjustable volume; 95 - 1000 µl

Notes on the order

Communication with the customer

Order no.

2021-11-12-14-10-58-DI

Test plan

4437.0 - EC-STD-ISO-PIpip-var-LHDiv

Order type

Individual test

Tester

D. Ihrig

Lot size

1

Random sample size

1 / 1

Order no.

Customer

Status

Released

OC-Number

End customer

Test characteristic

Test characteristic	Description	Target	Mean	Tol target	Tol OK	Standard devi	A% target	A% OK	CV% target	CV% OK	CV% OK	OS	CPK	Min CPK	CPK OK
1	2. IV	100µl	100.3 µl		OK	0 µl	8 %	0 %	OK	3 %	0 %	0 %	OK	0	OK
2	1. IV	500µl	500.4 µl		OK	0.88 µl	1.6 %	0.1 %	OK	0.6 %	0.18 %	0.18 %	OK	0	OK
3	NV	1000µl	1003.01 µl		OK	0 µl	0.8 %	0.3 %	OK	0.3 %	0 %	0 %	OK	0	OK

Device/calibration slip overview

Calibration slip type	Internal no.	Calibration slip no.	Serial number	Article code	Conducted tests	Status	Device status	Tester	Linked calibration slip
Individual test	13014	ISO-74		704780	30	Released		D. Ihrig	

Evaluation/release

Print internal log

Print certificates

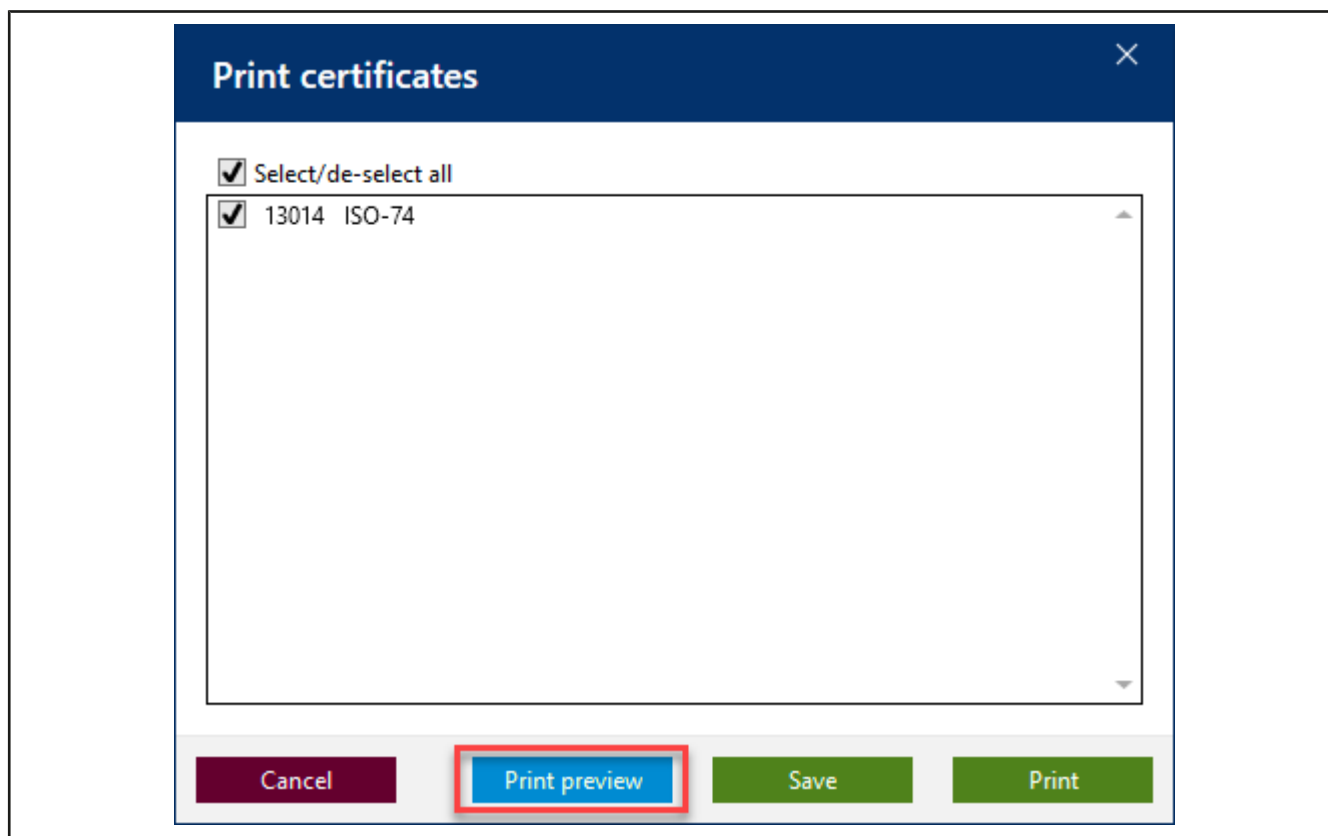
New calibration slip

Back


Back to the order overview

Save

g. Click on Print certificates.

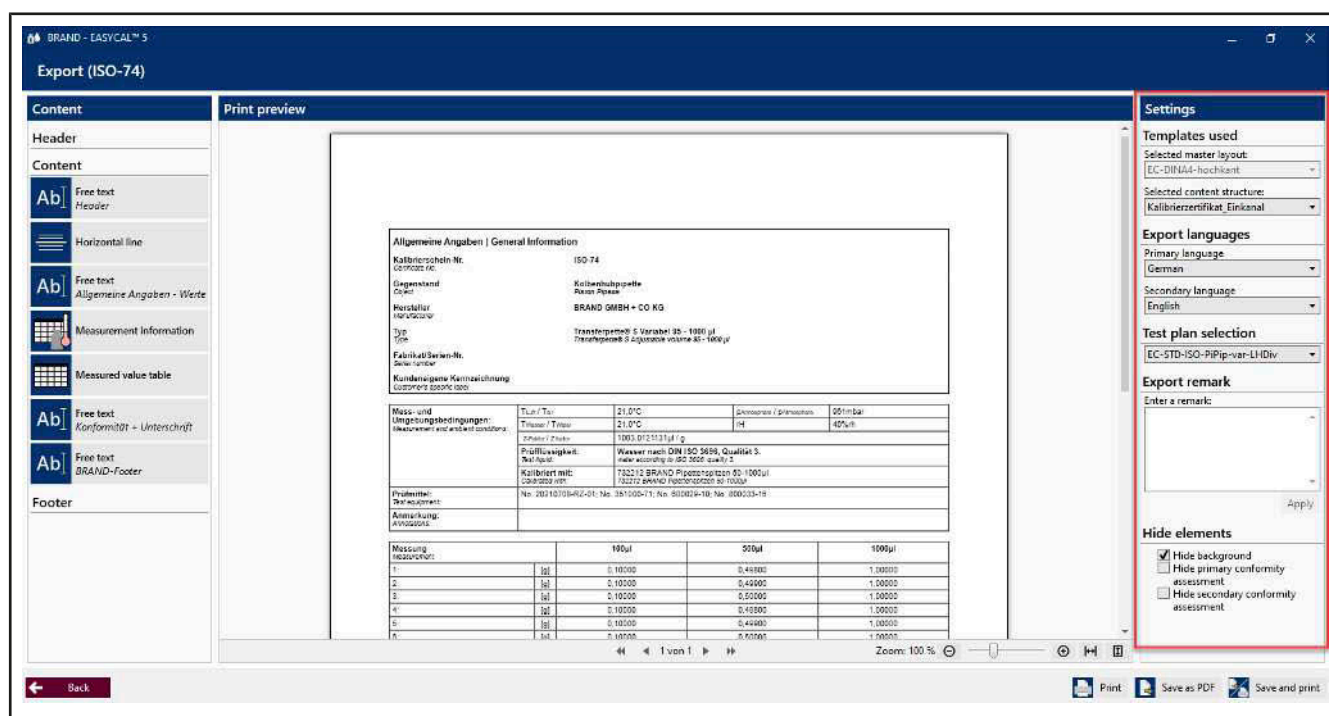
h. Click the check-box, and then click Print preview.

➤ The certificate editor opens in Print preview mode. You can still make settings in the right section and enter export annotations depending on the selected content structure.

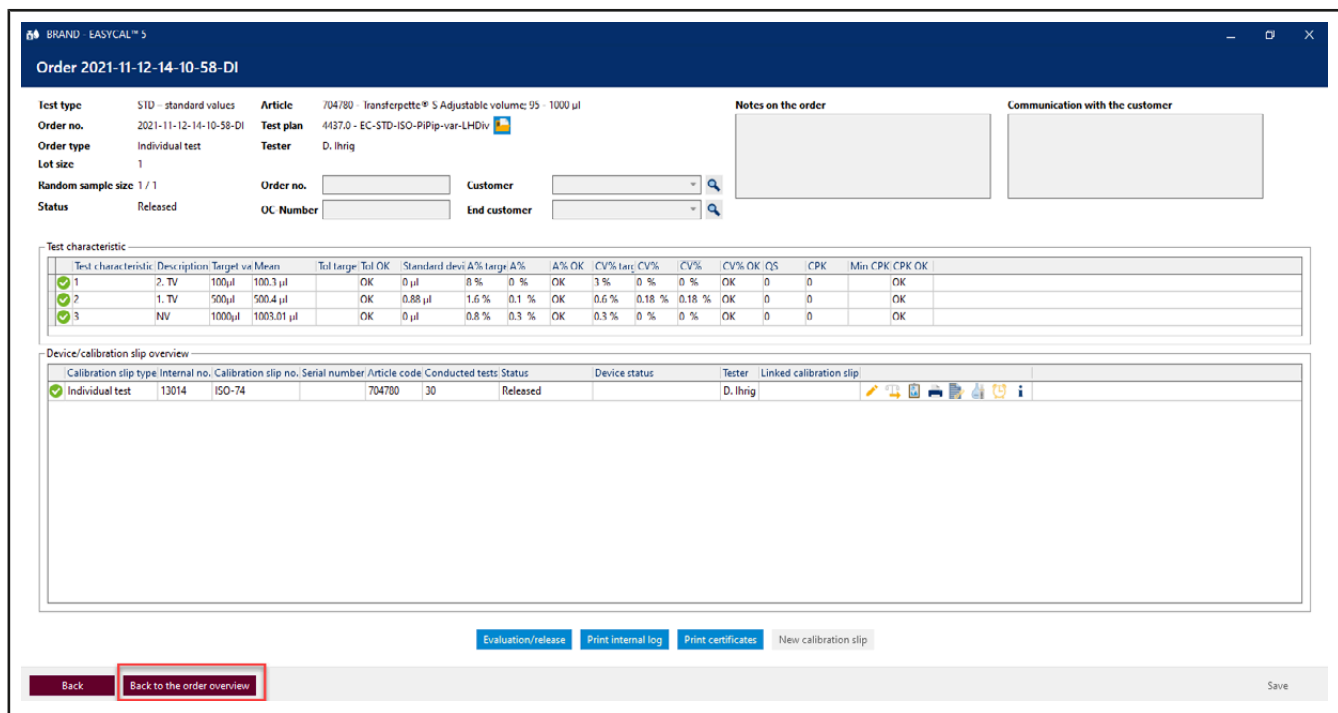
- i. Here you can also export the certificate as a PDF or print a hard copy. By clicking on Back, you leave the Print preview.

- j. Click on Back to the order overview

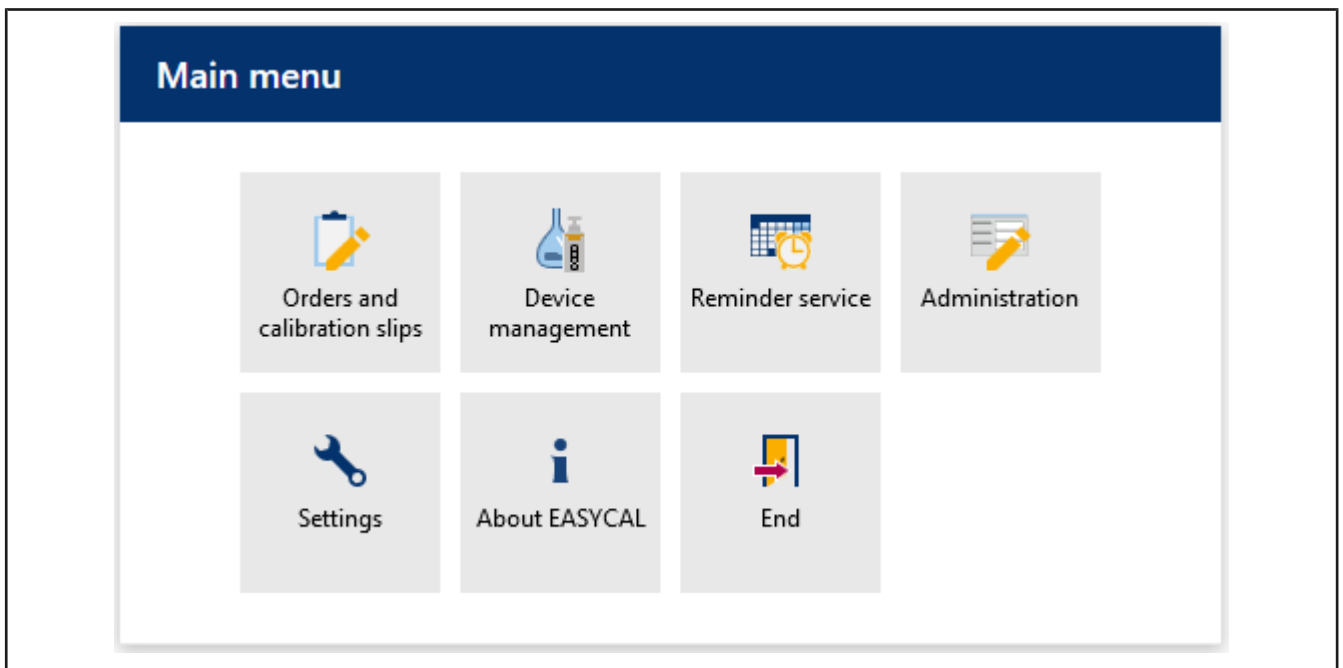- k. From here you can create additional jobs and start tests. By clicking on Back, you can return to the main menu.

➤ This is where this tutorial ends.

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BRAND GMBH + CO KG

Otto-Schott-Str. 25

97877 Wertheim (Germany)

T +49 9342 808 0

F +49 9342 808 98000

info@brand.de

www.brand.de

Do you need more operating manuals and translations?

Please refer to <http://www.brand.de/om> or use the following Quick

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Estado: Completado Número AB: Cliente final: