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Indien:

BRAND Scientific Equipment Pvt. Ltd.
303, 3rd Floor, 'C' Wing, Delphi
Hiranandani Business Park,
Powai

China:

BRAND (Shanghai) Trading Co., Ltd.
Guangqi Culture Plaza
Room 506, Building B
No. 2899, Xietu Road

Mumbai-400 076 (India)
T +91 22 42957790
F +91 22 42957791
info@brand.co.in
www.brand.co.in

Shanghai 200030 (P.R. China)
T +86 21 6422 2318
F +86 21 6422 2268
info@brand.com.cn
www.brand.cn.com

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Contents

1 Introduction	17
1.1 Scope of delivery	17
1.2 Terms of use	17
2 Safety regulations	17
3 Functional and operational components	19
4 Operation	20
4.1 Preparing the burette	20
4.2 Filling the burette	20
4.3 Deaerating the burette	20
4.4 Titrating	20
4.5 Post-titration	21
5 Cleaning	22
6 Recommendation	22
7 Order Information	23
7.1 Accessories	23
7.2 Spares	23
8 Warranty	26
9 Repairs	27
9.1 Sending for repair	27
10 Disposal	28

1 Introduction

1.1 Scope of delivery

BLAUBRAND® or SILBERBRAND model Pellet-type burette with PTFE valve stopcock or PTFE key; models with or without intermediate stopcock (PTFE key), BLAUBRAND® models with quality certificate.

1.2 Terms of use

- Carefully read the operating manual before using the device for the first time.
- The operating manual is part of the device and must be kept in an easily accessible place.
- Be sure to include the operating manual if you transfer possession of this device to a third party.
- You can find up-to-date versions of the operating manual on our website: www.brand.de.

1.2.1 Hazard levels

The following signal words identify possible hazards:

Signal word	Meaning
DANGER	Will lead to serious injury or death.
WARNING	May lead to serious injury or death.
CAUTION	May lead to minor or moderate injuries.
NOTICE	May lead to property damage.

1.2.2 Symbols

Symbol	Meaning
	Danger area

1.2.3 Format

Format	Meaning	Format	Meaning
1. Task	Indicates a task.	>	Indicates a condition.
a., b., c.	Indicates the individual steps of a task.	⇒	Indicates a result.

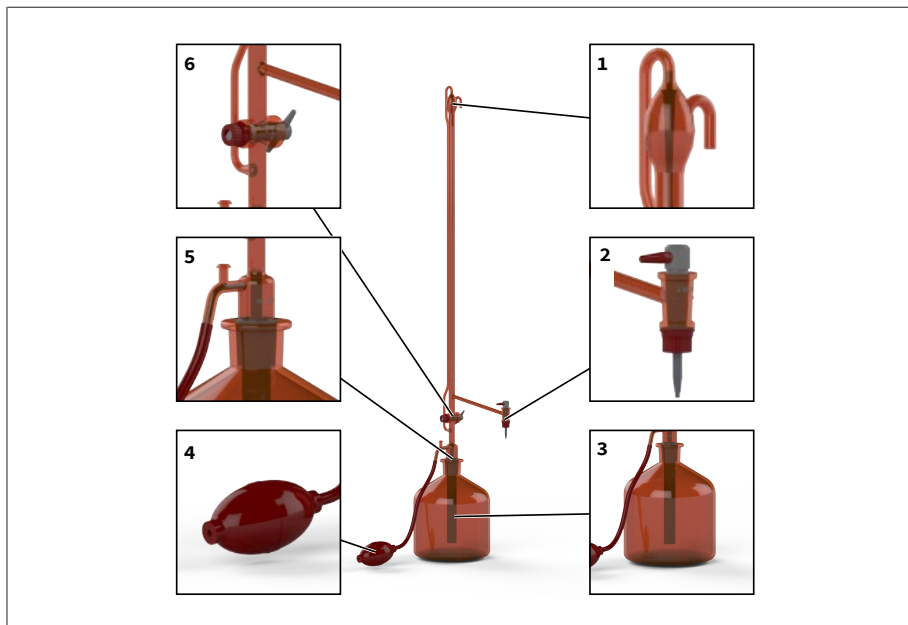
2 Safety regulations

Please read carefully!

The laboratory device can be used in combination with hazardous materials, work processes and equipment. However, the operating manual cannot cover all of the safety issues that may occur in doing so. It is the user's responsibility to ensure compliance with the safety and health regulations and to specify the corresponding restrictions before use.

1. Follow the general instructions for hazard prevention and the safety regulations. E.g., wear protective clothing, eye protection and protective gloves.
2. Follow the instructions given by the reagent manufacturer.
3. Every user must read this operating manual before using the instrument.
4. Always perform work in a manner that does not endanger users or other people.
5. Use the instrument only for titrating and only within the limits of its material resistance.
6. Never use force.
7. To unstick stuck stopcocks or ground joint connections, follow the instructions (see Cleaning, p. 22)
8. Use only original accessories and original replacement parts.

3 Functional and operational components

- | | |
|---|---|
| <p>1 Automatic zero point setting; located beside it is the overflow.</p> <p>3 Supply bottle*</p> <p>5 NS 29/32 ground joint connection, olive and vents</p> | <p>2 Burette stopcock, pictured here is a PTFE key. Other models have a valve stopcock.</p> <p>4 Rubber blowball*</p> <p>6 Intermediate stopcock; located behind it is the filling tube.</p> |
|---|---|

* This item is not included in the scope of delivery.

4 Operation

4.1 Preparing the burette

- a. Check whether the stopcocks are easy to turn.
- b. Lightly grease the ground joint connection to the supply bottle with stopcock grease, or use a PTFE ground joint sleeve.

4.2 Filling the burette

⚠ WARNING

Media leakage

In general, do not fill while the intermediate stopcock is open. If overfilled, media will escape through the overflow at the top of the burette.

- a. Place the burette onto the filled supply bottle.
 - b. Connect the rubber blowball to the side tubing connector.
 - c. Close the intermediate stopcock (model with intermediate stopcock).
 - d. Hold the vent closed with a finger. Squeeze the blowball to fill the burette.
 - e. When the liquid level is approx. 5 mm above the zero mark, lift your finger off the vent.
- ⇒ This ends the process.
- ⇒ The burette will automatically set to zero.

4.3 Deaerating the burette

- a. To remove trapped air, briefly open and close the burette stopcock.
- b. After the first filling operation, a small air bubble may remain on the burette stopcock.
- c. Hold the burette at an angle to remove the air bubble. Tap lightly with your finger against the area where the bubble is located.

4.4 Titrating

Preventing volume errors

- a. Begin titration from the zero point.
- ⇒ This prevents volume errors.
- b. Turn the valve stopcock gently when closing. Excessive force may lead to deformation of the valve spindle, resulting in an unwanted rise of the liquid column in the burette tube (reading error).

Observing wait times

NOTICE

If titration is carried out, the wait time can be omitted since the time required to reach the titration end point is normally longer than the specified wait time.

Class	Wait time in accordance with DIN EN ISO 385
AS	30 s
B	Not applicable

When calibrating Class A burettes, the wait time according to DIN EN ISO 385 must be adhered to.

4.5 Post-titration

On burettes with an intermediate stopcock, the content of the burette can be allowed to flow back into the supply bottle. It should be noted that liquid residue on the wall may partially evaporate.

Preventing volume errors

To prevent volume errors, fill and empty the burette one to two times before the next titration.

Storing the burette

Clean the burette if you will not be using it for several days. If it is not cleaned, stopcocks and ground joint connections may get stuck.

5 Cleaning

CAUTION

Risk of injury from broken glass!

Never use excessive force to unstick stuck stopcocks or ground joint connections.

Cleaning

If the burette does not discharge without dripping or if it sets the zero point too low, the burette must be cleaned. To clean, fill the burette with a mild alkaline cleaner and rinse it out with water several times after the prescribed soaking time.

Unsticking stuck parts

- a. Completely immerse stuck stopcocks or ground joint connections in a cleaning solution and allow to soak overnight.
- b. Carefully unstick stuck stopcocks or ground joint connections while wearing suitable safety gloves.
- c. If unsuccessful, send the instrument in for repair.

6 Recommendation

We recommend using a PTFE ground joint sleeve for the ground joint connection to the supply bottle, as these do not require grease for sealing.

7 Order Information

7.1 Accessories

Rubber blowball

Description	Order No.
Rubber blowball	23400

Ground joint sleeve

Description	Packaging unit	Order No.
PTFE ground joint sleeve	10 piece	51422

Ground joint socket

Description	Order No.
PTFE ground joint socket NS 29/32	51467

Joint clip NS 29/32

Description	Order No.
Stainless steel	55618
POM (red)	55646

Supply bottle 2000 ml, NS 29/32

Description	Order No.
Boro 3.3, clear	23310
Soda-lime glass, clear	126965
Boro 3.3, brown	23320
Soda-lime glass, amber	127065

7.2 Spares

7.2.1 Clear glass models

Burette stopcock, PTFE (10, 25, 50 ml)

Description	Order No.
Burette stopcock, PTFE (10, 25, 50 ml)	82170

Burette stopcock, PTFE with housing and tip

Description	Order No.
Burette stopcock, PTFE with housing and tip 10 ml, Boro 3.3, clear	82120
Burette stopcock, PTFE with housing and tip 25 ml, Boro 3.3, clear	82121
Burette stopcock, PTFE with housing and tip 50 ml, Boro 3.3, clear	82122

Intermediate stopcock, PTFE key 4NS/19

Description	Order No.
Intermediate stopcock, PTFE key 4 NS/19	81265

7.2.2 Amber glass models

Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip (25, 50 ml)

Description	Order No.
Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip (25, 50 ml)	81167

Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip and housing

Description	Order No.
Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip and housing (25, 50 ml, Boro 3.3 amber)	81927

Intermediate stopcock, PTFE key 4NS/14

Description	Order No.
Intermediate stopcock, PTFE key	81270

8 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the instrument or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

USA and Canada:

Find more warranty information on www.brandtech.com.

9 Repairs

9.1 Sending for repair

NOTICE

Transporting hazardous materials without approval is prohibited by law.

Clean the instrument thoroughly and decontaminate!

- When returning products, please enclose a general description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Only send the device without a battery installed.
- Shipment is at the risk and the cost of the sender.

Outside USA and Canada

Complete the “Declaration on Absence of Health Hazards” and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from www.brand.de.

Outside USA and Canada

Please clarify the requirements for the return delivery with BrandTech Scientific, Inc **before** sending the instrument in for service.

Return only cleaned and decontaminated instruments to the address provided with the Return Authorization Number. Place the Return Authorization number so that it is clearly visible on the outside of the package.

Contact addresses

Germany:

BRAND GMBH + CO KG
Otto-Schott-Str. 25
97877 Wertheim (Germany)
T +49 9342 808 0
F +49 9342 808 98000
info@brand.de
www.brand.de

USA and Canada:

BrandTech® Scientific, Inc.
11 Bokum Road
Essex, CT 06426-1506 (USA)
T +1-860-767 2562
F +1 - 860 - 767 2563
info@brandtech.com
www.brandtech.com

India:

BRAND Scientific Equipment Pvt. Ltd.
303, 3rd Floor, 'C' Wing, Delhi
Hiranandani Business Park,
Powai

China:

BRAND (Shanghai) Trading Co., Ltd.
Guangqi Culture Plaza
Room 506, Building B
No. 2899, Xietu Road

Mumbai-400 076 (India)
T +91 22 42957790
F +91 22 42957791
info@brand.co.in
www.brand.co.in

Shanghai 200030 (P.R. China)
T +86 21 6422 2318
F +86 21 6422 2268
info@brand.com.cn
www.brand.cn.com

10 Disposal

Before disposal, observe the relevant national disposal regulations and ensure that the product is disposed of properly.

Subject to technical changes, errors, and misprints.

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BrandTech[®] Scientific, Inc.
BRAND Scientific Equipment Pvt.Ltd.