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Table of Contents

1	Introduction	23
1.1	Use the instructions for use.....	23
1.2	Product markings.....	24
2	Safety	24
2.1	Function.....	25
2.2	Limitations of use.....	25
2.3	Operating exclusions	25
2.4	Storage Conditions	25
3	Functional and operating elements	26
4	Pipetting	27
5	Autoclaving	30
6	Cleaning	31
6.1	Replace filter and clean adapter	31
6.2	Replace valve system.....	34
6.3	Check tightness.....	34
7	Malfunction – what to do?	36
8	Ordering Information	36
9	Spare parts.....	36
10	Repairs.....	38
10.1	Sending for repair	38
11	Warranty	39
12	Disposal	39

1 Introduction

1.1 Use the instructions for use

- Carefully read the operating manual before using the device for the first time.
- Keep the user manual easily accessible. It is part of the device.
- Be sure to include the operating manual if you transfer possession of this device to a third party.
- You can find up-to-date versions of the operating manual on our website: www.brand.de.

1.1.1 Signal words and their meaning

Signal word	Meaning
⚠ WARNING or ⚠WARNING! ...	WARNING indicates a dangerous situation that, if not avoided, could result in death or serious injury.
⚠ CAUTION or ⚠CAUTION! ...	CAUTION indicates a hazardous situation that, if not avoided, could result in moderate or minor injury.
NOTICE or NOTICE ...	NOTE is used to address actions that are not related to physical injury. Example: Possible property damage.

1.1.2 Presentation of descriptions of actions

Format	Meaning
1. Task	Indicates a task.
a., b., c.	Indicates the individual steps of a task.
>	Indicates a prerequisite for a task.
⇒	Indicates a result of a completed task.

1.2 Product markings

Layout	Meaning
	General warning symbol
	Follow the operating instructions
XXZ	Batch number

2 Safety

Please read carefully!

The laboratory device macro can be used in combination with hazardous materials, work processes and equipment. However, the operating manual cannot cover all of the safety issues that may occur in doing so. It is the user's responsibility to ensure compliance with the safety and health regulations and to specify the corresponding restrictions before use.

1. Every user must read and observe this operating manual before using the device.
2. Follow the general hazard instructions and safety regulations (e.g., wear protective clothing, eye protection, and protective gloves).
3. Follow the instructions given by the reagent manufacturer.
4. Use the device only for pipetting liquids and only within the defined limits of use. Comply with the operating exclusions; see Operating exclusions, p. 25 . In case of doubt, contact the manufacturer or dealer.
5. Always perform work in a manner that does not endanger the user or other people. Avoid splattering. Use only suitable vessels.
6. Never use force.

7. Use only original accessories and original replacement parts. Do not make any technical modifications. Do not disassemble the device further than described in the user manual!
8. Always check that the device is in proper working condition before use. If the device malfunctions, stop pipetting immediately and refer to the section Malfunction – what to do?, p. 36 . If necessary, contact the manufacturer.

2.1 Function

The device serves as an aid for pipetting liquids with graduated and volumetric pipettes made of glass or plastic in the volume range from 0.1 ml to 200 ml and with a suction tube outer diameter < 9.2 mm. When the device is handled correctly, the pipetted liquid comes into contact with the pipette only.

2.2 Limitations of use

This device is intended for pipetting liquids within the following physical limitations:

- from 10 to 40°C (from 50 to 104°F) (device and reagent)
- vapor pressure up to 500 mbar. Absorb slowly above 300 mbar in order to prevent the liquid from boiling.
- Density up to 2.4 g/cm³

2.3 Operating exclusions

The device must not be used for liquids with vapors that can attack the materials silicone or PP.

The device is not suitable for use with Pasteur pipettes.


2.4 Storage Conditions

Store the unit and accessories in a cool and dry place in cleaned condition only.

Storage temperature from -20 to 50°C (-4 to 122°F).

Relative humidity: 5 % to 95 %.

3 Functional and operating elements

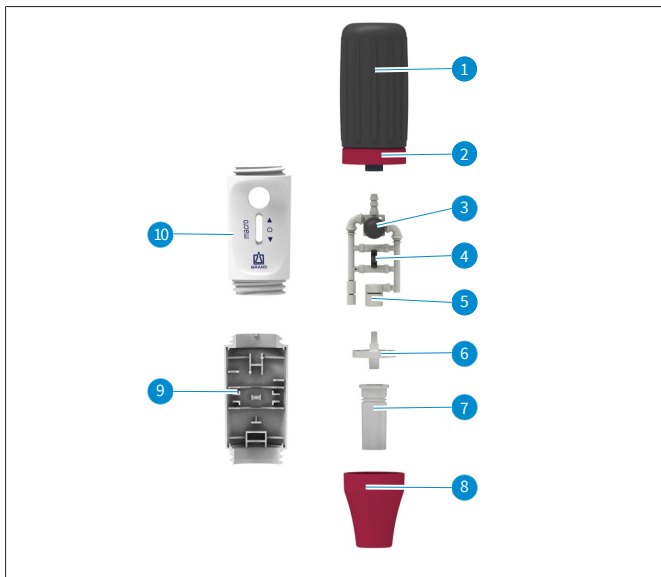1	Suction bellows (SI)	6	Membrane filter (PP/PTFE)
2	Screw ring (PP)	7	Adapter (SI)
3	Blow-out bubble (SI)	8	Adapter housing (PP)
4	Pipetting lever	9	Housing (PP)
5	Valve system (PP, SI)	10	Housing (PP)

Abbreviation	Material
PP	Polypropylene
PTFE	Polytetrafluoroethylene
SI	Silicone

4 Pipetting

1. Vent the suction bellows

1.

- a. Press the suction bellows together before attaching the pipette.

2. Attach the pipette

2.

- a. Always hold the pipette as close as possible to the upper end and push it carefully into the adapter until it is firmly in place.

⚠WARNING!
Risk of injury!


Never use force! Especially with thin pipettes, there is a risk of glass breakage!

Ensure that the pipette is firmly in place.

- b. Always hold the instrument with the pipette inserted vertically, tip down.

3. Fill the pipette

3.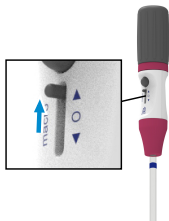

- a. Immerse pipette tip into the test liquid.
- b. Carefully push the pipetting lever upwards, and fill the pipette slightly above the desired mark.
The further the lever is moved upwards, the stronger the suction effect and the faster the meniscus rises!

NOTICE Make sure that no liquid gets into the device! The function of the membrane filter is impaired, and the suction capacity decreases. In this case, replace membrane filter; see Replace filter and clean adapter, p. 32 .



NOTICE With large pipettes (> 50 ml), the vacuum in the suction bellows is not sufficient to draw up the entire amount of liquid at once.
In this case:
Compress the suction bellows again and continue drawing up liquid.

4. Set the volume

- a. If necessary, wipe the tip of the pipette with a suitable lint-free cloth.
- b. Press the pipetting lever down gently until the meniscus is precisely adjusted.

5. Empty the pipette

5.

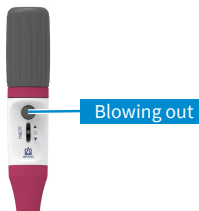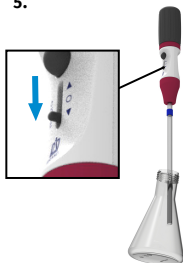- a. Hold collecting vessel at an angle. Place the pipette tip against the inner wall and press the pipetting lever downward.

When using pipettes with a waiting time (imprint e.g. 'Ex + 5 s'):

As soon as the meniscus in the pipette comes to a standstill, observe the waiting time indicated on the pipette (in this case, 5 s). Pull up the tip of the pipette a few millimeters along the wall of the vessel.

When using blow-out pipettes from other manufacturers:

As soon as the meniscus in the pipette tip has come to a standstill, press the small bellows once. Pull up the tip of the pipette a few millimeters along the wall of the vessel.

6. After pipetting

6.

- a. Grip the pipette as high up as possible, and pull it out of the adapter.

5 Autoclaving

NOTICE

Check the effectiveness of autoclaving yourself!

Maximum safety is achieved through vacuum sterilization. We recommend using sterilization bags.

- a. Clean the instrument thoroughly prior to autoclaving, see Cleaning, p. 31. Clean the housing parts with a damp cloth.
- b. Unscrew the suction bellows (1) prior to autoclaving.
- c. Pack the instrument, including the membrane filter, in a sterilization bag, observing any applicable packaging regulations.

NOTICE Not every membrane filter is suitable! Use only original accessories. The device comes

standard with filters of 3 µm pore width. This membrane filter can be autoclaved up to five times at 121°C.

- d. Autoclave the instrument. After autoclaving and before use, allow the instrument to cool down.

Temperature	121 °C
Pressure	2 bar
Holding time in autoclaves	15 min

6 Cleaning

Occasionally clean the housing parts with a damp cloth. When used properly, the instrument is maintenance-free.

6.1 Replace filter and clean adapter

NOTICE It is only necessary to replace and clean the filter if a pipette has been overfilled or suction performance is impaired.

⚠WARNING! Contamination and infection risk

Wear protective gloves and eye protection.

b.

- a. Pull out pipette.
- b. Unscrew the adapter housing.
- c. Remove defective membrane filter.
- d. Press the adapter upwards out of the adapter housing.

e.

- e. For cleaning, rinse the adapter carefully with a spray bottle filled with a suitable cleaning liquid (e.g., water). Then blow out and dry.

f.

- f. First press the new membrane filter (1) into the adapter (2) with the thicker connection facing downwards (the thinner, conical side facing upwards).
- g. Tighten the adapter housing.


6.2 Replace valve system

a.

- a. Unscrew suction bellows and adapter housing.

b.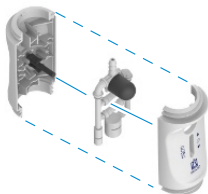

- b. Remove the upper part of the housing for replacement. Carefully remove and replace the valve system.
- c. Reassemble the unit in reverse order and check for leaks.

6.3 Check tightness

After adjusting the meniscus, no liquid should drip from the pipette.

- a. Adjust the meniscus.
- b. Hold the instrument vertically.
- c. If a drop forms after a short waiting time (30 s), follow the remedy under Malfunction - What to do?, p. 36

7 Malfunction – what to do?

Fault	Possible cause	What to do
Suction capacity impaired	Filter dirty or valve system defective	Replace filter, Replace filter and clean adapter, p. 31 .
Pipette drips	Filter not placed correctly or valve system defective	Insert filter correctly, or replace valve system; see Replace filter and clean adapter, p. 31 .
Pipette does not hold	Silicone adapter damaged	Unscrew the lower part of the device, and replace the silicone adapter
Suction bellows ventilates independently	Suction bellows or valve system defective	Replace upper screw connection with suction bellows or valve system

NOTICE A visual inspection of the device is often sufficient to detect a defective part. This can then be easily replaced with a spare part. The advantages of self-repair are obvious: cost and time savings.

8 Ordering Information

macro pipette aid

Description	Color*	Order No.
macro pipetting aid, including. Replacement membrane filter (3 µm) and operating instructions.	gray	26200
	Green	26201
	Blue	26202
	Magenta	26203

* The colors refer to the adapter housing and screw ring.

9 Spare parts

Adapter housing

Color*	Length (mm)	Material	Order No.
gray	49	PP	26220

Color*	Length (mm)	Material	Order No.
Green	49	PP	26223
Blue	49	PP	26221
Magenta	49	PP	26222

Suction bellows with screw ring

Color*	Order No.
gray	26225
Green	26228
Blue	26226
Magenta	26227

* The colors refer to the adapter housing and screw ring.

Replacement membrane filter (3 µm)

Packaging unit	Order No.
1 pc in blister packaging	26052

Replacement membrane filter (3 µm)

Packaging unit	Order No.
10 pc in blister packaging	26056

Silicone adapter

Length (mm)	Order No.
44	26146

Valve system

Description	Order No.
Valve system	26128

10 Repairs

10.1 Sending for repair

NOTICE

Transporting hazardous materials without approval is prohibited by law.

Clean the instrument thoroughly and decontaminate!

- When returning products, please enclose a general description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Shipment is at the risk and the cost of the sender.

Outside USA and Canada

Complete the “Declaration on Absence of Health Hazards” and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from www.brand.de.

Within USA and Canada

Please clarify the requirements for the return delivery with BrandTech Scientific, Inc **before** sending the instrument in for service.

Return only cleaned and decontaminated instruments to the address provided with the Return Authorization Number. Place the Return Authorization number so that it is clearly visible on the outside of the package.

Contact addresses

Germany:

BRAND GMBH + CO KG
Otto-Schott-Straße 25
97877 Wertheim (Germany)
T +49 9342 808 0

USA and Canada:

BrandTech® Scientific, Inc.
11 Bokum Road
Essex, CT 06426-1506 (USA)
T +1-860-767 2562

F +49 9342 808 98000
info@brand.de
www.brand.de

India:

BRAND Scientific Equipment Pvt. Ltd.
303, 3rd Floor, 'C' Wing, Delphi
Hiranandani Business Park,
Powai
Mumbai-400 076 (India)
T +91 22 42957790
F +91 22 42957791
info@brand.co.in
www.brand.co.in

F +1 - 860 - 767 2563
info@brandtech.com
www.brandtech.com

China:

BRAND (Shanghai) Trading Co., Ltd.
Rm 201-202, North Tower,
No. 199 Kaibin Rd, Xuhui District, Shanghai
Shanghai 200030 (P.R. China)
T +86 21 6422 2318
F +86 21 6422 2268
info@brand.com.cn
www.brand.cn.com

11 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the instrument or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

USA and Canada:

Find more warranty information on www.brandtech.com.

12 Disposal

Before disposal, observe the relevant national disposal regulations and ensure that the product is disposed of properly.

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