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# Contents

<b>1 Terms of use .....</b>	<b>26</b>
1.1    Hazard levels .....	26
1.2    Operating manual symbols .....	26
1.3    Format .....	26
<b>2 Safety .....</b>	<b>28</b>
2.1    Function.....	29
2.2    Operating exclusions .....	29
2.3    Storage Conditions .....	29
<b>3 Functional and operating elements .....</b>	<b>30</b>
<b>4 Pipetting .....</b>	<b>31</b>
<b>5 Cleaning.....</b>	<b>36</b>
5.1    Cleaning the suction system .....	36
<b>6 Storage .....</b>	<b>37</b>
<b>7 Replacing adapter and suction tube.....</b>	<b>38</b>
<b>8 Troubleshooting .....</b>	<b>39</b>
<b>9 Product markings .....</b>	<b>40</b>
<b>10 Order Information.....</b>	<b>41</b>
<b>11 Spares.....</b>	<b>42</b>
<b>12 Repairs.....</b>	<b>43</b>
12.1    Sending for repair .....	43
<b>13 Warranty .....</b>	<b>45</b>
<b>14 Disposal .....</b>	<b>46</b>

# 1 Terms of use

- Carefully read the operating manual before using the device for the first time.
- The operating manual is part of the device and must be kept in an easily accessible place.
- Be sure to include the operating manual if you transfer possession of this device to a third party.
- You can find up-to-date versions of the operating manual on our website: [www.brand.de](http://www.brand.de).

## 1.1 Hazard levels

The following signal words identify possible hazards:

Signal word	Meaning
DANGER	Will lead to serious injury or death.
WARNING	May lead to serious injury or death.
CAUTION	May lead to minor or moderate injuries.
NOTICE	May lead to property damage.

## 1.2 Operating manual symbols

Symbol	Meaning	Symbol	Meaning
	Danger area		Wear protective gloves
	Wear eye protection		

## 1.3 Format

Format	Meaning	Format	Meaning
<b>1. Task</b>	Indicates a task.	>	Indicates a condition.

<b>Format</b>	<b>Meaning</b>	<b>Format</b>	<b>Meaning</b>
a., b., c.	Indicates the individual steps of a task.	⇒	Indicates a result.

## 2 Safety

### Please read carefully!

The instrument micro-classic can be used in combination with hazardous materials, work processes and equipment. However, the operating manual cannot cover all of the safety issues that may occur in doing so. It is the user's responsibility to ensure compliance with the safety and health regulations and to specify the corresponding restrictions before use.

1. Every user must read and observe this user manual before using the device.
2. Follow the general hazard instructions and safety regulations (e.g., wear protective clothing, eye protection, and protective gloves).
3. Observe all specifications provided by reagent manufacturers.
4. Use the instrument only for pipetting liquids and only within the defined limits of use. If in doubt, contact the manufacturer or supplier.
5. Always perform work in a manner that does not endanger yourself or other people. Avoid splattering. Only use suitable vessels.
6. Always attach only unfilled pipettes to prevent contamination and infection risks.
7. Make sure that no liquid gets into the device.
8. Never use force.
9. Use only original accessories and original replacement parts. Do not make any technical modifications. Do not dismantle the instrument any further than is described in the operating manual!
10. Always check that the instrument is in proper working condition before use. If an instrument fault is signaled, stop pipetting immediately and follow the instructions in the section Troubleshooting, p. 39 . Contact the manufacturer, if necessary.

## 2.1 Function

The device serves as a pipetting aid. It reduces the risk of infection when pipetting liquids with disposable micropipettes with a ring mark, blood diluting pipettes and other pipettes with a shaft diameter < 5.5 mm and a volume of up to 1 ml.

When the device is handled correctly, the pipetted liquid comes into contact only with the pipette.

## 2.2 Operating exclusions

- Do not use the device with end-to-end micropipettes.

Liquid vapors may penetrate into the adapter (PVC) and into the suction tube (SI). Therefore, never use the device for liquids that corrode PVC or silicone.


## 2.3 Storage Conditions

Store the unit and accessories in a cool and dry place in cleaned condition only.

Storage temperature from -20 to 50°C (-4 to 122°F).

Relative humidity: 5 % to 95 %.

## 3 Functional and operating elements

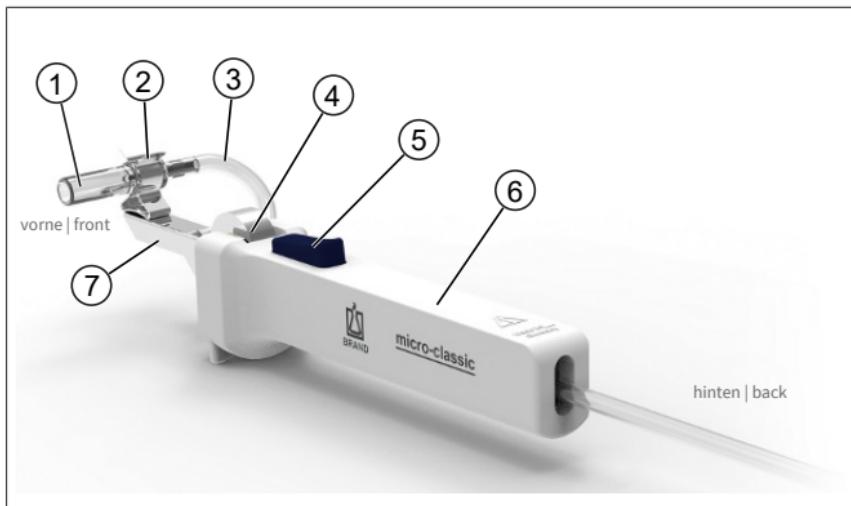<b>1</b>	Adapter	<b>5</b>	Venting button
<b>2</b>	Spring clip	<b>6</b>	Handle
<b>3</b>	Suction tube	<b>7</b>	Adapter support
<b>4</b>	Adjustment wheel		—

# 4 Pipetting

## Correct hand position before pipetting

- Hold the instrument horizontally.
- Push back the O-Ring from the venting button.

## Attaching the pipette


### ⚠ WARNING

Never attach filled pipettes:

- > Risk of contamination and infection.
- > Wear protective gloves.

a.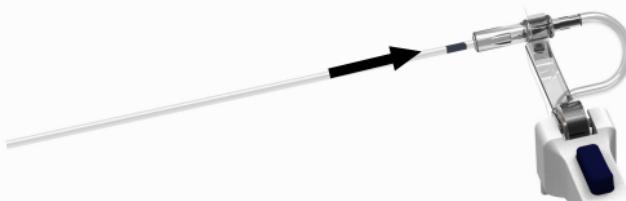


- Hold the pipette at the color code (short end up) to avoid glass breakage when fitting the pipette.

- b. Carefully insert the pipette into the adapter until a slight resistance can be felt. Never use force: Risk of damaging the adapter and pipette, hence a risk of injury.
- c. If the pipette breaks: Remove glass splinters by means of forceps. Check adapter for damages and, if necessary, replace adapter, see Replacing adapter and suction tube, p. 38.

## Aspirating liquid

b.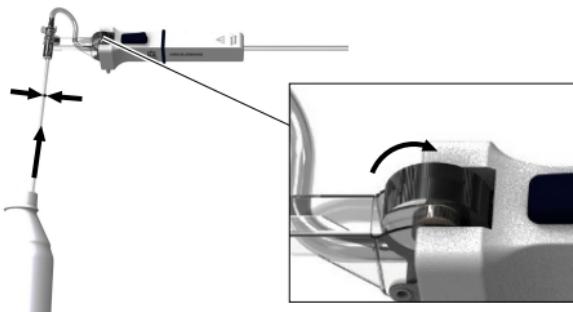

Allow the tube a length of approx. 100 mm between adapter and thumb wheel to keep the dead volume low.

- a. Immerse Pipette tip into the liquid.
- b. Slowly turn the thumb wheel towards the ejector key until the liquid has exactly reached the appropriate mark.

### NOTICE


Always ensure that the liquid does not contaminate adapter and suction tube. If they are contaminated, they must immediately be cleaned or disinfected or, if necessary, be replaced.

- c. Take the pipette out of the liquid.

- d. Carefully wipe the outside of the pipette with a lint-free cloth or swab. Take care not to touch the capillary orifice to avoid liquid being sucked out.

### Discharge liquid (pipettes calibrated to contain "In")


b.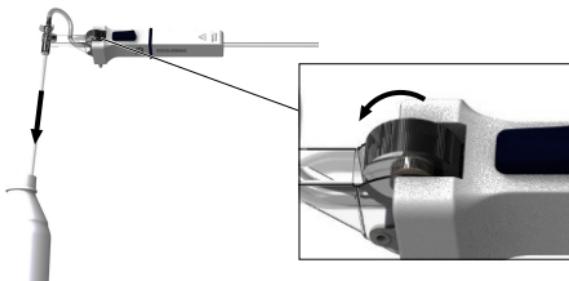

- Immerse the pipette into the diluting agent.
- Turn the thumb wheel towards the adapter until the liquid has been discharged.
- Rinse the pipette by drawing in and discharging the diluting agent: Carefully draw in the diluting agent up to the correct level of the liquid to be measured.
- Repeat this procedure at least 3 times.
- Lightly wipe the pipette against the wall of the vessel.

## Discharge liquid (pipettes calibrated to contain "Ex")

b.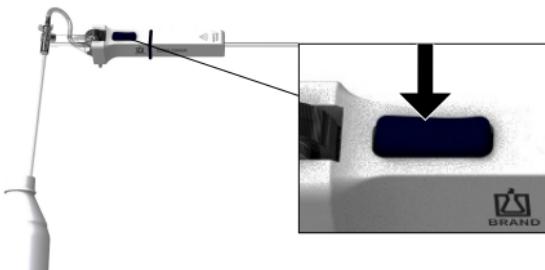

- a. Hold the instrument upright, and place the pipette tip against the wall of a vessel (inclined position of the vessel).
- b. Press the venting button until the liquid has run out.
- c. As soon as the meniscus comes to a standstill in the pipette tip, the waiting time (if applicable) starts.
- d. Then take the pipette out of the vessel by wiping it against the wall of the vessel. During this procedure the residue of liquid partly runs out.


### Removing the pipette

#### **WARNING**

Do not touch the contaminated area of the pipette. Wear protective gloves.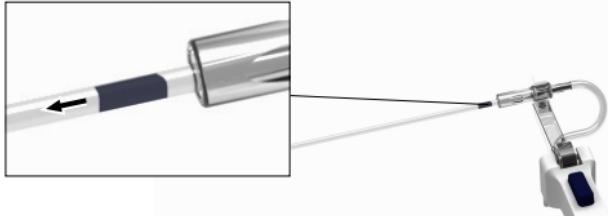

**a.**

- a. Grip the pipette just below the adapter.
- b. Carefully pull the pipette out of the adapter.
- c. Dispose the micropipettes according to the relevant provisions. Clean reusable micropipettes according to the instructions.

### Releasing the suction tube

After pipetting always release the suction tube by pushing the O-ring on the venting button.

# 5 Cleaning

Occasionally clean the housing parts with a damp cloth. When used properly, the instrument is maintenance-free.

If the adapter and the suction tube have been contaminated by liquid, these parts must be cleaned or disinfected immediately and replaced if necessary

## 5.1 Cleaning the suction system

### ⚠ WARNING

#### Risk of contamination and infection

Wear protective gloves and goggles.

### NOTICE

The adapter and suction tube are autoclavable at 121°C (2 bar) and a holding time of at least 15 minutes, in accordance with DIN EN 285. It is the user's responsibility to ensure effective autoclaving.

- a. Remove the adapter and tube from the device and place them in a cleaner or disinfectant solution.
- b. After this, allow to dry.
- c. Autoclave adapter and suction tube if necessary.
- d. Assemble the device, see. Replacing adapter and suction tube, p. 38
- e. Check that all controls are functioning properly.

## 6 Storage

Only store instrument with completely released suction tube, see  
Pipetting, p. 31 ‘Releasing the suction tube’

## 7 Replacing adapter and suction tube

- a. Pull off the adapter from the adapter support.
- b. Upon pressing the venting button the suction tube can be pull out forward.
- c. Insert the new suction tube with adapter between thumb wheel and roller while pressing the venting button.
- d. Clip the adapter into the adapter support. Simultaneously exert counterpressure at the bottom of the spring clip to avoid damage of the adapter support.

## 8 Troubleshooting

Problem	Possible cause	Corrective action
Pipette broken off in adapter.	Pipette was damaged or pushed into the adapter with too much force.	Remove glass fragments with tweezers; replace adapter with suction hose if necessary, see Replacing adapter and suction tube, p. 38 .
Instrument suction system not working/ pipette dripping.	Adapter or suction tube damaged.	Replace adapter with suction tube, see Replacing adapter and suction tube, p. 38 .
Liquid in adapter or suction tube.	Adjustment wheel turned too quickly.	Clean or disinfect adapter and suction tube and replace if necessary, see Cleaning, p. 36 .
Pipette not held in adapter.	Adapter is too moist or greasy.	Clean or disinfect adapter and allow to dry; replace if necessary, see Cleaning, p. 36 .

### NOTICE

A visual inspection of the device is often sufficient to detect a defective part. This can then be easily replaced with a spare part. The advantages of self-repair are obvious: Cost and time savings.

## 9 Product markings

Symbol or number	Meaning
	General warning symbol
	Follow the operating instructions
XXZXXXXX	Serial number

# 10 Order Information

## micro-classic

Description	Order No.
micro-classic complete with suction tube	259 00

## 11 Spares

### Spare adapter with suction tube

Description	Order No.
Spare adapter with suction tube	259 31

# 12 Repairs

## 12.1 Sending for repair

### NOTICE

Transporting of hazardous materials without a permit is a violation of federal law.

### Clean the instrument thoroughly and decontaminate!

- When returning products, please enclose a general description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Shipment is at the risk and the cost of the sender.

### Outside USA and Canada

Complete the “Declaration on Absence of Health Hazards” and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from [www.brand.de](http://www.brand.de).

### Outside USA and Canada

Please clarify the requirements for the return delivery with BrandTech Scientific, Inc **before** sending the instrument in for service.

Return only cleaned and decontaminated instruments to the address provided with the Return Authorization Number. Place the Return Authorization number so that it is clearly visible on the outside of the package.

### Contact addresses

**Germany:**

997782

**USA and Canada:**

Operating manual

**43**

BRAND GMBH + CO KG  
Otto-Schott-Straße 25  
97877 Wertheim (Germany)  
T +49 9342 808 0  
F +49 9342 808 98000  
info@brand.de  
www.brand.de

BrandTech® Scientific, Inc.  
11 Bokum Road  
Essex, CT 06426-1506 (USA)  
T +1-860-767 2562  
F +1 - 860 - 767 2563  
info@brandtech.com  
www.brandtech.com

**India:**

BRAND Scientific Equipment Pvt. Ltd.  
303, 3rd Floor, 'C' Wing, Delphi  
Hiranandani Business Park,  
Powai  
Mumbai-400 076 (India)  
T +91 22 42957790  
F +91 22 42957791  
info@brand.co.in  
www.brand.co.in

**China:**

BRAND (Shanghai) Trading Co., Ltd.  
Rm 201-202, North Tower,  
No. 199 Kaibin Rd, Xuhui District, Shanghai  
Shanghai 200030 (P.R. China)  
T +86 21 6422 2318  
F +86 21 6422 2268  
info@brand.com.cn  
www.brand.cn.com

## 13 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the device or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

### **USA and Canada:**

Find more warranty information on [www.brandtech.com](http://www.brandtech.com).

## 14 Disposal

Before disposal, observe the relevant national disposal regulations, and ensure that the product is disposed of properly.

Subject to technical changes, errors, and misprints.

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