

Dispenser selection chart

| Reagent | seripettor® | seripettor® pro |
|-----------------------------|-------------|--------------------|
| Acetaldehyde | | + |
| Acetic acid, 5% | + | + |
| Acetic acid, 96% | | + |
| Acetic acid (glacial), 100% | | + |
| Acetone | | + |
| Acetonitrile | | + |
| Acetophenone | + | |
| Acetylacetone | + | + |
| Acrylic acid | | + |
| Acrylonitrile | | + |
| Adipic acid | + | + |
| Agar (60 °C) | + | |
| Allyl alcohol | + | + |
| Aluminium chloride | + | + |
| Amino acids | + | + |
| Ammonia, 30% | + | + |
| Ammonium chloride | + | + |
| Ammonium fluoride | + | + |
| Ammonium sulfate | + | + |
| Amyl alcohol (Pentanol) | + | + |
| n-Amyl acetate | | + |
| Aniline | | + |
| Barium chloride | + | + |
| Benzaldehyde | | + |
| Benzyl alcohol | | + |
| Benzylamine | | + |
| Benzylchloride | | + |
| Boric acid, 10% | + | + |
| Butanediol | + | + |
| 1-Butanol | | + |
| Butylamine | | + |
| n-Butyl acetate | | + |
| Calcium carbonate | + | + |
| Calcium chloride | + | + |

| Reagent | seripettor® | seripettor® pro |
|----------------------------|-------------|--------------------|
| Calcium hydroxide | + | + |
| Calcium hypochlorite | | + |
| Chloroacetaldehyde, 45% | | + |
| Chloroacetic acid | | + |
| Chromic acid, 50% | | + |
| Copper sulfate | + | + |
| Cumene (Isopropyl benzene) | | + |
| Diethylene glycol | + | + |
| Dimethyl sulfoxide (DMSO) | | + |
| Dimethylaniline | | + |
| Ethanol | + | + |
| Formaldehyde, 40% | + | + |
| Formamide | + | + |
| Formic acid, 100% | | + |
| Glycerol | + | + |
| Glycol (Ethylene glycol) | + | + |
| Glycolic acid, 50% | + | + |
| Hexanoic acid | + | + |
| Hexanol | | + |
| Hydriodic acid | + | + |
| Hydrobromic acid | | + |
| Hydrochloric acid, 37% | | + |
| Hydrogen peroxide, 35% | + | |
| Isoamyl alcohol | | + |
| Isobutanol | + | + |
| Isopropanol (2-Propanol) | + | + |
| Lactic acid | + | + |
| Methanol | + | + |
| Methyl benzoate | | + |
| Methyl ethyl ketone | | + |
| Methyl propyl ketone | | + |
| Mineral oil (Engine oil) | | + |
| Monochloroacetic acid | | + |

| Reagent | seripettor® | seripettor® pro |
|---|-------------|--------------------|
| Nitric acid, 10% | | + |
| Oxalic acid | + | + |
| Perchloric acid | | + |
| Phenol | | + |
| Phosphoric acid, 85% | | + |
| Piperidine | | + |
| Potassium chloride | + | + |
| Potassium dichromate | + | + |
| Potassium hydroxide | + | + |
| Potassium hydroxide in ethanol | + | + |
| Potassium permanganate | + | + |
| Propionic acid | + | + |
| Propylene glycol (Propanediol) | + | + |
| Pyridine | | + |
| Pyruvic acid | + | + |
| Salicylaldehyde | | + |
| Salicylic acid | + | + |
| Silver acetate | + | + |
| Silver nitrate | + | + |
| Sodium acetate | + | + |
| Sodium chloride | + | + |
| Sodium dichromate | + | + |
| Sodium fluoride | + | + |
| Sodium hydroxide, 30% | + | + |
| Sodium hypochlorite 20% (active chlorine approx. 10%) | | + |
| Sulfuric acid, 10% | + | + |
| Tartaric acid | | + |
| Urea | + | + |
| Zinc chloride, 10% | + | + |
| Zinc sulfate, 10% | + | + |

Note: seripettor® and seripettor® pro are not suitable for HF.

The above recommendations reflect testing completed prior to publication. Always follow instructions in the operating manual of the instrument as well as the reagent manufacturer's specifications. In addition to these chemicals, a variety of organic and inorganic saline solutions (e.g., biological buffers), biological detergents and media for cell culture can be dispensed. Should you require information on chemicals not listed, please feel free to contact BRAND. Status as of: 0614/8

Additional information on the seripettor® and seripettor® pro (operating manual, etc.) can be found at www.brand.de

BRAND®, seripettor® and seripettor® pro are trademarks of BRAND GMBH + CO KG, Germany.

Our technical literature is intended to inform and advise our customers. However, the validity of general empirical values, and of results obtained under test conditions, for specific applications depends on many factors beyond our control. Please appreciate, therefore, that no claims can be derived from our advice. The user is responsible for checking the appropriateness of the product for any particular application.

Subject to technical modification without notice. Errors excepted.

BRAND GMBH + CO KG · P.O. Box 11 55 · 97861 Wertheim · Germany
Phone: +49 9342 808-0 · Fax: +49 9342 808-98000 · E-Mail: info@brand.de · Internet: www.brand.de

