HandyStep®

BRAND GMBH + CO KG
Postfach 11 55
97861 Wertheim · Germany
Tel: +49 9342 808-0
Fax: +49 9342 808-236
E-Mail: info@brand.de
Internet: www.brand.de

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Operating Manual

Please read carefully before use, and follow all operating and safety instructions!

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Safety Instructions

Please read the following carefully!

This instrument may sometimes be used with hazardous materials, operations, and equipment. It is beyond the scope of this manual to address all of the potential safety risks associated with its use in such applications. It is the responsibility of the user of this pipette to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Every user must read and understand this operating manual prior to using the instrument and observe these instructions during use.
2. Observe general instructions for hazard prevention and safety instructions; e.g. wear protective clothing, goggles and gloves. When working with infectious or other hazardous samples, all appropriate regulations and precautions must be followed.
3. Observe all safety precautions provided by reagent manufacturers.

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Purpose

The HandyStep® is a repetitive pipette for quick and simple repetitive dispensing of liquids. After selecting an PD-Tip from a range of sizes and setting the stroke length, volumes from 2 µl to 5 ml can be dispensed with the highest precision and accuracy within the following limitations.

Limitations of use

- +15 °C to +40 °C (59 °F to 104 °F)
(instrument and reagents
- other temperatures on request)
- vapor pressure up to 500 mbar
- viscosity:
 - 20 mPa s with 50 ml PD-Tips
 - 260 mPa s with 5 ml PD-Tips
 - 977 mPa s with 1,25 ml PD-Tips

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Operating exclusions

During proper operation, the dispensed sample only comes in contact with the tip and not with the HandyStep®.

The user has to ensure the compatibility of the instrument with the intended application.

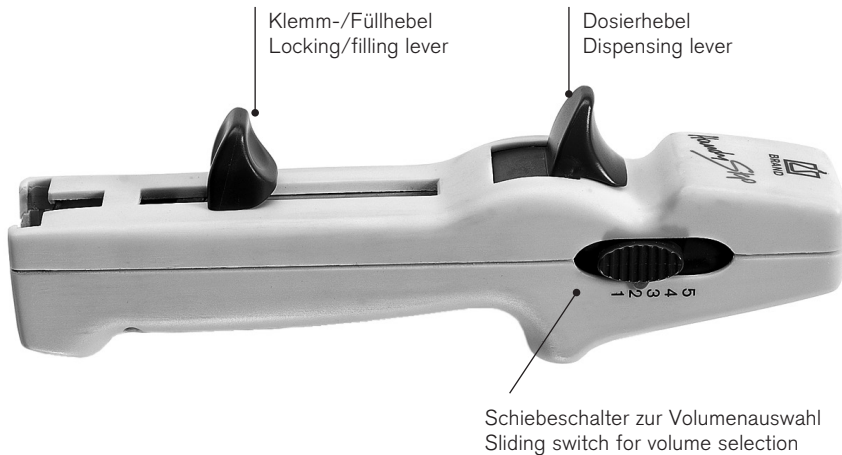
Never dispense samples which attack polypropylene, polyethylene or polycarbonate (housing).

Note:

Instrument and tips are not autoclavable (121 °C / 250 °F). See sterile PD-Tips. PD-Tips are disposable. For disinfection you may use commercially available disinfecting solutions as long as they do not attack PC/PBT.

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

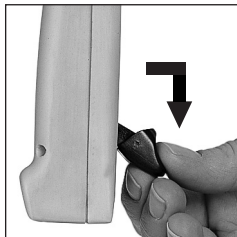
Operation

Mounting the tip

1. Push the locking/filling lever to the lower stop.
2. Pull out the locking/filling lever slightly and swing it out.
3. Insert the PD-Tip.
4. Swing the locking/filling lever back to lock the tip.

Note:

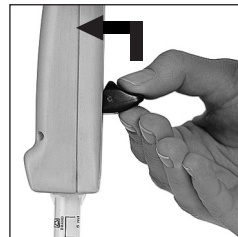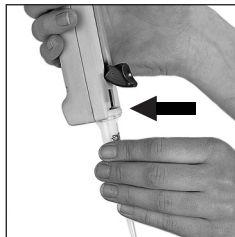To mount 25 ml and 50 ml PD-Tips, use the adapter supplied with the PD-Tips. The reusable adapter twists on and off and can be removed after use.

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Dispense volumes for the HandyStep® are shown in Table 1 below. The table is a matrix of all sizes of PD-Tips and all stroke settings.

The dispensed volume is obtained by multiplying the smallest volume unit (printed on the PD-Tip) by the stroke setting.

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Hinweis:

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Filling the PD-Tip

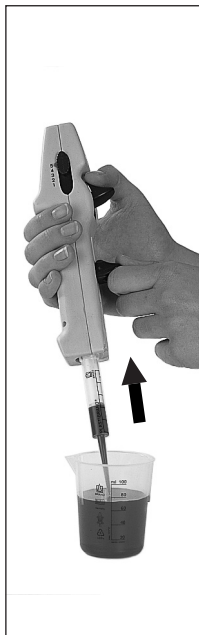
1. Immerse the orifice of the PD-Tip to a depth of 3 - 10 mm into the liquid.
2. Raise the locking/filling lever slowly to prevent cavitation.

Discard the first dispensing step.

Any small air bubble near the piston can be ignored since a locking mechanism prevents the residual stroke from being dispensed.

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Repetitive dispensing

1. Verify the volume setting.
Wipe off excess sample on the PD-Tip around the orifice with a lint-free tissue.
2. Place the orifice of the PD-Tip against the inner wall of the vessel.
3. Dispense the liquid by pressing down the dispensing lever completely each time.
4. Take care to dispense smoothly and evenly.

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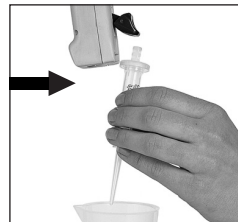
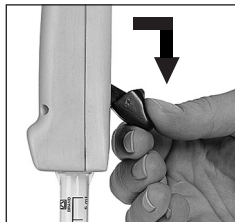
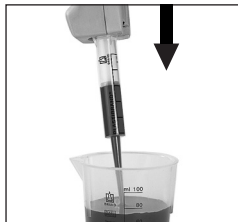
Removing the PD-Tip

1. Hold the HandyStep® over a vessel.
2. Empty the PD-Tip by pushing the locking/filling lever to its lower stop.
3. Pull out the locking/filling lever slightly and swing it out.
4. Remove the PD-Tip.

The adapter required for 25 ml and 50 ml PD-Tip is reusable.

Note:

The tightness of the seal of the PD-Tips is not guaranteed for repeated use with high viscosity media.

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Troubleshooting

Trouble	Possible cause	Corrective action
PD-Tip cannot be inserted	Locking/filling lever is not pushed all the way down and is not swung out The PD-Tip piston is not fully inserted	Push the locking/filling lever to its lower stop and swing it out Push the piston of the PD-Tip fully into the cylinder.
Locking/filling lever cannot be pushed up	Locking/filling lever is not fully pushed in	Make sure locking/filling lever is fully pushed in
HandyStep® dispenses unevenly	Sliding switch is not locked into position	Lock sliding switch into position
PD-Tip drips	PD-Tip seal worn Tip loosely mounted	Replace PD-Tip Push tip on firmly

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Maintenance

The HandyStep® is calibrated at the factory. It may be cleaned with water or isopropyl alcohol.

The HandyStep® contains no user-servicable parts. Never attempt to disassemble the HandyStep®.

Technical data

Measuring results were obtained by using original PD-Tips from BRAND.

Accuracy table HandyStep® with PD-Tips from BRAND, DE-M marking

PD-Tip size	Volume range (µl)	A' ≤ ± %			CV' ≤ %		
		Stroke setting			Stroke setting		
		1**	3**	5**	1**	3**	5**
0.1 ml	2 - 10	4.0	2.4	1.6	6.0	3.0	2.0
0.5 ml	10 - 50	2.5	1.5	1.0	2.5	1.5	1.0
1.0 ml	20 - 100	2.5	1.5	1.0	2.0	1.2	0.8
1.25 ml	25 - 125	2.5	1.4	0.9	2.0	1.1	0.7
2.5 ml	50 - 250	1.8	1.1	0.7	1.5	0.9	0.6
5.0 ml	100 - 500	1.8	1.1	0.7	1.5	0.9	0.7
10.0 ml	200 - 1000	1.8	1.1	0.7	2.0	1.2	0.8
12.5 ml	250 - 1250	1.8	1.1	0.8	3.2	2.0	1.4
25.0 ml	500 - 2500	1.5	0.9	0.6	3.0	1.5	1.0
50.0 ml	1000 - 5000	1.5	0.8	0.5	5.0	1.8	1.2

A' = Accuracy. CV' = Coefficient of variation

** 1 ≅ 2%. 3 ≅ 6%. 5 ≅ 10% of nominal volume

Error limits are based on the maximum volumes for each PD-Tip size obtained at equal temperature (20 °C) of instrument, tip, ambience and dist. H₂O, and with smooth, jerk-free operation.

The HandyStep® can also be used with other compatible tips. The nominal volume is the maximum volume printed on the PD-Tip.

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* incl. 1 adapter, ** sterile/free of endotoxin – single wrapped,

*** PD-Tips 25 ml: non sterile pack of 50 / sterile pack of 25

PD-Tip Set

20 each:

0.5, 1.25, 2.5, 5.0, 12.5 ml

Cat. No. 7023 68

PD-Tips in BIO-CERT®
quality (sterile, free of
endotoxins, DNA, RNase and
ATP) available on request.

Adapter for 25 and 50 ml PD-Tips, PP, autoclavable

Pack of	non-sterile Cat. No.	pack of	sterile** Cat. No.
10	7023 98	5	7023 99

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Return for repair

Important:

For safety reasons, only clean and decontaminated instruments will be accepted for inspection or repair.

- Thoroughly clean and decontaminate the instrument.
- Complete the "Declaration on Absence of Health Hazards" (ask your supplier or the manufacturer for the form. The form can also be downloaded from **www.brand.de**) and send the instrument to the manufacturer or supplier and describe:
 - The nature of the problem
 - What samples have been used

Return shipment has to be effected at the risk and the cost of the sender.

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Disposal

For the disposal of instruments and tips, please observe the relevant national disposal regulations.

Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operation or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original parts have been used.

Subject to technical modification without notice. Errors excepted.