cell-culture™

Absauggerät/Aspiration Controller

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Operating Manual

Please read carefully before use, and follow all operating and safety instructions!

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Contents

	Page
Safety Instructions	4
Applications	6
Operating exclusions	6
Components	7
Before Aspiration	8
Correct Hand Position	8
Aspiration	8
Attaching the Pipette/Pipette Tip	8
Aspiration of Liquid	9
After Aspiration	10
Removing the Pipette/Pipette Tip	10
Releasing the Suction Tube	10
Cleaning/Autoclaving	11
Storage	12
Instructions for Repair	13
Replacing Adapter and Suction Tube	13
Trouble Shooting	15
Return for Repair	16
Contact addresses	18
Warranty	19
Disposal	19
Ordering Data	20

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Please read carefully before use!

Safety Instructions

This instrument may sometimes be used with hazardous materials, operations, and equipment. It is beyond the scope of this manual to address all of the potential safety risks associated with its use in such applications. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Every user must read and understand this operating manual prior to using the instrument and observe these instructions during use.
2. Follow general instructions for hazard prevention and safety instructions; e.g., wear protective clothing, eye protection and gloves. When working with infectious or other hazardous samples, all appropriate regulations and precautions must be followed.
3. Observe the reagent manufacturers' information.

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Applications

This instrument allows aspiration of excess liquids.

The instrument is applied with 3 different adapters to hold pasteur pipettes, capillaries, or pipette tips (100 µl). For easy fitting of pipette/pipette tips the adapter can be rotated.

For aspiration of liquids, the instrument must be connected to a collection bottle and a vacuum pump.

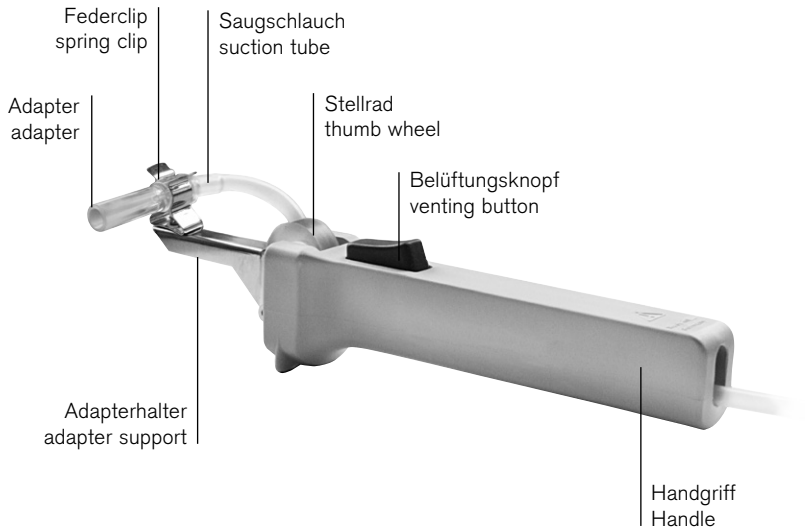
Operating exclusions

The user has to ensure the compatibility of the instrument with the intended application.

Vapours of liquids may come into contact with the adapter (SI, PVC or PP) and suction tube (SI). Therefore never use the instrument for liquids which attack PVC, PP or SI.

The cell-culture™ Aspiration Controller (K) is not autoclavable. The included tube and the adapters are autoclavable.

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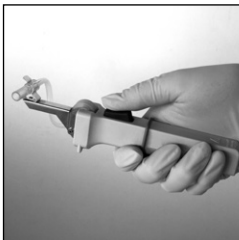
Before Aspiration

Correct hand position

Hold the instrument horizontally as shown.

Push back the O-ring from the exhaust button.

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Attaching the Pipette/Pipette Tip

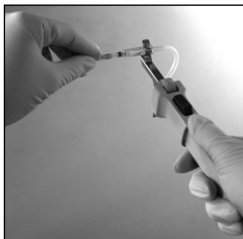
Warning!

Never insert filled pipettes due to risk of contamination and infection.

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
Note:

- Hold the pasteur/capillary pipette at the short end to avoid glass breakage when fitting the pipette.
- Carefully insert the pasteur/capillary pipette into the adapter until a slight resistance can be felt. Never use force.
- If the pasteur/capillary pipette breaks: Remove glass splinters with forceps. never use bare hands. Check adapter for damage and, if necessary, replace adapter. (► "Instructions for repair", p. 13).

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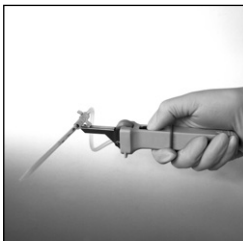Aspiration of Liquid

- Press the control button until the liquid has been aspirated.
- Release the control button to stop aspiration.

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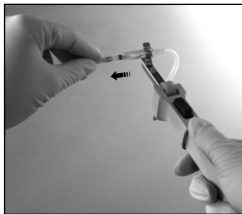
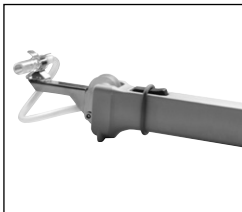
Removing the Pipette/Pipette Tip

Warning!

Never touch the contaminated area of the pipette/pipette tip. Always wear protective gloves.

- Gently grip the pipette just below the adapter.
- Carefully pull it out of the adapter and dispose of properly, observing all applicable laws of regulations.

Releasing the Suction Tube

After aspiration always release the suction tube by pushing the O-ring on the exhaust button.

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Cleaning/Autoclaving

Occasionally wipe the instrument with a damp cloth especially if it comes in contact with sample. When properly used, the instrument is maintenance-free.

Warning!

Risk of contamination and infection. Wear protective gloves and eye protection.

- Remove adapter and suction tube and put all parts in a cleansing or disinfection bath.
- Allow for complete drying.
- If necessary, autoclave adapter and suction tube.

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Adapter and suction tube are autoclavable at 121 °C (250 °F), 2 bar absolute (30 psi), according to DIN EN 285. It is the user's responsibility to ensure effective autoclaving.

- Reassemble instrument (► "Instructions for repair", page 13).
- Check unit for correct function of all operating components.

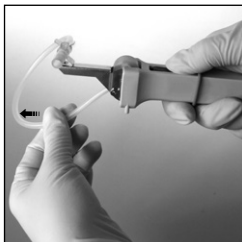
Storing

Only store instrument with completely released suction tube (► "Release suction tube", page 10).

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Instructions for Repair

Replacing Adapter and Suction Tube

- Gently remove the adapter from the adapter support.
- While pressing the control button the suction tube can be pulled out.
- While pressing the control button insert the new suction tube with adapter between thumb wheel and roller.
- Clip the adapter into the adapter support.
Gently exert counterpressure at the bottom of the spring clip to avoid damage of the adapter support.

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Trouble Shooting

Trouble	Possible Cause	Action to be taken
Broken pipette in the adapter.	The pipette was damaged/forced into the adapter.	Remove the glass splinters by means of forceps, replace adapter and suction tube, if necessary, ► page 13.
The instrument does not aspirate liquid/pipette drips.	Adapter or suction tube damaged.	Replace adapter and suction tube, ► page 13.
Pipette not held properly in the adapter.	Adapter is greasy.	Clean and disinfect adapter and suction tube, and allow to dry completely. Replace if necessary, ► page 11.

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Return for Repair

Attention:

Transporting of hazardous materials without a permit is a violation of federal law.

- Clean and decontaminate the instrument carefully.
- It is essential always to include an exact description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Shipment is at the risk and the cost of the sender.

Outside the U.S. and Canada:

- Complete the "Declaration on Absence of Health Hazards" and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from www.brand.de.

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In the U.S. and Canada:

- Contact BrandTech Scientific, Inc. and obtain authorization for the return **before** sending your instrument for service.
- Return only cleaned and decontaminated instruments, with the Return Authorization Number prominently displayed on the outside of the package to the address provided with the Return Authorization Number.

Kontaktadressen/Contact addresses

BRAND GMBH + CO KG

Otto-Schott-Straße 25
97877 Wertheim (Germany)
Tel.: +49 9342 808-0
Fax: +49 9342 808-98000
E-Mail: info@brand.de
www.brand.de

India:

BRAND Scientific Equipment Pvt. Ltd.
303, 3rd Floor, 'C' Wing, Delphi
Hiranandani Business Park, Powai
Mumbai - 400 076 (India)
Tel.: +91 22 42957790
Fax: +91 22 42957791
E-Mail: info@brand.co.in
www.brand.co.in

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Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operation or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual.

We are not liable for damage resulting from any actions not described in the operating manual or if non-original spare parts or components have been used.

U.S. and Canada:
Information for warranty please see www.brandtech.com.

Disposal

For the disposal of instruments, please observe the relevant national disposal regulations.

Subject to technical modification without notice. Errors excepted.

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Best.-Nr. 259 62

Ordering Data

cell-culture (Aspiration Controller),

complete with suction tube
and 3 adapters

Cat. No. 259 50

Spare parts

Adapter (SI, PVC) for pasteur pipettes,

pack of 1

Cat. No. 259 60

Adapter (PVC) for capillaries,

pack of 1

Cat. No. 259 33

Adapter (PP) for pipette tips (100 µl),

pack of 1

Cat. No. 259 61

Suction tube (SI), 2 m,

pack of 1

Cat. No. 259 62