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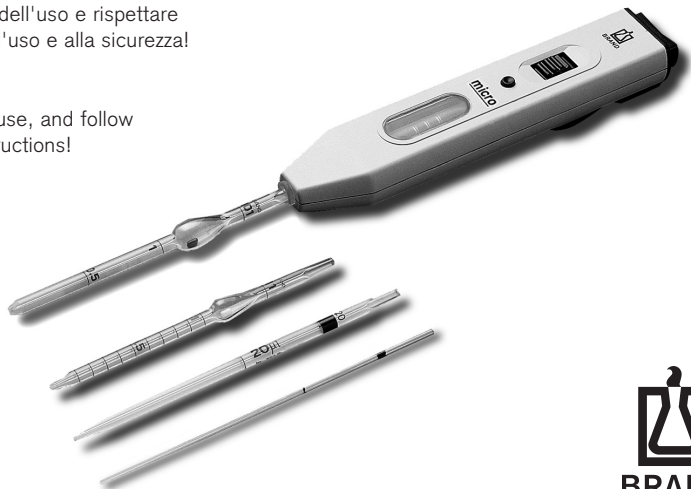
Operating Manual

Please read carefully before use, and follow all operations and safety instructions!

BRAND GMBH + CO KG
Otto-Schott-Str. 25
97877 Wertheim
Germany
Tel.: +49 9342 808-0
Fax: +49 9342 808-236
E-Mail: info@brand.de
Internet: www.brand.de

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Read carefully before use!

Safety Instructions

This instrument may sometimes be used with hazardous materials, operations, and equipment. It is beyond the scope of this manual to address all of the potential safety risks associated with its use in such applications. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

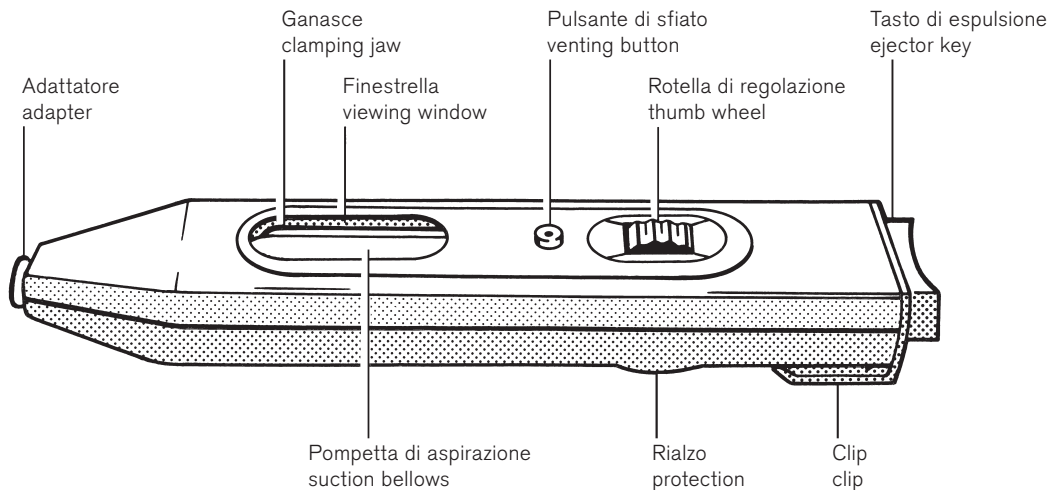
1. Every user must read and understand this operating manual before operation.
2. Follow general instructions for hazard prevention and safety instructions; e.g., wear protective clothing, eye protection and gloves.
3. Observe all specifications provided by reagent manufacturers.
4. Use the instrument only for pipetting liquids, with strict regard to the defined operating limitations. Observe operating exclusions (s. page 8! If in doubt, contact the manufacturer or supplier.

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5. Always use the instrument in such a way that neither the user nor any other person is endangered. Avoid splashes. Only use suitable vessels.
6. Only fit empty pipettes to avoid the risk of contamination and infection.
7. Take care that the liquid does not penetrate into the instrument.
8. Never use force on the instrument!
9. Use only original manufacturer's accessories and spare parts. Do not attempt to make any technical alterations. Do not dismantle the instrument any further than is described in the operating manual!
10. Always check the instrument for visible damage before use. If there is a sign of a potential malfunction, immediately stop pipetting. Consult the 'Troubleshooting' section of this manual (see page 28), and contact the manufacturer if needed.

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Application

This instrument provides for pipetting liquids, and reduces the risk of infection, when pipetting liquids with **ring-mark disposable micropipettes, blood diluting pipettes, and other pipettes of a shaft diameter < 5.5 mm and a volume of up to 1 ml.**

When the instrument is correctly used, the pipetted liquid comes only into contact with the pipette.

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Operating Limitations

Attention:

This instrument must not be used with end-to-end micropipettes.

Vapors of liquids may penetrate into the suction system. Therefore never use the instrument for liquids which attack silicone.

Note:

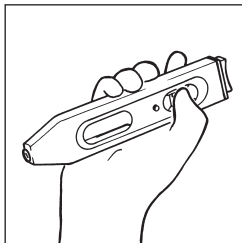
This instrument is autoclavable at 121 °C (250 °F), 2 bar absolute according to DIN EN 285. It is the user's responsibility to ensure effective autoclaving.

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Before Pipetting

Correct hand position

Hold the instrument horizontally as shown in figure. The projection at the back of the housing is placed between index and middle fingers.

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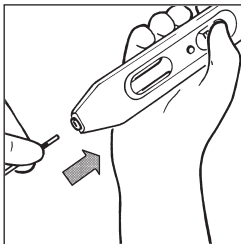
Attaching the pipette

Warning:

Never fit filled pipettes: Risk of contamination and infection.

- Hold the pipette at the color code (short end up) to avoid glass breakage when fitting the pipette.
- Carefully insert the pipette into the adapter until a slight resistance can be felt. Never use force: Risk of damaging the instrument and pipette, hence a risk of injuries.
- If the pipette breaks: Remove the glass splinters by means of forceps. Inspect the adapter for damage. If necessary, replace it (➤ "Instructions for repair", page 20).

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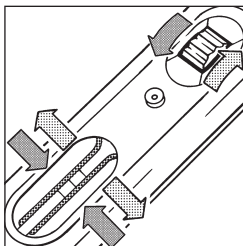
Adjusting the suction volume

Micropipettes $\leq 5 \mu\text{l}$

Turn the thumb wheel towards the ejector key until the clamping jaws are completely open.

Micropipettes $> 5 \mu\text{l}$ up to $100 \mu\text{l}$

Turn the thumb wheel towards the window until the clamping jaws compress the suction bellows to about half of the volume (see figure).

Pipettes $> 100 \mu\text{l}$

Turn the thumb wheel towards the window until a slight resistance can be felt. The clamping jaws completely compress the suction bellows.

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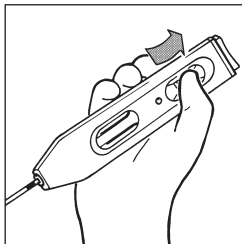
Drawing up liquid

Attention:

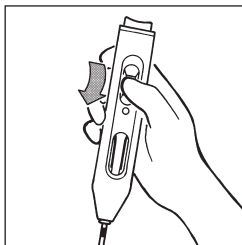
Always ensure that the liquid does not contaminate the suction system. If it is contaminated, it must immediately be cleaned or disinfected or, if necessary, be replaced.

- Hold the instrument at an angle.
- Immerse the pipette into the liquid.
- **Slowly** turn the thumb wheel towards the ejector key until the liquid has exactly reached the appropriate mark. (The micropipettes $\leq 5 \mu\text{l}$ fill up automatically. Therefore turn the thumb wheel slowly towards the viewing window to stop the suction process in time at the appropriate mark.)
- Take the pipette out of the liquid.
- Carefully wipe the outside of the pipette with a lint-free cloth or swab.

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Discharging liquid

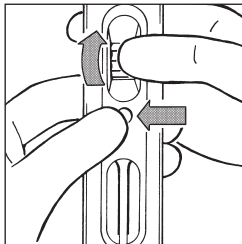
Pipettes calibrated to contain "In"

- Immerse the pipette into the diluting agent.
- Turn the thumb wheel towards the viewing window until the liquid has been discharged.
- Rinse the pipette by drawing in and discharging the diluting agent: Carefully draw in the diluting agent up to the correct level of the liquid to be measured. Repeat this procedure at least three times.
- Lightly wipe the pipette against the wall of the vessel.

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
Note:

If accidentally the suction bellows have been compressed too much when discharging the pipette and a small residue of liquid remains in the pipette after rinsing:

- Press the venting button, and simultaneously turn the thumb wheel slightly towards the ejector key.
- Release the venting button, and turn the thumb wheel towards the window until the residue of liquid has been discharged.

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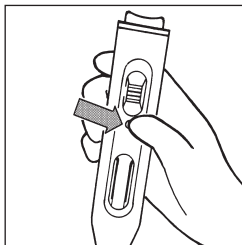Pipettes calibrated to deliver "Ex"

- Hold the instrument upright, and place the pipette tip against the wall of a vessel (inclined position of the vessel).
- Press the venting button until the liquid has run out.
- As soon as the meniscus comes to a standstill in the pipette tip, the waiting time (if applicable) starts.
- Then take the pipette out of the vessel by wiping it against the wall of the vessel. During this procedure the residue of liquid partly runs out.

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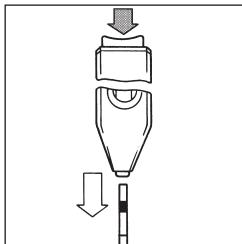
Removing the pipette

Warning:

Do not touch the contaminated area of the pipette. Wear protective gloves.

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
Disposable micropipettes up to 50 μ l

- Hold the instrument upright over a vessel (pipette down).
- Press the ejector key.
- Dispose the removed pipette according to the relevant provisions.

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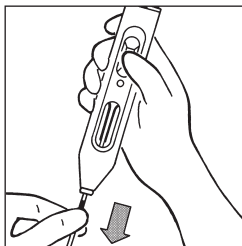Disposable micropipettes > 50 µl and all reusable pipettes

These pipettes cannot be ejected.

- Grip the pipette just below the adapter.
- Carefully pull it out of the adapter.
- Dispose the micropipettes according to the relevant provisions.

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Releasing the suction bellows

After pipetting always release the suction bellows. For this purpose turn the thumb wheel towards the ejector key until the clamping jaws are completely open.

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Cleaning / Autoclaving

Occasionally wipe the instrument with a damp cloth. When properly used, the instrument is maintenance free.

If liquid has contaminated the suction system, these parts must immediately be cleaned or disinfected or, if necessary, be replaced (➤ "Instructions for repair", page 20).

Warning!

Risk of contamination and infection.
Wear protective gloves and eye protection.

- Remove the suction system and put all parts in a cleansing or disinfecting bath.
- Then allow to dry completely.
- Autoclave the instrument, if necessary.

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Note:

The instrument is autoclavable at 121 °C (250 °F), 2 bar absolute (30 psi) with a holding time of at least 15 minutes according to DIN EN 285. It is the user's responsibility to ensure effective autoclaving.

- Reassemble the instrument
(➤ "Instructions for repair", page 20).
- Check that all operating elements work properly.

Storing

When storing the instrument, the suction bellows must be completely released (➤ "Releasing the suction bellows", page 17).

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Instructions for Repair

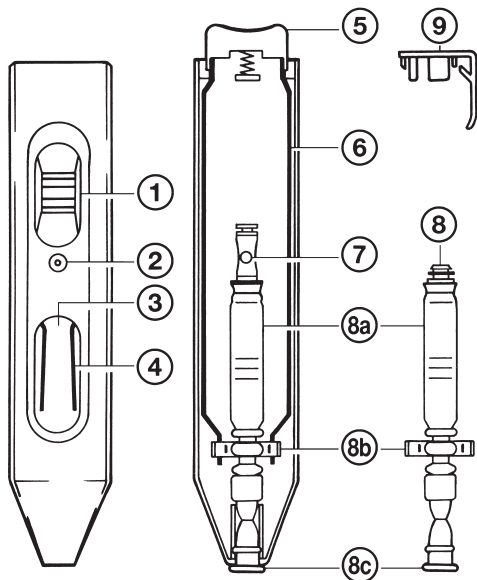
Replacing the suction system / Dismantling the instrument

Warning!

Do not touch the contaminated area of the pipette. Wear eye protection and protective gloves.

Never remove glass fragments of broken pipettes by hand but always use a forcep.

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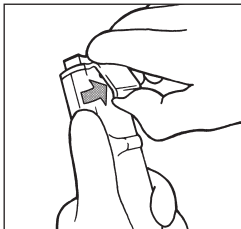
Back part of the instrument

- 5 ejector key
- 6 side rails
- 7 valve
- 8 suction system
- 8a suction bellows
- 8b plastic ridge
- 8c adapter
- 9 clip

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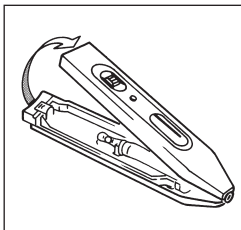
Dismantling the instrument

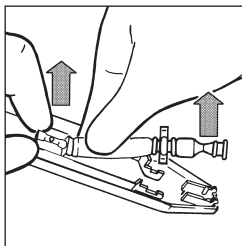
(➤ "Components", page 21)

- Turn the thumb wheel ① towards the ejector key until a slight resistance can be felt. Then turn the thumb wheel with 2 rotary finger motions towards the window.

- Carefully lift the clip ⑨ until the pin is released from the recess. Press the clip with its cover-ring upwards and off.

- Open the instrument (➤ see figure), and put the front part aside.

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
- Remove the suction system ⑧ with the valve ⑦ from the support.
- Pull off the suction bellows ⑧a from the valve ⑦.
- If the suction system is just dirty: Put it into a cleansing or disinfecting bath and, if necessary, autoclave it (121 °C). It is the user's responsibility to ensure effective autoclaving.
- If the suction system remains dirty, or if it is damaged: dispose and replace it by a new one.

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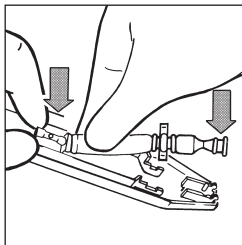
Reassembling the instrument

(➤ "Components", page 21)


- Insert the cleaned or new suction system:
 - Slide the suction system ⑧ on the valve ⑦ and insert it into the support.
 - Insert the plastic ridge ⑧b into the two grooves of the side rails ⑥.
 - Press the adapter ⑧c into the back part of the instrument.

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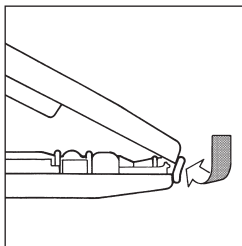- Place the front part of the instrument behind the outer bulge of the adapter, and attach it to the hooks of the back part of the instrument.

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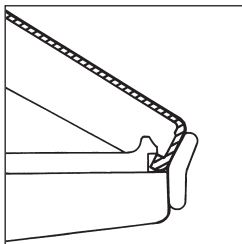- Press both parts of the instrument slowly and uniformly together.
- Hold the instrument loosely, and shortly press down the ejector key.
- Slide on the clip, and make it snap in.
- Check that all operating elements work properly.

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Return for Repair

Attention:

For safety reasons, instruments returned for repair must be clean and decontaminated.

- Therefore: Clean and decontaminate the instrument carefully.
- Complete the "Declaration on Absence of Health Hazards" (ask your supplier or manufacturer for the form. The form can also be downloaded from **www.brand.de**) and send the instrument to the manufacturer or supplier and describe:
 - the nature of the problem
 - the pipetted liquids.

Return shipment has to be effected at the risk and the cost of the sender.

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Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original parts have been used.

Subject to technical modification without notice. Errors excepted.

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Troubleshooting

Trouble	Possible Cause	Corrective action
Broken pipette in the adapter	Pipette was damaged/ forced into the adapter	Remove the glass splinters by means of forceps; if necessary, replace adapter (➤ p. 10)
The instrument does not draw in liquid/pipette drips	Suction system damaged	Replace the suction system (➤ p. 20-24)
Liquid in suction system	Thumb wheel was turned too quickly	Clean or disinfect the suction system or, if necessary, replace it (➤ p. 18, 19)
Ejector key does not work	Ejector key not properly positioned	Remove the clip. Hold the instrument loosely, and quickly press down the ejector key. Slide on the clip (➤ p. 22, 23, 24)
Pipette is ejected from the adapter	Adapter is wet or greasy	Clean or disinfect suction system and allow to dry completely. If necessary, replace suction system (➤ p. 18-19, 20-24)

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