# micro-classic (Controlador de Pipetagem/Pipette Controller)

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## Operating Manual

Please read carefully before use, and follow all operating and safety instructions!

BRAND GMBH + CO KG  
Otto-Schott-Str. 25  
97861 Wertheim  
Germany  
Tel.: +49 9342 808-0  
Fax: +49 9342 808-98000  
E-Mail: [info@brand.de](mailto:info@brand.de)

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## Contents

	Page
Safety Instructions	4
Components	6
Application	7
Operating Limitations	7
Before Pipetting	8
Correct Hand Position	8
Pipetting	8
Attaching the Pipette	8
Drawing up Liquid	9
Discharging Liquid	11
After Pipetting	13
Removing the Pipette	13
Releasing the Suction Tube	13
Cleaning/Autoclaving	14
Storage	15
Replacing Adapter and Suction Tube	16
Trouble Shooting	17
Return for Repair	19
Warranty	20
Ordering Data	21

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## Safety Instructions

This instrument may sometimes be used with hazardous materials, operations, and equipment. It is beyond the scope of this manual to address all of the potential safety risks associated with its use in such applications. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

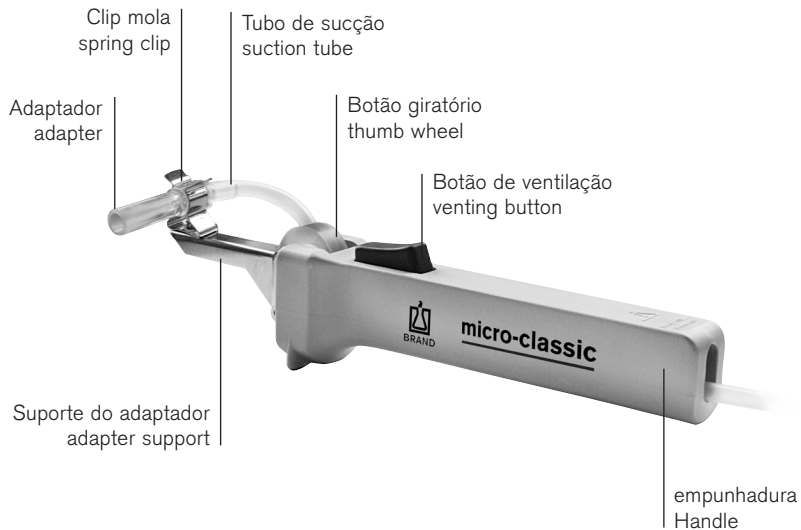
### **Please read the following carefully!**

1. Every user must read and understand this operating manual before operation.
2. Follow general instructions for hazard prevention and safety instructions; e.g., wear protective clothing, eye protection and gloves.
3. Observe all specifications provided by reagent manufacturers.
4. Use the instrument only for pipetting liquids, with strict regard to the defined operating limitations. If in doubt, contact the manufacturer or supplier.

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## Application

This instrument provides for pipetting liquids, and reduces the risk of infection, when pipetting liquids with **ring-mark disposable micropipettes, blood diluting pipettes, and other pipettes of a shaft diameter < 5.5 mm and a volume of up to 1 ml.**

When the instrument is correctly used, the pipetted liquid comes only into contact with the pipette.

## Operating Limitations

### Attention!

This instrument must not be used with end-to-end micropipettes.

Vapours of liquids may penetrate into the adapter (PVC) and into the suction tube (SI). Therefore never use the instrument with liquids that are incompatible with PVC or SI.

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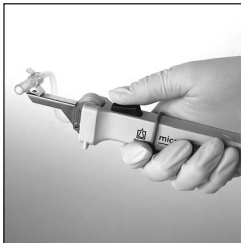
## Before Pipetting

### Correct Hand Position

Hold the instrument horizontally as illustrated.

Push back the O-Ring from the venting button.

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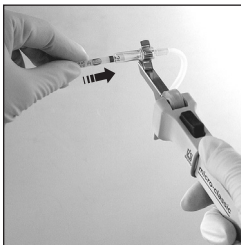
### Attaching the Pipette

#### **Warning!**

Never insert filled pipettes:  
Risk of contamination and infection.  
Wear protective gloves.

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- Hold the pipette at the colour code (short end up) to avoid glass breakage when fitting the pipette.
- Carefully insert the pipette into the adapter until a slight resistance can be felt. Never use force: Risk of damaging the adapter and pipette, hence a risk of injury.
- If the pipette breaks:  
Remove glass splinters by means of forceps. Check adapter for damages and, if necessary, replace adapter. (► Instructions for repair, page 16).

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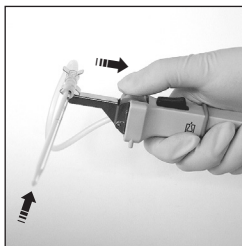
## Drawing up Liquid

Allow the tube a length of approx. 100 mm between adapter and thumb wheel to keep the dead volume low.

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- Immerse the pipette into the liquid.
- Slowly turn the thumb wheel towards the ejector key until the liquid has exactly reached the appropriate mark.

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### Attention!

Always ensure that the liquid does not contaminate adapter and suction tube. If they are contaminated, they must immediately be cleaned or disinfected or, if necessary, be replaced.

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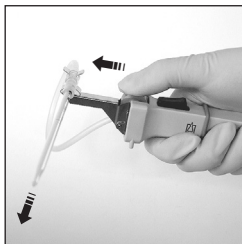
- Take the pipette out of the liquid.
- Carefully wipe the outside of the pipette with a lint-free cloth or swab. Take care not to touch the capillary orifice to avoid liquid being sucked out.

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
## Discharging Liquid

### Pipettes calibrated to contain "In"

- Immerse the pipette into the diluting agent.
- Turn the thumb wheel towards the adapter until the liquid has been discharged.
- Rinse the pipette by drawing in and discharging the diluting agent: Carefully draw in the diluting agent up to the correct level of the liquid to be measured. Repeat this procedure at least three times.
- Lightly wipe the pipette against the wall of the vessel.

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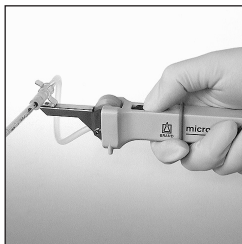### Pipettes calibrated to deliver "Ex"

- Hold the instrument upright, and place the pipette tip against the wall of a vessel (inclined position of the vessel).
- Press the venting button until the liquid has run out.
- As soon as the meniscus comes to a standstill in the pipette tip, the waiting time (if applicable) starts.
- Then take the pipette out of the vessel by wiping it against the wall of the vessel. During this procedure the residue of liquid partly runs out.

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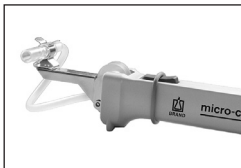
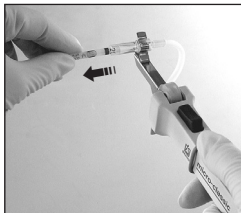
### Removing the Pipette

#### Warning!

Do not touch the contaminated area of the pipette. Wear protective gloves.

- Grip the pipette just below the adapter.
- Carefully pull it out of the adapter.
- Dispose the micropipettes according to the relevant provisions.

### Releasing the Suction Tube

After pipetting always release the suction tube by pushing the O-ring on the venting button.

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## Cleaning/Autoclaving

Occasionally wipe the instrument with a damp cloth. When properly used, the instrument is maintenance-free.

If liquid has contaminated adapter and suction tube, these parts must immediately be cleaned or disinfected and, if necessary, be replaced (► Instructions for repair, page 16).

### **Warning!**

Risk of contamination and infection. Wear protective gloves and eye protection.

- Remove adapter and suction tube and put all parts in a cleansing or disinfection bath.
- Allow for complete drying.
- If necessary, autoclave adapter and suction tube.

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Adapter and suction tube are autoclavable at 121 °C (250 °F), 2 bar absolute (30 psi) with a holding time of at least 15 minutes according to DIN EN 285. It is the user's responsibility to ensure effective autoclaving.

- Reassemble instrument (► Instructions for repair, page 16).
- Check correct function of all operating elements.

## Storage

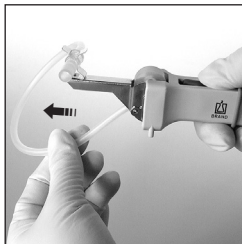
Only store instrument with completely released suction tube (► "Release suction tube", page 13).

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## Replacing Adapter and Suction Tube

- Pull off the adapter from the adapter support.
- Upon pressing the venting button the suction tube can be pull out forward.
- Insert the new suction tube with adapter between thumb wheel and roller while pressing the venting button.
- Clip the adapter into the adapter support.

Simultaneously exert counterpressure at the bottom of the spring clip to avoid damage of the adapter support.

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## Trouble Shooting

<b>Trouble</b>	<b>Possible Cause</b>	<b>Action to be taken</b>
Broken pipette in the adapter.	The pipette was damaged/forced into the adapter.	Remove the glass splinters by means of forceps, replace adapter and suction tube, if necessary, ► page 16.
The instrument does not draw in liquid/pipette drips.	Adapter or suction tube damaged.	Replace adapter and suction tube, ► page 16.
Liquid in the adapter and the suction tube.	The thumb wheel was turned too quickly.	Clean or disinfect adapter and/or suction tube, if necessary, replace, ► page 14.
Pipette not held properly in the adapter.	Adapter is wet or greasy.	Clean or disinfect adapter and suction tube, and allow to dry completely. If necessary, replace, ► page 14.

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## Return for Repair

### Attention!

For safety reasons only cleaned/decontaminated instruments can be checked/repared.

- Therefore: Clean and decontaminate the instrument carefully!
- Complete the "Declaration on Absence of Health Hazards" (ask your supplier or the manufacturer for the form. The form can also be downloaded on **www.brand.de**) and send the instrument to the manufacturer or supplier and describe:
  - the nature of the problem
  - the pipetted liquids

Return shipment has to be effected at the risk and cost of the sender.

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## **Warranty**

We shall not be liable for the consequences of improper handling, use, servicing, operation or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original spare parts or components have been used.

Subject to technical modification without notice.  
Errors excepted.

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